

# NIS-Elements

## Manual

Publication date 08.07.2020

v. 5.30.00

No part of this publication may be reproduced or transmitted except with the written permission of Laboratory Imaging s.r.o. Information within this publication is subject to change without notice. Changes, technical inaccuracies and typographical errors will be corrected in subsequent editions.


Laboratory Imaging s.r.o.  
Za Drahou 171/17  
102 00 Praha 10  
Czech Republic



---

# Table of Contents

1. Startup .....	1
1.1. Command Line Startup Options .....	1
1.2. Fixed Grabber Startup .....	2
2. Installation and Settings .....	5
2.1. Installation and Updates .....	5
2.1.1. The Installation DVD-ROM Content .....	5
2.1.2. NIS-Elements Installation Steps .....	5
2.1.3. Command Line Installation Options .....	9
2.1.4. Additional Module/Device Installation .....	9
2.1.5. Sample Database Installation .....	10
2.1.6. Software Copy Protection .....	10
2.1.7. Device Updates .....	10
2.1.8. Fixes .....	11
2.1.9. Installing the Database Module on 64-bit Systems .....	11
2.2. Additional Modules Available in NIS-Elements D .....	11
2.2.1. Acquisition and Analysis Modules .....	11
2.2.2. Industrial Modules .....	11
2.2.3. Device Modules .....	12
2.2.4. Other .....	13
2.3. User Rights .....	13
2.3.1. NIS-Elements authentication .....	14
2.3.2. Creating a Shared Layout .....	14
2.3.3. Closing Another Instance of NIS-Elements .....	14
2.3.4. User Rights Options .....	15
2.3.5. Users Tab Options .....	16
2.3.6. Groups Tab Options .....	17
2.3.7. Privileges Tab Options .....	18
2.4. NIS-Elements Preferences .....	19
2.4.1. Adjusting Program Preferences .....	19
2.4.2. General .....	19
2.5. Hardware licences .....	21
2.5.1. What is a hardware licence? .....	21
2.5.2. How do I find out the content of my USB key? .....	22
2.5.3. Is there an update for my software? .....	22
2.5.4. What is SUA? .....	22
3. User Interface .....	25
3.1. Advanced Layout .....	25
3.1.1. Main Window Components .....	25
3.1.2. Image Window .....	28
3.1.3. Arranging User Interface .....	30
3.2. Modifying Layouts .....	31
3.2.1. Layout Manager .....	31
3.3. Modifying Toolbars .....	33
3.4. Modifying Menus .....	35
3.5. Running a Macro Upon Layout Change .....	36
3.6. Appearance Options .....	36
3.7. Simplified User Interface .....	37
3.7.1. Behavior .....	37
3.7.2. Toolbar Buttons .....	38
3.7.3. Camera Control Panel .....	38

3.7.4. Save Control Panel .....	38
3.7.5. Preview Control Panel .....	39
4. Cameras & Devices .....	41
4.1. Basic Workflows .....	41
4.1.1. Startup Hardware Configuration Selection .....	41
4.1.2. Dual, Triple, Quad Camera Modes .....	41
4.1.3. Using a <i>Multiview</i> Device .....	41
4.1.4. Optical Configurations .....	42
4.1.5. Objectives .....	44
4.1.6. Connecting a Device to NIS-Elements .....	49
4.1.7. What are “Logical Devices”? .....	50
4.1.8. XY Stages and Z Drives Tips .....	52
4.1.9. Camera Settings .....	53
4.1.10. Controlling Illumination Devices .....	59
4.2. Visual Device Manager .....	59
4.2.1. Connecting Cameras .....	61
4.2.2. Connecting Devices .....	62
4.2.3. Lightpath Scheme  .....	63
4.3. Legacy Cameras and Devices .....	65
5. Image Acquisition .....	67
5.1. Introduction to Image Acquisition .....	67
5.2. Camera ROI .....	68
5.3. Shading Correction .....	68
5.3.1. Simple Shading Correction .....	69
5.3.2. Shading Correction per Optical Configuration .....	70
5.3.3. Removing Dust from Camera Chip Using Shading Correction .....	70
5.3.4. Shading Correction Panel .....	71
5.4. About ND Acquisition .....	73
5.4.1. Introduction to ND Acquisition .....	73
5.4.2. Common ND Experiment Options .....	73
5.4.3. ND Sequence Options .....	74
5.5. Time-lapse Acquisition .....	75
5.5.1. Special Options .....	75
5.6. Multi-point Acquisition .....	76
5.6.1. Point by Point (Manual) Multi-Point .....	76
5.6.2. Well Plate (Rectangular) Multi-Point .....	77
5.6.3. Large Image (Covering) Multi-Point .....	77
5.6.4. Random Multi-Point .....	77
5.6.5. To Change a Single Z Coordinate .....	78
5.6.6. Special Options .....	78
5.7. Z-series Acquisition .....	79
6. Displaying Images .....	83
6.1. Open, Save, Close Images .....	83
6.1.1. Opening Image Files .....	83
6.1.2. Saving Image Files .....	84
6.1.3. Closing Images .....	85
6.1.4. Supported Image Formats .....	85
6.1.5. Floating Point Images .....	86
6.2. Image Layers .....	86
6.2.1. Introduction to Image Layers .....	86
6.2.2. Image Types .....	87
6.2.3. Channel Coloring .....	88

6.2.4. Assigning Colors to Channels .....	89
6.2.5. Copying Channels by Drag and Drop .....	89
6.3. ND Dimensions .....	90
6.3.1. Navigation in ND2 Files .....	90
6.3.2. ND Views .....	92
6.4. Large Images .....	94
6.4.1. Opening Files in Progressive Mode .....	94
6.4.2. Splitting Large Images .....	94
6.4.3. Locating XY Positions Between Images .....	95
6.4.4. Methods Used for Stitching Large Images .....	95
6.5. LUTs - Non-destructive Image Enhancement .....	96
6.5.1. Introduction to Look-Up Tables (LUTs) .....	96
6.5.2. LUTs on RGB Images .....	97
6.5.3. LUTs on Monochromatic Images .....	99
6.5.4. LUTs on Multichannel Images .....	101
6.6. Organizer .....	101
6.6.1. About Organizer .....	101
6.6.2. Image Filter .....	102
6.6.3. Operations with Images within Organizer .....	103
6.6.4. Resizing the Organizer Panes .....	105
6.6.5. Reporting .....	105
7. Image Analysis .....	107
7.1. Preparation .....	107
7.1.1. Histogram .....	107
7.1.2. Image processing .....	109
7.2. Image Segmentation .....	110
7.2.1. Thresholding .....	110
7.3. Binary Layers .....	113
7.3.1. Introduction to Binary Layers .....	113
7.3.2. Managing Binary Layers .....	113
7.3.3. Binary Editor .....	115
7.3.4. Mathematical Morphology Basics .....	116
7.4. Regions of Interest - ROIs .....	120
7.4.1. Introduction to ROIs .....	120
8. Measurement .....	123
8.1. Basic Workflows .....	123
8.1.1. Calibration .....	123
8.1.2. Rough Measurement .....	124
8.1.3. Object Count .....	125
8.1.4. Measurement Options .....	130
8.1.5. Measurement on Graph .....	131
8.1.6. Pixel Classifier .....	132
8.1.7. Exporting Results .....	134
8.2. Manual Measurement .....	136
8.2.1. Basic Workflow .....	136
8.2.2. Working with the Measurement and Annotation Objects .....	137
8.2.3. Measurement Tools .....	138
8.3. Automated Measurement .....	142
8.4. Measurement Features .....	144
9. Creating Reports .....	159
9.1. Report Generator .....	159
9.2. Report Objects .....	159
9.2.1. Aligning Objects .....	160

---

9.2.2. Dynamic Data .....	161
9.3. Report Templates .....	161
10. Macros .....	163
10.1. Creating Macros .....	163
10.2. Running a Macro .....	165
10.3. Interactive Advanced Macro (API) .....	165
10.4. Macro Language Syntax .....	166
10.4.1. Variable types .....	166
10.4.2. Structures and Unions .....	166
10.4.3. Arrays .....	166
10.4.4. Local and Global Variables .....	166
10.4.5. Statements .....	167
10.4.6. Directives .....	168
10.4.7. Operators .....	170
10.4.8. Priority of Operators (!) .....	171
10.4.9. C-like functions .....	172
10.5. Controlling Cameras by Macro .....	173
10.6. Macro Preferences .....	173
11. Movies .....	175
11.1. Capturing AVI Movie .....	175
11.2. Save ND2 as AVI .....	175
11.3. Creating a Movie from EDF Surface View .....	175
11.3.1. Creating a Movie Step by Step .....	175
11.3.2. Movie Maker Controls .....	176
11.3.3. Directors Mode .....	177
11.4. About Video Compression .....	177
Index .....	179

# 1. Startup

## 1.1 Command Line Startup Options

See also 2.1.3 Command Line Installation Options.

When starting *NIS-Elements* from the command line (or when editing the desktop shortcut properties), you may append some switches with parameters to the main command and thereby modify the startup behavior.

---

### Note

*Some switches do not have parameters.*

*Example 1.1. Command line switches syntax*

"*NIS-Elements* executable" *switch#1 parameter#1 switch#2 parameter#2* etc.

For example, to run *NIS-Elements* and open the *starting\_image.jp2* at straight away, run.:

```
"c:\Program Files\NIS-Elements D\nis_d.exe" -f "C:\Images\starting_image.jp2"
```

### **Startup Switches**

**-?** Displays a help screen with the description of switches (basically this page).

**-c "Command"** The application runs the specified macro *Command*.

**-cam "Camera Name"** The specified Camera will be used. When using this option you must also specify the *Grabber Name* and the *HW Unit Name*. To find the camera name, search the application log file for "Camera Name". It is also possible to use only a part of the Camera Name, but the part of the name must be unique among all the available cameras.

**-cw "Command Wait"** The application runs the specified macro command and waits until it is completed

**-f "Filename"** The application tries to open the image pointed to by the *Filename* parameter.

**-g "Grabber Name"** The image grabber (driver) of the specified *Grabber Name* will be used and no camera-selection window will show at startup. Set the parameter value to "last" and the most recently used grabber/camera will be loaded. As the grabber name, you must use internal name of the grabber, not the one you see in the startup dialog window.

Either use one of the following names or search the application log file for the phrase "Grabber Name" (see 1.2 Fixed Grabber Startup).

**-gn** No grabber (camera) driver will be loaded. Image acquisition will not be possible.

**-h "HW Unit Name"** The specified HW Unit will be used. When using this option you must also specify the Grabber Name. Two HW Units (e.g. two DS-U2 cameras) can be run by one grabber/driver. The available HW Units are listed in the *Acquire > Select [Camera Name]* window. Set the parameter value to "select" and the selection box will be shown. When using this option you must also use the **-g** switch. To find the HW Unit Name, please search through the application log file for "HW Unit Name".

**-i "Config Name"** A user can specify the configuration to be used by *NIS-Elements*. The default configuration is saved in the "C:\ProgramData\Laboratory Imaging\Platform" folder. If you want to create a custom configuration, use the **-i** switch, e.g.: `nis_ar.exe -i "My_Configuration"`. The configuration of this instance will be saved to the "C:\ProgramData\Laboratory Imaging\My\_Configuration" folder.

This is useful when using two cameras on one microscope for example. You can create a separate program shortcut which uses the **-i** switch to load a different configuration.

**-l “Language”** The application will run in the specified language (if available). Use three-letter language codes according to the ISO 639 [[http://en.wikipedia.org/wiki/List\\_of\\_ISO\\_639-2\\_codes](http://en.wikipedia.org/wiki/List_of_ISO_639-2_codes)] standard.

**-m “Macro File”** The application will execute the macro file (\*.mac) pointed to by the *Macro File* parameter.

**-mw “Macro File”** The application will execute the macro file (\*.mac) pointed to by the *Macro File* parameter and wait until it is completed.

**-p “Command” [“Param”]** The application window will be placed according to the *Command*. Possible values are:

**left, right, top, bottom** The window is placed on the corresponding monitor screen.

**minimize** The software window is minimized to the task bar on startup.

**monitor N** The window is placed on the *N*th monitor.

**rectangle (x0,y0,x1,y1)** The window is placed in the specified rectangle (in workspace coordinates). The coordinates must be in parenthesis without any spaces.

**-passive** The application will not connect to any device on startup (*NIS-Elements* will start with blank Device Manager). Camera selection is not affected by this option.

**-q** A new instance of the application will run. By default, only one instance of *NIS-Elements* is allowed to run.

**-s “Settings File”** The application will use the specified *Settings File* to load and save its settings. The settings file can be created from MS Windows Start menu by the *NIS Settings Utility*.

**-var, -vbr, -vd, -vf** For testing purposes, you can switch the application to run as if you have a different version (AR/BR/D/Freeware) installed. However, you have to have the corresponding licence written in your key (Even if you normally run NIS-AR, the -vd switch will not work to run NIS-D unless you have also the NIS-D licence).

#### Note

*Mind that the set of functions you will get once the software is run depends on the setup file used for installation (for example, some functions of NIS-D are not included in the NIS-AR setup so these will not appear in the software run by: `nis_ar.exe -vd`*

## 1.2 Fixed Grabber Startup

Each time *NIS-Elements* is run or the grabber / camera / HW unit is changed, four parameter values are written to the application log file (C:\ProgramData\Laboratory Imaging\Platform\Logfiles):

- Grabber Name
- HW Unit Name
- HW Unit Connection String
- Camera Name

Let's say we want to set the last used configuration for every *NIS-Elements* start:

1. Lookup the parameters in the most recent log file. E.g.:

**Grabber Name** CLxGrabberDriverSim

**HW Unit Name** SimGrabber

**HW Unit Connection String** (empty)

**Camera Name** Sim\_Camera\_Color



2. Modify the shortcut, E.g.:

```
"c:\Program Files\NIS-Elements\nis_d.exe" -g CLxGrabberDriverSim -h SimGrabber  
-cam Sim_Camera_Color
```

3. Double-click the shortcut. *NIS-Elements* will start without asking for grabber selection



## 2. Installation and Settings

### 2.1 Installation and Updates

#### 2.1.1 The Installation DVD-ROM Content

- *NIS-Elements* software setup file
- Drivers and utilities for the HASP Key
- Drivers for selected cameras
- User's Guide in the PDF file format
- Sample image database
- Sample ND2 image sequences

#### 2.1.2 NIS-Elements Installation Steps

##### Caution

---

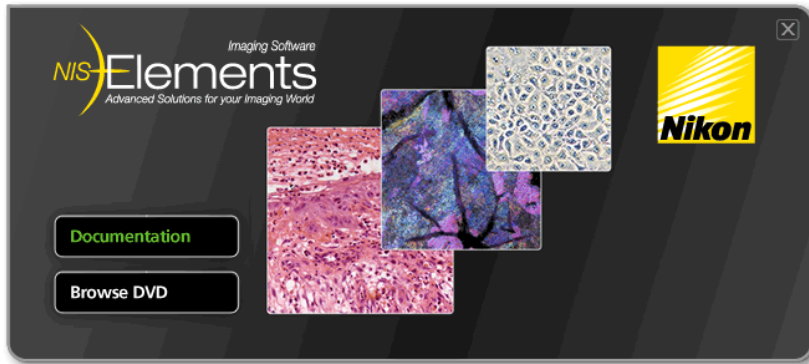
*You have to possess the administrator rights to your computer to be able to install NIS-Elements successfully.*

##### **Quick Guide**

- Insert the installation DVD in the DVD-ROM drive. A window automatically appears.
- Install the selected *NIS-Elements* software version, additional modules and device drivers.
- Plug the provided HASP key into the USB port of your PC.
- Run *NIS-Elements*.

## **Step by Step**


- 1) When the installation DVD is inserted, a selection window appears automatically.



Select the software package to be installed. Select the one you have got the licence for and which is properly coded in your HASP key. The installation wizard welcome-window appears. Click *Next* to continue.

- 2) **Install Local Options**

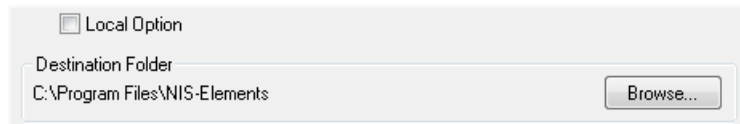*Figure 2.2. Local Options installation*




Select whether to install *Local Option* or not. Define the folder where *NIS-Elements* should be installed. We recommend to use the predefined directory. If you want to change the directory anyway, press the *Browse* button and select a new one. Otherwise click *Next*.

### Note

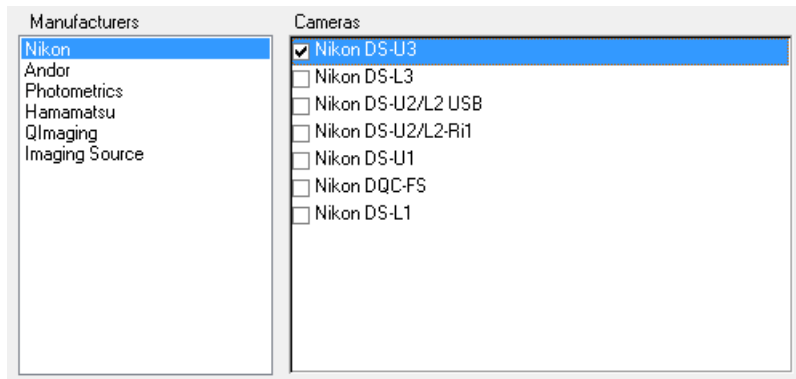*The Local Option installation provides some advanced features which did not pass the quality assurance procedure yet. We recommend to wait until they are released officially.*

3)




Now, select the cameras which will be used with *NIS-Elements*.

Note

*When a camera is installed, CPU Power Management C-State is automatically disabled to secure stable camera operation. Please restart your computer after changing the C-State settings. C-state is automatically restored to its original condition after uninstalling NIS-Elements.*

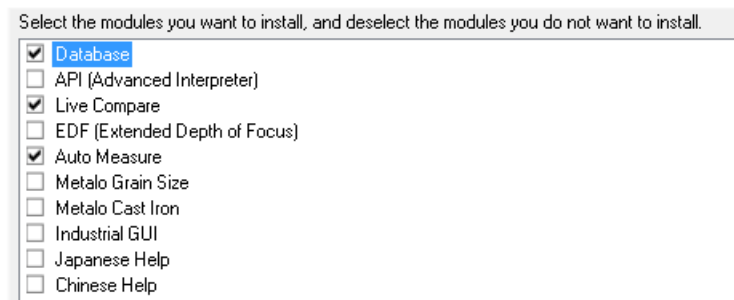4)




If your licence contains some additional modules besides the *NIS-Elements* core software, please select them in this window.

Note

*Any module selected will be installed along with NIS-Elements automatically. However, you might not be licensed to use it. The module will run after you get the corresponding code registered in your HASP key.*

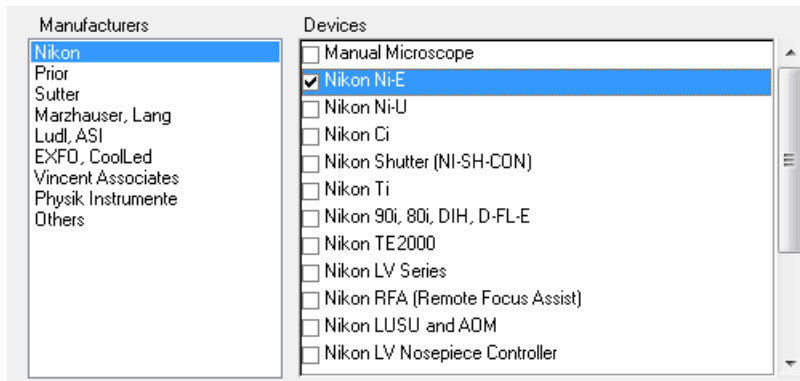5)



Select the devices which will be used with *NIS-Elements*. Finish the installation by clicking the *Install* button.

Note

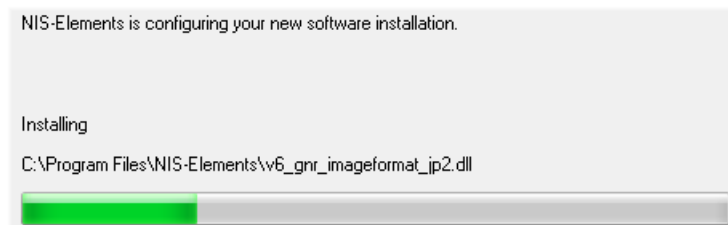
*NIS-Elements* may not be connected to devices after installation depending on the PC status at installation. If so, execute the *Modify installation* command of the Windows Start menu (see the next step) to perform *Repair* that corrects the previously installed status.

Warning

For Windows 7 users: Some devices do not work correctly after recovery from Sleep mode. Please turn off the Windows 7 Sleep Mode to prevent possible problems:


1. Run the *Start > Control Panel* command on the main Windows toolbar.
2. Make sure the control panel is displayed in the *Category mode* (*View by* in the top-right corner).
3. Click *Hardware and Sound*
4. In the *Power Options* section, click *Change when the computer sleeps*
5. Set the *Put the computer to sleep* option to *Never*
6. Confirm the settings by the *Save changes* button and close the *Control Panel* window.

6)



The setup creates a new program group in the *Start* menu containing the following items: *NIS-Elements* application shortcut, the HASP key information shortcut, the *Modify Installation* shortcut (for adding hardware drivers, modules, etc.), the *Uninstall* procedure, and the Send Info Tool. A shortcut to *NIS-Elements* is created on the desktop too. These changes affect all user profiles of your local Windows operating system.

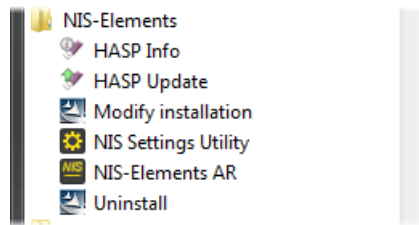Figure 2.7. Start menu



#### Note

Clicking the *Uninstall* command deletes all installed files from disk, and removes the *NIS-Elements* program group from the *Start* menu as well as it removes the desktop icon.

### 2.1.3 Command Line Installation Options

See 1.1 Command Line Startup Options for details about how to use command line switches. The following switches are to be used with the *NIS-Elements* installation file to modify behavior of the installation wizard.

**-all** This parameter unhides options in the setup which are hidden by default. It concerns mainly drivers of devices which were formerly supported but are no longer available.

**-new** If this command line parameter is appended, the setup checks if any other setup has already been installed. If so, the setup restarts and runs in an independent mode. The update mode is disabled. This enables to install two or more builds together on one operating system.

**-xp** This parameter enables the user to install *NIS-Elements* on Windows XP even though this operating system is NOT supported officially.

### 2.1.4 Additional Module/Device Installation

You may need to install a device or an additional module after the *NIS-Elements* main system installation.

- Go to [Start menu > Programs > *NIS-Elements*] program group.
- Select the *Modify Installation* command.
- Follow the installation wizard instructions. Select the check-boxes by the items you would like to add.
- Finish the installation.

## 2.1.5 Sample Database Installation

If you chose to install the Sample Database, a new subdirectory “Databases” is created inside the *NIS-Elements* installation directory (e.g. C:\Program files\*NIS-Elements* \Databases\...). The “Sample\_Database.mdb” file is copied to there along with database images (stored in subdirectories). An administrator username/password to access this database is set to:

- Username: “sa”
- Password: “sa”

## 2.1.6 Software Copy Protection

*NIS-Elements* is delivered with a hardware key (also called HL = hardware licence or HASP).

Figure 2.8. Hardware key



The key contains information about the software licence and allows users to run the corresponding software. A warning message is displayed when user starts *NIS-Elements* with an incorrect hardware key. Please do not connect the USB hardware key until after the software installation is finished. The utility called *HASPinfo* is installed to the installation directory. It enables the user to view information about the software licence and is accessible via the *Help > HASP Info* menu command. See also 2.5 Hardware licences.

## 2.1.7 Device Updates

The main goal of the Device Updates setup is to add new devices/cameras or to fix problems that are localized in the drivers only and do not affect any other *NIS-Elements* functionality.

Device Update setup requires installation of the corresponding *NIS-Elements* version. The setup contains only .dll files supporting a new device or some corrections and upgrades of already existing drivers. This update form keeps *NIS-Elements* stable and the update file size small.

Device Update numbering is consecutive and appropriate to its *NIS-Elements* version until a new full setup of *NIS-Elements* is released (Major version, Service Pack or HotFix) - then the numbering starts from the beginning.

Device Updates are cumulative and contain all the changes from the previous updates. Also full *NIS-Elements* setups (Major versions, Service Packs and HotFixes) will contain all the changes from the previously released Device Updates.



## 2.1.8 Fixes

Fixes represent driver changes which serve for testing purposes or for correcting problems occurring only under special conditions. Fixes are packed into a zip file containing modified driver files. Fixes are usually user requested, therefore they are not considered as Device Update candidates.

## 2.1.9 Installing the Database Module on 64-bit Systems

The *NIS-Elements* installation file contains 64bit MDB drivers which improve the speed of the database module on 64bit systems. However, 32bit version of MS Office 2010 cannot be used along with these 64bit drivers. You can install these drivers, if:

- you run 64bit version of MS Office 2010
- you use other version of MS Office such as Office 2013
- you do not use MS Office at all

Find the drivers in the *NIS-Elements* folder, typically: "C:\Program Files\NIS-Elements\Drivers\Database\_MDB\"

## 2.2 Additional Modules Available in NIS-Elements D

### 2.2.1 Acquisition and Analysis Modules

**EDF Module (Extended Depth of Focus)** The EDF module allows you to combine an existing Z-stack of images into one focused image by picking the focused regions from each frame and the pieces together.

**High Dynamic Range** This module brings the functionality of capturing and creating HDR images within *NIS-Elements*.

**INTADV (Interactive Advanced Macro - API)** This module dramatically extends the macro programming capabilities of *NIS-Elements*. It contains a set of user-interactive macro functions and the ability to import functions from external DLLs.

**LCOMP (Live Comparisons)** This module provides several comparison modes which can help to identify and visualize differences or similarities between two images.

**MEAS (Automatic Measurement)** This module extends the measurement abilities of *NIS-Elements* by adding number of automatic processing features. Particularly, it enables users to threshold the image and process and measure the binary layer. See also 8.3 Automated Measurement.

### 2.2.2 Industrial Modules

**Concrete** (*requires: Local Option*)

This module is for automated and semi-automated measurement of air void characteristics of hardened concrete.

**Filters Particle Analysis** (*requires: 64bit NIS-Elements*) This industrial module is aimed at analysis of filters according to the ISO 16232 standard.

**Industrial GUI** After the *Industrial GUI* module is installed, the *View > Simple Control* command is added to the application. The command activates a simple layout intended to be used in industry imaging applications.

See 3.7 Simplified User Interface

**Layer Thickness** This module adds the *Applications > Layer Thickness Measurement > Measure Layers* command to the main menu. The module can perform different types of layer thickness measurement.

**Measurement Sequencer** With this module, you can:

- Create custom measurement sequences.
- Install predefined measurement sequences suitable to measure welds of different shapes.
- Install predefined measurement sequences suitable for hardness measurement (Vickers, Brinell, Knoop)

**Metalog - Cast Iron Analysis** After the module installation, the *Applications > Metallography > Cast Iron* command appears in the main menu. Run this command to display the cast iron measurement layout.

**Metalog - Grain Size Analysis** Adds the *Applications > Metallography > Grain Size* command to the main menu. The command displays a layout dedicated to grain size measurement.

## 2.2.3 Device Modules

**Digital Mirror Device** (*requires: Local Option*) DMD support. E.g.: Nikon Ti-Lapp, Andor Mosaic 3, Mightex Polychrome 400.

**Dual Camera Support** Support for dual camera drivers (two mono cameras behaving as a single dual-channel camera). Specific camera models are supported.

**Monochromator** (*requires: Local Option*) Thirdparty monochromator support. E.g.: Till Photonics Polychrome V.

**Revolution DSD** (*requires: Local Option*) Andor Revolution DSD confocal support.

**SFC** (*requires: Local Option*) Swept field confocal support.

**Slide Loader** (*requires: Local Option*) This module enables *NIS-Elements* connect and control a slide loader.

**Splitter DV** *Dual View* and *Quad View* support.

**Stage** This module enables to connect and control motorized XY stages. All non-Nikon stages require this module.

**Stage Incubator** This module enables to connect and control gas and temperature-controlling chambers.

**Structured Light** (*requires: Local Option*) Support of structured-light-based devices (OptiGrid, ViCo).

**Sutter** Third party filter-wheels support.

**TTL/Analog IO** This module enables *NIS-Elements* to send and receive TTL and analog signals via a DAQ board.

**Wavelength Switcher** This module enables *NIS-Elements* connect and control wavelength switchers.

**Well Plate Loader** This module enables *NIS-Elements* connect and control a well plate loader.

**X-CITE** Third party illumination device support.

**Z Drive** This module enables connect and control motorized Z drives.

See also 4.1.6 Connecting a Device to *NIS-Elements* .

## 2.2.4 Other

**64bit NIS-Elements** Some functions and applications run only under the 64bit version of *NIS-Elements*. Unlike with 32bit applications, the size of memory available to a single application is not restricted on 64bit.

**DBASE (Database)** If the database module was installed, the system can create and control image databases. A database can help you to efficiently organize your image archives and manage additional information related to the images.

**Local Option** When Local Option is installed (see Step 2), *NIS-Elements* will provide some advanced features which did not pass the quality assurance procedure yet. We recommend to wait until they are released officially.

**MS Excel** *NIS-Elements* supports export of results straight to an MS Excel sheet (see 8.1.7 Exporting Results). MS Excel is not part of the *NIS-Elements* and must be purchased and installed separately.

## 2.3 User Rights

User management is very useful when a single workstation is shared by number of users. Some user accounts may have administrator rights to *NIS-Elements* while other's privileges are quite restricted. A per-person system may be established or user accounts may be shared. The following principles are utilized within *NIS-Elements*:

**MS Windows accounts support** *NIS-Elements* can assign privileges to MS Windows accounts. Whoever is logged in the operating system and runs *NIS-Elements* obtains a default set of privileges. System administrator can then further restrict or extend the user's rights.

**NIS-Elements password protection** It may be not comfortable enough to log the current user off from MS Windows when a user change is required in *NIS-Elements*. For such case a completely independent list of user accounts, which is not connected with MS Windows users, can be created. Then just restart of *NIS-Elements* is needed to switch the user.

**Private vs. Shared storage** Every user-created item (setting) important in the work-flow can be protected from unwanted change in two levels. First, the creator of the item (optical configuration, layout, objective, etc.) can save it as *private*. Unless its status is changed to *shared*, no other user will even see this item within the lists throughout the application. The second level of protection concerning all *shared* items is provided by assigning users to *Groups* and *Privileges* groups.

**Groups** Every user is a member of a group. The group enumerates shared items (Hardware Configurations, Objectives, Layouts, etc.) which will be visible for its members. *NIS-Elements* administrators can create any number of groups and select the items to be visible.

**Privileges** There are basic groups of privileges which allows users or restrict them to perform certain actions. This concerns for example modifying macros, reports, selecting cameras, etc. There are four levels of privileges by default, but further can be created by an administrator:

- Admin
- Common
- Guest
- Default

### 2.3.1 NIS-Elements authentication

If you have chosen “*NIS-Elements* (Password protected)” user authentication method, a login dialog window appears at every start-up of *NIS-Elements*. Enter a user name and a password for existing user account. If the option *Allow Create New User on Startup* is enabled, you can create a new user account without needing an administrator. This account is assigned to the *Default* group of users and *Default* group of privileges automatically.

### 2.3.2 Creating a Shared Layout

Please read the following example on user management. Step by step, we describe the creation of a shared layout which can not be modified by members of the *Common* and *Guest* groups.

- 1) Run the *View > Layout > Layout Manager* command. Set at least one of the defined layouts as *Shared*. (see 3.1.3 Arranging User Interface)
- 2) In the left column of the window, select *User Rights*.
- 3) Make sure you are logged as a user with the *Privileges to Modify user rights*. If not, use the  button to log in under a different account which has the privilege.

#### 4) Set layout visibility

Select the *Groups* tab. Here you can see at least one group of items called *Default*. Either select one a group or create a new one by the *New* button and choose *Layout* within the list of items. All layouts set as *shared* will be listed on the right side within the *Accessible items* field. Make sure the check boxes next to the layouts which you allow to be used by the group are selected.

#### 5) Disable layout modification

Select the *Privileges* tab. Select the group of privileges which should have not the permission to modify the selected shared layouts. In the *Privileges* field on the right, make sure you de-select the *Modify Shared Layouts* option.

#### 6) Apply the new policies


Select the *Users* tab. In the list of users, select the ones which you would like to apply the policy to (users can be selected one by one or by multi-selection with the  or  key pressed). Then, assign the appropriate *Group* and *Privileges* by selecting them from a pull-down menu on the right side of the window.

- 7) Confirm the changes by the  button.

### 2.3.3 Closing Another Instance of NIS-Elements

Only one instance of *NIS-Elements* can be running on one computer. If a second user starts *NIS-Elements* when another instance of the program is already running, the following dialog appears.

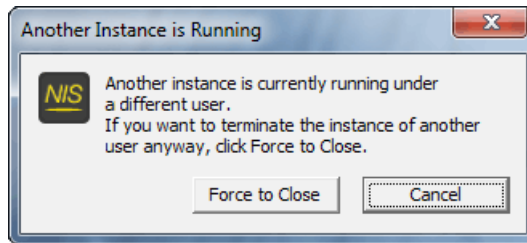Figure 2.9. Another instance running



The user can press the *Force to Close* button to close the running instance. This option is only available if there is no acquisition experiment in progress.


### 2.3.4 User Rights Options

**User Authentication** The *Windows* authentication utilizes Windows accounts to automatically log in to *NIS-Elements*. The *NIS-Elements (Password protected)* authentication utilizes application user accounts and passwords at the start-up of *NIS-Elements*. Check the *Allow Create New User on Startup* option to enable creating of a new user in the *Login* dialog window.

**Current user** Displays currently logged in user.

**Log As** Opens the *Login As* dialog which enables you to login as a different user e.g. an administrator with more user rights for the purpose of changing *User rights* settings.

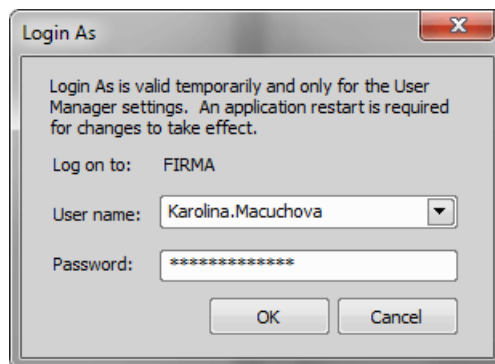Figure 2.10. Login As



Select the user name you want to log-in as from the drop-down menu. Enter a valid password and press **OK**. The new user identity is available only while the User manager is opened.


**Statistics** Edit path and file name of a database file for saving statistics.

**Export/Import user rights** These buttons enable anyone to save/load complete settings of user accounts and their privileges to/from an external XML file. Standard *Open* and *Save As* windows appear. Having the data exported may be useful when copying the settings to other computers.

## 2.3.5 Users Tab Options

**New** Enables to create new *NIS-Elements* user account. You have to possess the privilege Modify User Rights.

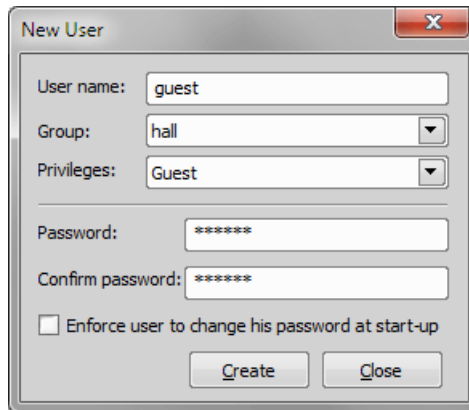Figure 2.11. Creating new user



Enter a new user name, assign him to one of the existing Groups and group of Privileges and set the password. Check the *Enforce user to change his password at start-up* option to make the user change his password at his first login into the *NIS-Elements*.

### **User name and password properties**

Length of the user name and password is limited to 50 characters. You can use any combination of characters, symbols or numbers (except the Windows ones: ", ;, \*, ?, ", \, /, <, >, |").

**Remove** Removes selected user account. You can select multiple accounts and delete them at once.


**Copy To** Choose a user from the pull down menu whose group and group of privileges will be changed according to settings of a selected user.

**Duplicate** You can easily create a new user by duplicating the existing. This preserves all account settings under a different name. If you have selected to duplicate more than one user, enter a name prefix, which is added in front of the original name for all selected duplicated accounts. User which was created as a duplicate has an empty password and is forced to change his password at next application startup.

**User Accounts List** This table displays all existing user accounts. You can sort them according to any of four columns. The first column displays type of the account. The second one shows user names. The third one displays assigned group of users. The fourth one displays assigned group of privileges. Select one or more users whose settings needs modification.

**Set password** To change password of a selected *NIS-Elements* user account, press the *Set password* button. You have to possess the Privilege Modify User Rights. Otherwise you can change only password of the current user. Following window appears:

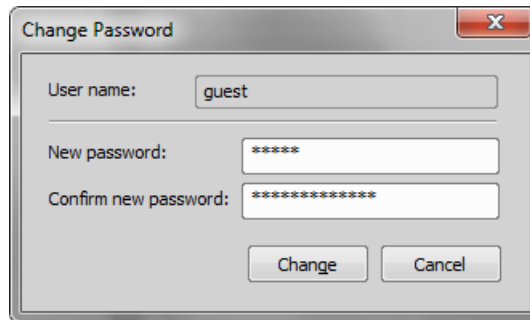Figure 2.12. Reset of password




Name of the user is displayed. Enter and confirm new password. When done, press the *Change* button.

**Enforce user to change his password at start-up** Check this option to make the user change his password at his first login into the *NIS-Elements*.

**Enumerate All Windows Users** This button starts copying MS Windows accounts to *NIS-Elements*. The accounts are assigned to user groups (admin, common, guest) according to their Windows permission settings.

## 2.3.6 Groups Tab Options

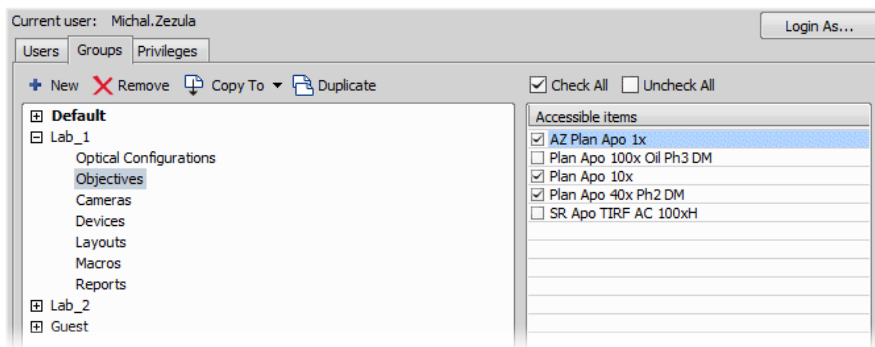Figure 2.13. Groups tab



**New** Enables to create new group of users. You have to possess the privilege Modify User Rights. Enter a name of new group.

**Remove** Removes selected groups. You can select multiple groups and delete them at once.

**Copy To** Choose a group from the pull down menu which settings will be changed according to selected group.

**Duplicate** You can easily create a new group of users by duplicating the existing. This preserves all group settings under a different name.

**List of Groups** All defined groups are listed in this window. Each group contains subsets. If you click on a subset, its items appear in the *Accessible items* list.

**Accessible items** This list shows items of a selected subset. Every item has its own check box. If the check box is checked, all users assigned to the corresponding group will see the item. Otherwise they will not see the item.

**Optical Configuration** Displays all shared optical configurations.

**Objectives** Displays all shared objectives.

**Cameras** Displays all installed cameras. From the pull down menu below select a camera which is used as default for users without the privilege Select Camera.

**Devices** Displays all installed devices.

**Layouts** Displays all shared layouts.


**Macros** Displays all shared macro commands. Check a macro to make the content of it visible to this Group.

**Reports** Displays all shared reports and report templates.

**Sequencer Definitions** Displays all shared *Measurement Sequencer* definitions.

## 2.3.7 Privileges Tab Options

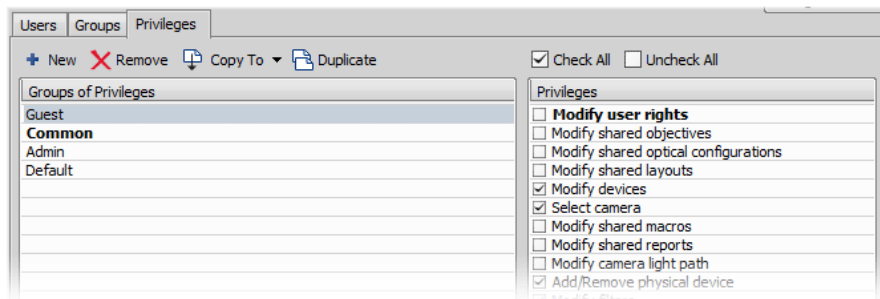Figure 2.14. Privileges tab



**New** Enables to create new group of privileges. You have to possess the privilege Modify User Rights. Enter a name of new group.

**Remove** Removes selected groups. You can select multiple groups and delete them at once.

**Copy To** Choose a group from the pull down menu which settings will be changed according to selected group.

**Duplicate** You can easily create a new group of privileges by duplicating the existing. This preserves all group settings under a different name. If you have selected to duplicate more than one group of privileges, enter a name prefix, which is added in front of the original name for all selected duplicated accounts.

**Groups of Privileges** A list of all groups of privileges is displayed.

**Privileges** A list of privileges is displayed. Every item has its own check box. If the check box is checked, it grants all users assigned to the corresponding group relevant privilege to access, change, modify, etc.



---



 Note
 

---

Users that do not have the privilege “Modify Shared Optical Configuration” can temporarily change the brightness settings (e.g. exposure time, gain, etc.) in a shared optical configuration and use such adjusted optical configuration for example in Multichannel acquisition. But all such modifications will not be stored and they will disappear when the application is restarted.

## 2.4 NIS-Elements Preferences


### 2.4.1 Adjusting Program Preferences

- 1) Run the *Edit > Options*  command. The *Options* dialog window appears.
- 2) Select a tab which contains requested options. Options are sorted into several groups:
  - General** Options concerning basic image operations. See 2.4.2 General.
  - Appearance** Options concerning the graphical user interface. See 3.6 Appearance Options.
  - Open Next** Options concerning the *File > Open/Save Next > Open Next* menu command.
  - Save Next** Options for the *File > Open/Save Next > Save Next*  command.
  - Macro** Configures key shortcuts to macros and sets macros to run automatically on startup See 10.6 Macro Preferences.
  - Measurement** See 8.1.4 Measurement Options for more information.
  - Data export** See 8.1.7.2 Data Export Options for more information.
  - User rights** See 2.3 User Rights for more information.
  - Layout Manager** See 3.2.1 Layout Manager for more information.
- 3) Make any changes you need in preferences and use the following buttons to manage them:
  - Defaults for this Page** Restores default settings of the currently displayed options tab.
  - OK** Confirms and saves changes made to the preferences. The dialog window is closed afterwards.
  - Cancel** Cancels all changes made to the preferences. The dialog window is closed afterwards.
  - Apply** Applies changes made in preferences but keeps the *Options* dialog window open. You can apply changes individually on each tab.
  - Help** Displays the relevant help page.

### 2.4.2 General

#### Documents and History

---


**Use fix path for images** The defined directory is always used when using the *File > Open*  or *File > Save As* command.

**Allow only one opened document** If checked, only one document at a time can be opened.

**Limit number of opened documents** When this options is checked, the user is limited to a single opened image at one time.

**Capture always creates a new document** If checked, every acquisition creates a new image.

**Use last LUTs on image open** When a new image is opened, this option turns the look up tables automatically on, and copies the settings from the current or the last opened image. Some image formats (JP2, ND2) can contain the LUTs settings. If such image is opened, the saved LUTs are loaded instead of the most recently used.

**Use AutoLUTs on image open for images without LUTs information** This option keeps the AutoLUTs feature of the *File > Open*  window always ON, and it applies AutoLUTs to the image after opening. This option is *ignored* if LUTs settings are saved in the image, or if the *Use last LUTs on image open* option is applied.

**Show mapping dialog in organizer after drag and drop** Displays the *Mapping* window every time an image is inserted to a database table. It enables you to check/modify mapping of the image info values to the database table fields.

**Open new ND views to new window** Displays each newly generated ND view in a separate window.

**Enable authentication** Authentication verifies each image frame and detects changes in image data made after acquisition. Modifications are indicated in the *Image Properties* dialog. See *File > Image Properties*.

---

#### Note


*Authentication works with images acquired by the Capture button, the ND Acquisition dialog(s) or using the JOBS module.*

**Show image info window on save** Displays the *File > Image Properties* window every time an image is saved via the *Save As* command.

**Detect sequence on image open** When you select to open an image that is a part of image sequence such as 001.JP2, 002.JP2..., it is recognized automatically and you are offered to convert it to an ND2 file.

**Show binary layers' contours** Check this item to display contours of binary layers.

**Use zero based time scale for ND documents** Select this option to ensure that the first frame of a time sequence will always start at 0.0s.

**State of saturation indicator after start** Select what to do with the saturation indicator settings (  ) after *NIS-Elements* is restarted. Turn it ON, OFF, or remember the last setting.

## Rotation, Flips and Shifts

**Apply to overlaid binary layer** If checked, in overlay mode both binary and color images are rotated, shifted and flipped.

## Optical Configuration

**Save changes to optical configuration automatically** Make changes to the current optical configuration without warning? If selected, some setting changes are written to the optical configuration immediately and the exclamation mark would not appear by the OC name. Select the settings to be changed without warning:

**Exposure Only** Camera exposure time only.

**All Camera Settings** All camera settings such as exposure, gain, resolution, etc.

**All Brightness Settings** All Camera Settings plus selected settings of illumination-related devices such as laser power, light power, etc.

**Select corresponding optical configuration when filter or nosepiece changed** Enabled if “Unselect” in the option below is chosen.

Allows *NIS-Elements* to automatically change the *optical configuration* based on the filter or nosepiece selected on the microscope.

#### When Optical Configuration setting changed

**Keep Selected** The Opt. configuration name is marked by the “\*” sign and the change is offered to be saved.

**Unselect** Every change of Optical configuration setting results in its deselecting and switching to user settings.

**Save all changes** All changes regarding the Opt. configuration settings are updated automatically according to the current state.

**Temp** *NIS-Elements* uses the Temp directory for storing temporary data when there is not enough RAM available. You can redirect the system to use a directory on another hard disk, which may speed up the whole system.

**Defaults for this Page** Restores the default setting for General Options.

## 2.5 Hardware licences

### 2.5.1 What is a hardware licence?

Together with the *NIS-Elements* software package you received a hardware USB key - so called hardware licence (HL) or HASP. This key contains information about your software licences. If the USB key is not inserted into the computer USB port or if the key does not contain the appropriate licence *NIS-Elements* cannot be executed.

Figure 2.15. Hardware USB key




---

#### Note

Formerly the USB key was called “HASP key” often abbreviated to “HASP”.

## 2.5.2 How do I find out the content of my USB key?

To verify the contents of the licence recorded on your USB key, service tool called *HASP Info* is installed together with *NIS-Elements*.

1. Insert the USB key into your PC.
2. Go to the *Start* menu of your operation system and search for the installed *NIS-Elements* folder.
3. Click on *HASP Info* inside this folder. Window containing information about your key (modules which can be executed using this key) is shown. See also 2.2 Additional Modules Available in *NIS-Elements D*.

### Note

Displayed information can be saved into a text file (\*.txt) using the **Save info** button. Sending this text file (e.g. to the technical support) is the easiest way of conveying the full information about your USB key. This can be useful for verifying whether a specific function is available with your current licence.

## 2.5.3 Is there an update for my software?

New updates of *NIS-Elements* are released to serve as a tool for:

- expanding the software functionality
- adding the support of new devices (cameras, microscopes, etc.)
- bug fixes

You can easily find out whether an update is ready for your *NIS-Elements* version:

- If the computer running *NIS-Elements* is connected to the Internet, execute *Help > Check Version Update* from the program main menu. A web page with information about your USB key is automatically loaded inside your browser.
- Frequently the laboratory computes are not connected to the Internet. In such case write down your licence number (USB key), find a computer having Internet connection, go to this web page: <http://www.nis-elements.cz/version> and fill in the form using your licence number.

**Version** List of released program versions compatible with the USB key.

**Release date** Release date of the particular version.

**Available for free** Is the software update available for free? If so, "Yes" is shown in the column.

**Available after purchasing** For updates not available for your original licence, "Software Upgrade Agreement" (SUA) has to be ordered.

## 2.5.4 What is SUA?

SUA - "Software Upgrade Agreement" is a time limited agreement enabling the owner of the *NIS-Elements* licence to update the software during the next 12 months for free. The business model works as follows:

- After buying the licence to *NIS-Elements* it needs to be activated. Once activated the activation date is recorded into the manufacturer's database.


- From this activation date you get a 13 months period during which all updates (even updates to new software versions) are available for free.
- After the expiry of the 13 month period from the activation date, all new updates are available only after buying a SUA. 1× SUA = free updates for the next 12 months.



## 3. User Interface

### 3.1 Advanced Layout

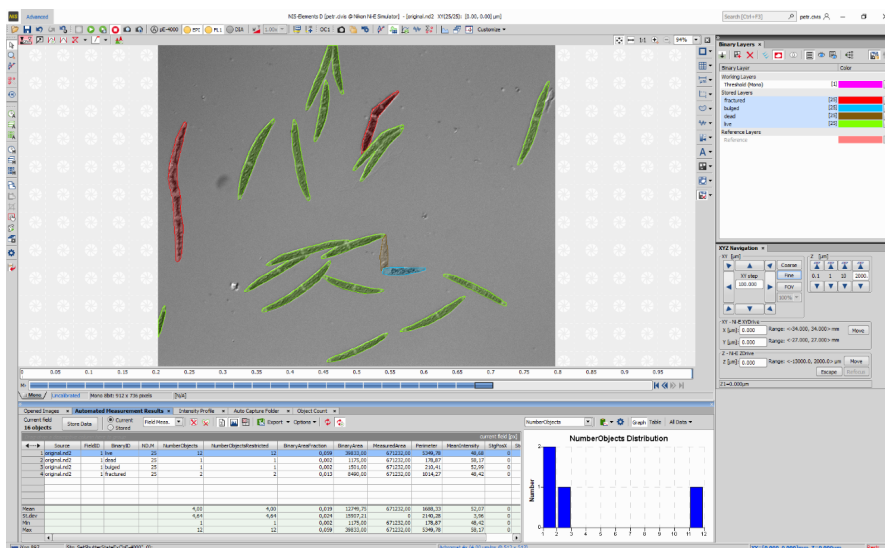Figure 3.1. NIS-Elements Advanced Layout



Context menu over the *Advanced* tab enables to load original settings of the selected predefined layout (*Load Default*), lock the selected layout so that its components cannot be moved or closed (*Lock Layout*), *Save Current Layout As* a new layout or open the *Layout Manager* (3.2 Modifying Layouts).

### 3.1.1 Main Window Components

#### 3.1.1.1 Main Menu

All *NIS-Elements* functions are accessible from the main menu at the top of the screen. Menu commands are grouped according to their purpose.


#### 3.1.1.2 Toolbars

There is a default set of toolbars, each toolbar containing number of buttons. There is also one fully customizable toolbar - the main left toolbar - to which any button can be added. Every button or whole toolbar can be hidden by the user. Please see 3.1.3 Arranging User Interface for further details.

### 3.1.1.3 Status Bar

The status bar at the bottom of the screen displays the following information:

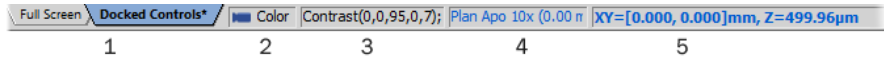Figure 3.2. The application status bar



1. This part of the status bar displays available layouts.

#### Note

*The layout Tabs may be hidden when the Show Layout Tabs option in the 3.1.3 Arranging User Interface window is deselected.*

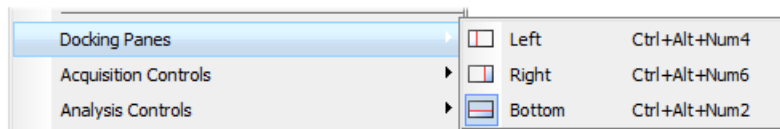
2. This status bar section displays the type of the currently selected camera.
3. Here you can get information about the most recently performed command. The FPS / Exposure / Focus info is shown in case of live image. The black bar indicates the focus rate. Longer black bar represents more of the image in focus.
4. This section show the name of the current objective.
5. Current coordinates of XY (Z) stage are shown in this part of the status bar.

### 3.1.1.4 Docking Panes

Docking panes are square spaces inside the application window, where you can place (dock) any of the control panels. There is one docking pane available at the Right, Bottom, and Left side of the application screen.

#### **To Display a Docking Pane**

1. Go to the *View > Docking Panes* sub-menu and select the pane you would like to display.



#### Note

*The Docking Panes sub-menu can be also displayed by right-clicking into the empty application screen.*

2. The docking pane appears, either empty or with some window(s) docked inside.
3. Repeat this procedure to display more docking panes.

#### **Handling Control Panels**


Various control panels can be displayed docked within the docking panes or they can be floating. See the following picture:



Figure 3.4. The Docked Control Panel Caption



To handle the control panels (CPs), you can:

**Open recently closed CP** Locate the  button on the main left toolbar. When you click it, the list of recently closed CPs appears. Pick one to display it again.

**Add CP to a docking pane** Right click inside a docking pane (3) to display the context menu. Select the control panel to be displayed. If the window is already opened somewhere else (in another docking pane or floating), it closes and moves to the new destination.

**Close CP** Click the cross button (2) in the right top corner of the tab.

**Drag CP** Drag any CP by the tab and drop it somewhere. If you drop it by the edge of a docking pane, it will create another column of this pane. If you drop it over the caption of another CP, it will be docked in the same pane as a new tab. If dropped somewhere else, the CP will be floating.

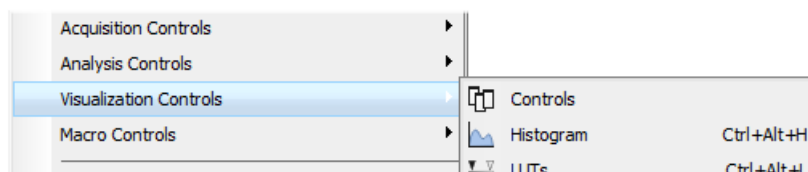
A color frame appears when you place the mouse cursor dragging a CP over the edge of a docking pane or a caption of another CP. It indicates that if you drop it, its placement will be handled automatically.

**Minimize the docking pane** Click the arrows in the top left corner. The pane minimizes to a stripe by the edge of the screen. It can be restored to its original position by double clicking this stripe or by clicking the arrows again.


**Close the docking pane** Click the cross button (4) in the docking pane caption. Or you can right click the pane and unselect the *Docking View* option.

**Dock/undock CP** To dock (and undock) a CP, double click its tab.

**Display CP** Another way to display a CP is to go to the *View* menu and select the desired control panel. After that, the CP appears on the screen - floating or docked. Positions of the windows are being saved by the system so that each control panel appears in the same position as it was before it being hidden. The controls are sorted to several groups.



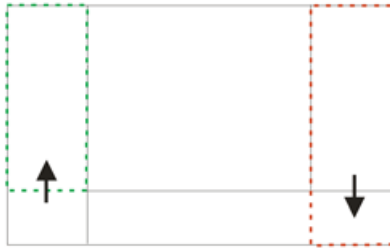
**Shrink/Expand dockers** Having more docking panes opened, a situation where there is not enough room for the control panels can occur. In such case, as you can see in our example, the *Shrink bottom* (green color) or *Expand bottom* (red color) commands shall be used.




1. Right click the pane you would like to shrink/expand. A context menu appears.
2. Select p.e. the *Expand /Shrink right* command. When one of the panes shrinks, the neighboring pane expands to the emptied corner and vice-versa.

### 3.1.2 Image Window

Tools affecting the appearance of the current image are gathered within the image window toolbars (the top image toolbar and the right image toolbar). There are the following buttons by default:


 **Enable LUTs (L)** Applies LUTs to the image. See 6.5 LUTs - Non-destructive Image Enhancement.



 **Keep Auto Scale LUTs** Applies the AutoScale command to the image continuously.

 **Auto Scale** Performs automatic setting of LUTs.

 **Reset LUTs** Discards the LUTs settings.

 **Show LUTs window** Opens the window with LUTs.


 **Pixel Saturation Indication (Ctrl+Shift+S)** Turns on/off pixel saturation indication without setting on/off LUTs. Select the highlighting color from the nearby pull-down menu for *Oversaturated pixels* and/or *Undersaturated pixels*. See 6.2.3 Channel Coloring.

 **Open Attachment** Opens a file previously attached and saved with the ND2 file. To attach any file to an image, right-click inside the image, select *Add Attachment* and choose a file to be attached. Maximum size of the attached file is 64 MB. Context menu over the image also enables to *Open Attachment...* (same functionality as the  button), *Save Attachment As...* (saves the attached file separately) or *Remove Attachment* (attachment is removed from the ND2 file).
























 **Switch to Main View**,  **Show Slices View**,  **Show Volume View**,  **Show Tiled View**,  **Show Maximum Intensity Projection**,  **Show Maximum Intensity Projection** When working with the ND images, these functions become available.

For more information see 6.3.2 ND Views.

**Enable Dynamic Threshold** Activates the dynamic threshold in the image.

 **Mouse XY** Requires a motorized stage. If turned on, the mouse cursor on a live image changes and you can control the XY stage by dragging the live image.

 **Create Focused Image** Creates an EDF focused image.

-  **Create EDF Focused Document** Creates a new simple focused image from an ND file.
-  **Show EDF Z-profile** Opens the EDF Z-profile panel with the Z-profile graph of the image. See *View > Analysis Controls > EDF Z-Profile* .
-  **Show EDF Z-map** Creates a new image displaying the EDF Z-map. Opens the Z-map setting window, where you can choose the Z-map method. Press OK and the new focused image is created. See *Applications > EDF > Show Z-Map Image*.
-  **Show EDF 3D Surface View** Opens the 3D surface view of the image. See *Applications > EDF > Show Surface View* .
-  **Show EDF Anaglyph** Enables/disables Anaglyph 3D view. New options appear in the toolbar - you can define the position of the first frame (Lowest/Highest), *Style* (the color combination used for anaglyph image), *Effect* (Descend into the Screen / Rise from the Screen) and *Z-zoom* (define the percentage).
-  **Refresh ND View** Refreshes the ND view.
-  **Fit to screen** Adjusts zoom to view the whole image within the *NIS-Elements* screen.
-  **Best Fit** Adjusts zoom to fit the image window in one direction and fill the available area with the image.
- 1:1 **1:1 Zoom** Adjust zoom so that one pixel of the image matches one pixel of monitor.
-  **Zoom In** Increases magnification of the image.
-  **Zoom Out** Decreases magnification of the image.
-  **Show Probe** This button activates the probe. The probe affects histograms, auto exposure and auto white balance functions.
-  **Show Grid** Displays the grid for indicative measurements.
-  **Show Scale** Displays the image scale.
-  **Show Frame** Displays and applies the measurement frame. (*requires: MEAS (Automatic Measurement)*)
-  **Turn ROI On/Off** Displays the Measurement Region Of Interest. (*requires: MEAS (Automatic Measurement)*)
-  **Show Profile** Displays the *Measure > Intensity Profile*  control panel. It allows you to specify a linear section in the image of which the pixel intensities graph will be created.
-  **View LUT Intensity** Displays the scale of intensities used inside the image. It works on monochromatic images or a single image channel.
- A Show Annotations (A)** Displays the vector layer which typically consists of annotation objects (text labels, arrows) and measurement objects.
-  **View Binary** Displays the binary layer of the image. (*requires: MEAS (Automatic Measurement)*)
-  **View Color** Displays the color layer of the image.
-  **View Overlay** Displays the color layer and the binary layer in overlay. (*requires: MEAS (Automatic Measurement)*)

---

**Tip**

Right click the icons to invoke a context menu where properties of each tool can be modified.

## Channel Tabs

Figure 3.7. Channel tabs of an RGB image




Channel tabs at the bottom left corner of the image window enable switching between image channels. You can also edit their properties using commands available via a context-menu. See also 6.2 Image Layers.

## Status Bar

The status bar at the bottom of the image window displays the following information:

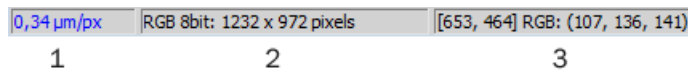Figure 3.8. Status bar of the image window



1. The first field of the image window status bar Image displays the calibration. See also 8.1.1 Calibration, 8.1.1.3 Units.
2. Image bit depth (8bit, 12bit, 16bit, etc.) followed by Image size. You can change the displayed units from the context menu.
3. Pixel coordinates of the mouse cursor along with channel intensities, Binary layer value (0 or 1) and the Color mode (RGB, Monochromatic, etc.).

### 3.1.3 Arranging User Interface

Having a well organized application layout can help you make the work with *NIS-Elements* very effective. There are the following options on customizing the appearance of *NIS-Elements*:

**Custom window placement** All control panels (Camera Settings, Measurement, Histogram, LUTs, etc.) can be arranged inside or outside of the main application window.

**Compact window or multiple windows** The control panels as well as toolbars can be floating or docked on sides of the application screen.

**Multiple monitor support** The *NIS-Elements* window can be stretched to occupy two monitors. When you switch from different application, *NIS-Elements* will be activated on both monitors.

**Customized toolbars** Toolbar buttons may be added and removed from toolbars. See 3.3 Modifying Toolbars.

**Maximizing the Image Area** You can hide some of the GUI elements which are displayed by default:

- The channel tabs and the layout tabs may be hidden to save some screen-space. Display the 3.6 Appearance Options window and de-select the *Show Channel Tabs* and the *Show Layout Tabs* options.
- Image controls and the image status bar may be hidden. Use the *Common for all images, auto hide bottom toolbar* option of the 3.6 Appearance Options window.
- When an image is displayed in great magnification, scroll bars automatically appear by the sides of the image window. You can hide them by de-selecting the *Show Scrollbars* context menu option.

## 3.2 Modifying Layouts

A layout in the context of *NIS-Elements* is a set of options describing the arrangement of control panels, tool bars and menu items.


### To Create a New Layout

1. Select the *Advanced* layout.
2. Right click the tab and select *Save Current Layout As*.
3. Write the new layout name and confirm it by .
4. A new layout is created.
5. Modify the layout so that it suits your concept of work. You can add or remove control panels, tool bar buttons etc.
6. Right-click the tab of the new layout and select *Save* from the context menu.

### 3.2.1 Layout Manager

Run *View > Layout > Layout Manager* to display the Layout Manager. The list of currently available layouts is placed on the left side of the layout manager. Each layout may contain information about controls, toolbars, menu, and commands to be performed when switching between layouts.

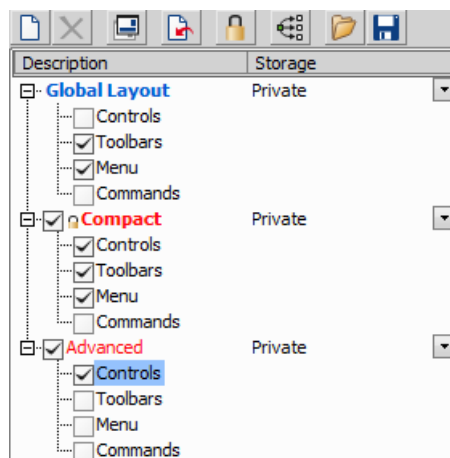Figure 3.9. List of Layouts



### Modifying the Layout Settings

1. Select items within the *Global Layout* list which you want to be shared by all layouts.

2. The check marks on the left of the layout names indicate the layout visibility. Select the ones you want to display in the application status bar.
3. If an item within some particular layout is not selected, it means you do not want to customize it and the global settings - if selected within the *Global Layout* - will be used. If the item is not selected within the Global Layout either, the settings of the most recently active layout will be used.
4. Set whether the layout is *Private* or *Shared*. If set *Private*, it will not be visible for other users. See 2.3 User Rights.
5. Customize each item according to your needs (see below).

---

#### Note

*When upgrading to NIS-Elements version 5.30, only the modified layouts from the older version are loaded into the new Layout Manager.*

### **Making a Layout Default for All Users**

1. Create the layout you would like to propagate to all *NIS-Elements* users.
2. Right-click the command name and select the *Propagate this layout to* command (you can choose either to All Users, or to a specific Group).
3. Restart *NIS-Elements*.

---

#### Note

*No apparent change will be visible in the Layout Manager after propagating the layout. However, after the software is restarted, the layout will be present.*

---

#### Caution

*Users who have the privilege to Modify shared layouts can remove the layout via a context-menu, but this action will also affect all users!*


### **Layout Manager Tools**


**Propagate this layout to all users** This context-menu command makes the layout default to all users of the software. Whenever the software is started the layout will be activated.


 **Create Layout** Adds a new layout to the list of layouts.


 **Remove Layout** Deletes the selected layout. The first two layouts cannot be deleted.


 **Set Layout as Active** Makes the selected layout active.

 **Load Default** Loads original settings of the selected pre-defined layout (*Full Screen*, *Docked Controls*, *Measurement*) so that it looks just like after the program installation.

 **Lock Layout** Select a layout component and click this button to lock these elements so that they cannot be moved or closed. To unlock these components, click the button again or right-click the locked layout tab at the bottom and click *Unlock Layout*.

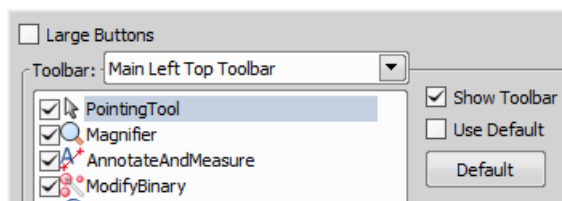
 **Apply to All** Changes made on the currently edited layout are applied to all other layouts using this button. Global layout is not affected however changes made on the global layout are applied to other layouts.

 **Import Layouts** Enables to load a previously saved set of layouts from an XML file. When you try to import a layout with already existing name, you will be prompted to choose whether: the imported layout replaces the existing colliding layout, or is not imported, or is imported and renamed.

 **Export Layouts** The settings of layouts can be saved to an external XML file. Use this *Export* button. In the windows that appears, define the destination file name and check which layouts will be included in the exported xml file.

## 3.3 Modifying Toolbars


### Hiding Toolbar Buttons



1. Display Layout Manager by the *View > Layout > Layout Manager* command.
2. Select the *Toolbars* item within the list of layouts. The right-side part of the window changes.
3. Choose one of the toolbars which you would like to modify from the *Toolbar* pull-down menu.
4. If selected, de-select the *Use Default* option on the right.
5. Any button of the toolbar may be hidden by de-selecting it. No buttons can be added to any of the toolbars except the *Main Left Toolbar*.
6. The whole toolbar can be hidden by de-selecting the *Show Toolbar* check box.
7. There are two sizes of buttons available. Select the *Large Buttons* option to use the larger one. This setting is shared by all toolbars.

### Adding Buttons to the Left Toolbar

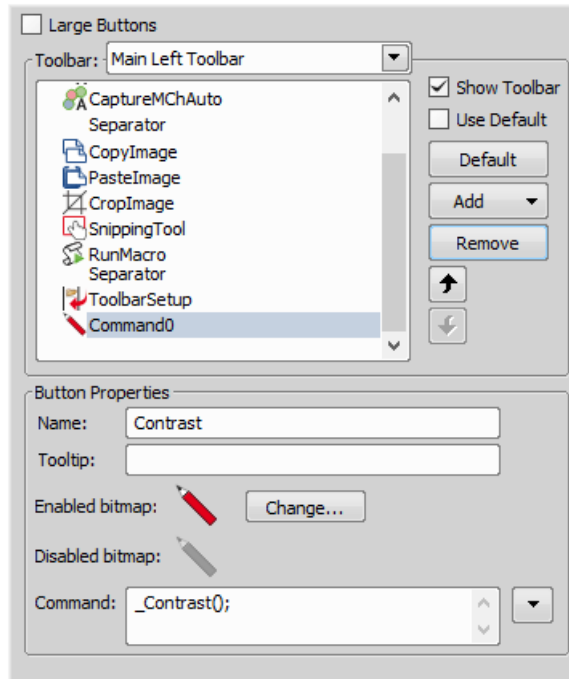Custom user buttons can be added to the main left toolbar. You can define your own buttons which run single macro functions or execute macros. Select the *Main Left Toolbar* from the pull-down menu.



Let's say that we very often use the *Image > Contrast*  command. It is useful to add a shortcut button to the toolbar.

1. Press the *Add* button, and choose *Command* from the pull-down menu:
2. A new command - *Command0* - is added to the list.
3. Now, assign a macro command: Open the pull-down menu on the right side of the *Command* edit box and click *Command List*.
4. A list of commands appears. Choose `_Contrast()`.
5. Confirm the selection by .

#### Note

*It is possible to assign a sequence of commands to a single button by repeating this procedure.*

If you are not satisfied with the default icon, you can change it by pressing the  button. A window for selecting the icon appears. You can select the image from the *NIS-Elements* icon set or load some other from any file containing icons (ico, dll). You can define another icon for the command in a disabled state, too.

It is handy to define a tool-tip (a text that appears when the mouse cursor is placed over the icon) for your command. Simply write the text into the *Tooltip* box. You can change the position of the command in the toolbar using the arrow buttons. The  button discards your changes and inserts the default set of commands to the toolbar.

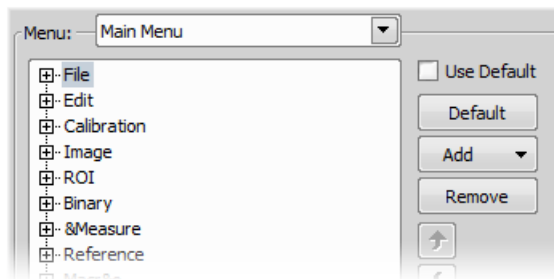


## 3.4 Modifying Menu


The *Main Menu* and some of the *Context Menus* within the application window may be modified. Items of the context menus can be hidden by de-selecting them similarly to the toolbars. The main menu can be modified as follows:

### Modifying the Main Menu

1. Display Layout Manager by the *View > Layout > Layout Manager* command.
2. Select the *Main Menu* in the topmost pull-down menu.



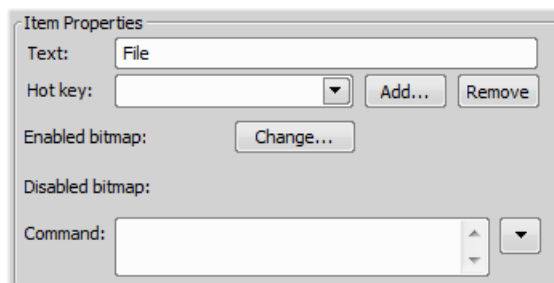3. Any item may be added to the main menu - a *Separator*, a *Menu Command*, a sub-menu (*Menu Popup*), and even a new menu (*Main Menu Popup*) - by the *Add* button.
4. Select the existing menu item under which you would like to place the new item.
5. Click the *Add* button and select the item to be added from the pull-down menu.
6. Edit the *Item Properties*.



**Text** This is the text which appears in the pull-down menu. You may add “&” before any letter - such letter will be considered a keyboard shortcut when browsing the menu.

**Hot key** One or more hot key shortcuts may be assigned to the command. Just press the *Add* button and press the key combination. Press *Remove* to remove the selected hot key.

**Enabled/Disabled bitmap, Command** These fields serve for assigning a bitmap image and a macro function to the menu command. It works the same way as when modifying the Main Left Toolbar (described above).


## Note

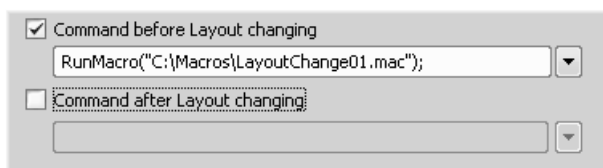The *Default* button discards all changes and loads the main menu original configuration. The *Remove* button deletes the selected item. The arrow buttons move the selected item up/down. The *Use Default* check box, when selected, applies the default settings to the menu.

## 3.5 Running a Macro Upon Layout Change


A macro command or a macro can be run upon layout change.

1. Display the layout manager by the *View > Layout > Layout Manager* command.
2. Select the layout by displaying of which the macro function will run and select the *Commands* check box.
3. The following box appears on the right side:



4. Select the timing. The “before” option will run the command when you click on the layout tab, but before actually changing the layout. The “after” option runs the command right after the layout is changed.
5. When the field is enabled, type the command or insert it via the pull-down menu.

## 3.6 Appearance Options

General appearance adjustments can be made in the *Options* window. Run the *Edit > Options*  command and switch to the *Appearance* tab.

**Background** Background of the main screen may use default tiles or custom color.

**Color scheme** There are the following color schemes predefined within the application: Light scheme, Dark scheme, Black scheme.

**Language** Select the language to be used in GUI. The language pack is a part of the *Local Option* feature set.

**Prompts on Image Save dialog** Actual words used on buttons of the window which appears if an image has been modified and is about to be closed.

**The Z value displayed in status bar** Having two Z devices, select whether to display Z1, Z2 or both on the main status bar.

**Close ND Acquisition window after Run** The *View > Acquisition Controls > Capture Timelapse* window will be closed automatically upon acquisition start.

**Image window toolbars** You can decide whether the top and side image toolbar is integrated into the main toolbar (*Common for all Images*) or whether each image has its own set of toolbars (*Visible for each image*). The most space-saving option is to use shared toolbars and hide the image status bar and the ND2 control bar (6.3.1.1

Control Bar) automatically (*Common for all images, auto hide bottom toolbar*). It will show up only when the mouse cursor rolls over the bottom part of the image.

In the *Right Toolbar* section you can select the appearance of buttons having more sub-options. See 3.1.2 Image Window .


**Common for all images, auto hide bottom toolbar** This option hides the image status bar and the ND2 control bar (6.3.1.1 Control Bar) automatically. It is displayed only when the mouse cursor rolls over the bottom part of the image picture.



**Show channel tabs** Show channel names and colors in the image window status bar.

**Show Layout tabs** Show layout tabs in the main status bar.

**Show Task Bar** If unselected, the Windows taskbar will be hidden.

**Default vertical docker on the right/left side** Select the preferred side for the default vertical docking pane.

**Keep text size while zooming** When selected, text annotations will not be zoomed with the image. See *View > Analysis Controls > Annotations and Measurements*  .

**Allow zoom factors lower than best fit** When selected, the image can be resized to be smaller than displayed using the *View > Zoom*  > *Fit to Screen*  command.

**Lock camera magnification** If you are used to switch between two camera modes (resolutions), this option ensures that the scene observed does not change its size or position.

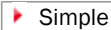

**Keep picture window aspect** If checked, the image window respects the image size while zooming.

**Initial Zoom** You can select the zoom factor of the newly opened images. The options are Best Fit, 200, 100, 50, 25%.

**Initial zoom 100% for Live and Capture** Whenever a new Live or Captured window is opened, the zoom will be automatically set to 100%.

**Default Volume View Rendering Engine** Switch between the Old/New engine used in the Volume Viewer.

## 3.7 Simplified User Interface

After the *Industrial GUI* module is installed, a simple layout intended to be used in industry imaging applications is added to the application. Switch to the layout by the  **Simple** button in the top right corner. Return to the main layout by the same button (which is changed to  **NIS-D**).

### Note

*This button corresponds to the View > Simple Control menu choice.*

*A Nikon-manufactured digital camera (DS-U1/U2/U3, DXM1200C, 1QM) must be connected to NIS-Elements in order to display the View > Simple Control command.*



### 3.7.1 Behavior


- The “Live” and “Captured” images can be switched between the main image area and the preview window on the right.
- The number of control panels available within the docking panes is reduced.


- The main toolbar is reduced, some buttons are added. See 3.7.2 Toolbar Buttons.


### 3.7.2 Toolbar Buttons

The following toolbar buttons are added to the layout:




 **Print** This button sends the current image (the large one) to the default printer. Invoke the nearby pull-down menu to set printing properties. See *File > Print* .

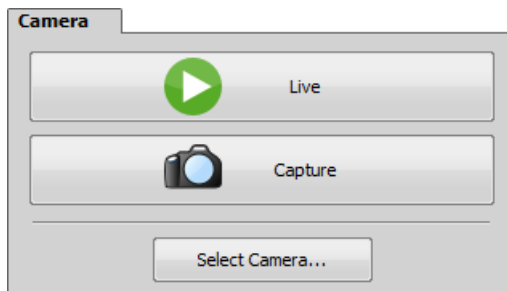
 **Change View** This button swaps the Live and the Captured images between the main image window and the preview window in the right-side docking pane.

 **Show Graticule** This button has been moved here from the right image toolbar of the main view and provides the same functionality. See 3.1.1 Main Window Components.

 **Show Scale** This button has been moved here from the right image toolbar of the main view and provides the same functionality. See 3.1.1 Main Window Components.


### 3.7.3 Camera Control Panel

The  *Live* button displays live image in the picture window. The second button either captures a single image after clicking the  *Capture* or captures a single image and saves it to the *Auto Capture Folder* automatically after clicking  *Auto Capture* (check the *Auto Capture (Auto Save)* option in the *Save* control panel).




### 3.7.4 Save Control Panel

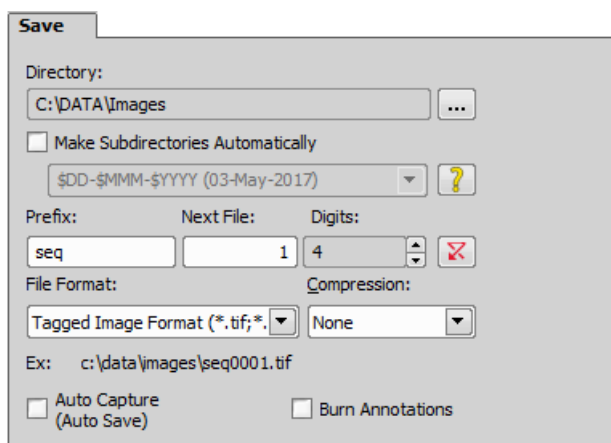This control panel enables the user to specify image saving options. Most of the settings only applies when the *Auto Capture (Auto Save)* option is checked.



**Directory** Press the ... button and select the *Auto Capture Folder* where images captured by the  **Auto Capture** button will be saved.

**Make Subdirectories Automatically** The images can be sorted to subdirectories automatically based on the time of acquisition. Select this option and use place-holders from the drop-down menu to specify the right format (click  to display the place-holder legend). E.g.: “\$YYYY\$MM\$DD” would create a new subdirectory for each day named like this: “20150103”.

**Prefix, Next File, Digits** Type the prefix and numbering style intended for the files being saved automatically. The *Next File* field 1/ informs you and enables you to change the number which will be used for the next automatically saved image.

**File Format, Compression** Select the file format to save the images in. Some formats enable you to set the *Compression* parameter. It is recommended to use either “none” or “lossless” in order to preserve good image quality.

**Ex.** The path and file name of the next image to be saved is displayed here.


**Auto Capture (Auto Save)** Check this option and a new image will be created and saved to the *Auto Capture Folder* each time you press the *Auto Capture* button.


**Burn Annotations** This option determines whether the annotations created by the tools of the *Annotations and Measurements* window will be included in the saved file.

#### Warning

*The separate annotation layer will be merged with the image data.*

### 3.7.5 Preview Control Panel

The Live or Captured image is displayed within this window. Use the  **Change View** button of the main toolbar to select which one.

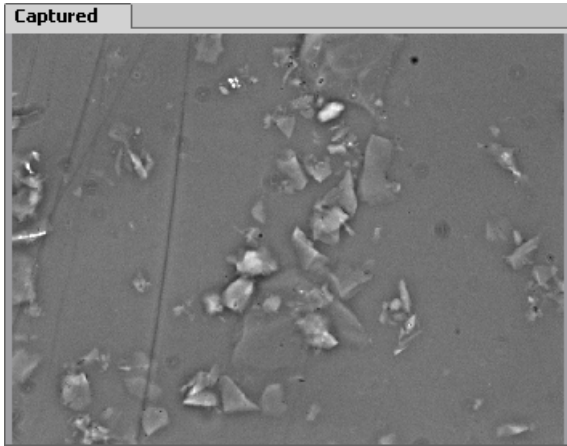



## 4. Cameras & Devices

### 4.1 Basic Workflows


#### 4.1.1 Startup Hardware Configuration Selection

If you are running *NIS-Elements* for the first time, you will be asked to select a microscope and define the first hardware configuration. Click **Select Microscope** to open the *New Setup* window. Select the microscope connected to your computer and click **Create Setup**. The *Device Manager* window opens. Use this manager to add all cameras and devices connected to the microscope, then create hardware configurations and light paths. Please see 4.2 Visual Device Manager .

You will be asked to select a hardware configuration every time you launch *NIS-Elements*. Only the cameras and devices present in the selected configuration are connected. You can switch to another hardware configuration later in *Devices > Device Manager*  by switching the tabs or using the *Devices > Select HW Configuration* command. If you need to create a new hardware configuration without loading an existing one, please select *Offline Analysis* during the startup of *NIS-Elements* and then create a new configuration in the *Devices > Device Manager*  (see 4.2 Visual Device Manager ).



#### 4.1.2 Dual, Triple, Quad Camera Modes

Some camera models may be used in a dual, triple or quad mode. The software combines their grayscale images into one multi-channel image.

1. Run the *Devices > Device Manager*  command.
2. Add the *Dualcamera* device and two or more cameras to the configuration. Arrange them to match the physical setup.
3. Right-click a camera and select the *Make Dual Camera with* command. A dashed line will connect the two selected cameras.
4. Repeat the previous step for all of your cameras. For a triple camera setup use two *Dualcamera* devices connected in a row, for a quad setup use three.

#### 4.1.3 Using a *Multiview* Device

A *Multiview* device (dual, triple or quad) enables you to divide the camera chip into halves/thirds/quarters and observe the chip parts in an overlay as a multichannel image.

1. Run the *Devices > Device Manager*  command
2. Add your multi-view device to the configuration by the **Devices**  button. If you cannot find your multi view brand in the list, use the generic *MultiView* device.
3. Arrange the configuration to match the physical setup.

4. Right-click the multi view device and select *Configure Multiview* and select whether it is a *Dual View*, *Triple View* or *Quad View*.

5. Click .

#### Note


- *Bypassing the multi view feature is done by switching to the light path with the “MV(Out)” postfix.*
- *To run a multi view mode on a dual camera, connect two “Multiview” devices to one “Dualcamera” device, display a context menu over a camera and select *Make Dual Camera with*.*

## 4.1.4 Optical Configurations

### 4.1.4.1 Introduction to Optical Configurations

Typically, a laboratory computer image analysis system consists of a computer, a camera, and a microscope equipped with certain accessories (objectives, filters, shutters, illumination, rotary changers, etc). Most of the mentioned microscopic hardware can be motorized and therefore can be controlled via *NIS-Elements*. In addition, it is possible to integrate single settings of all these devices into one compact set called *Optical Configuration*. It is recommended to create several optical configurations containing particular devices settings. Then a single click can completely change the current hardware configuration.

### 4.1.4.2 Creating New Optical Configuration

- 1) Please check that all the devices (microscopes, cameras, etc.) which you want to associate with the new optical configuration are properly attached to the system and working.
- 2) Choose the *Calibration > New Optical Configuration*  command. In the window which appears, adjust the settings of the devices to match the intended state which will be saved to the optical configuration.
- 3) Type the name of the new optical configuration to the *Name* field. Use a short descriptive name, the name is used on the button in the main toolbar when you select the *Show on toolbar* option.
- 4) In the left column, select which device settings to associate with the Optical configuration. Only the devices related to the currently selected light path are listed here.

**Camera setting** A list of the current camera properties appears on the right. It is being updated dynamically.

**Use Stored ROI** If you want to include the camera ROI settings, select this option. Status of the camera ROI (ON/OFF) and its size will be set when switching to this configuration.

**Channel setup** These settings determine how channels of newly captured images will be named and what color will be assigned to them. The available properties depends on the current camera setup (color/mono). In the mono camera mode, you can either assign the name, the emission length and color to the channel *Manually*, or leave this task to *NIS-Elements* (the *Automatically* option). In such case, information of the light path (emission wavelength) will be used to determine the channel name and color.



**Microscope setting** If there is more than one shutter available and you would like to associate a shutter with the optical configuration, select which one is the *Active Shutter* from the pull-down menu. Select which parts of the microscope shall be included in the configuration by checking them in the *Used devices* dialog box.


Note

*Active shutters remember their aperture setting and display it in the Microscope setting section.*


**Objective** An objective mounted to a motorized nosepiece can be included in the configuration. Select the objective from the pull-down menu. Objectives which are currently assigned to any position of the nosepiece are listed. See 4.1.5 Objectives.

Note

*The objective must be assigned to a nosepiece position via the microscope control pad or the nosepiece control panel beforehand.*

- 5) If some of the device settings still need to be adjusted, click the **Camera & Devices Controls** button and select the appropriate control panel from the pull-down menu. Adjust the settings within the control panel, the optical configuration will be updated automatically.
- 6) Click *Finish* to save the new optical configuration and to close the window.
- 7) You can create more optical configurations by repeating the procedure. The optical configurations are saved to registry immediately. A backup of optical configurations can be made by running the *Calibration > Optical Configurations*  command and clicking the *Export* button.

#### 4.1.4.3 Managing Optical Configurations

To display the optical configurations manager window, run the *Calibration > Optical Configurations*  command. You can make the following actions from the window:

- Create, duplicate, rename, delete, copy settings and switch between optical configurations.
- Modify optical configuration properties.
- Import and export optical configurations to/from an XML file.

#### 4.1.4.4 Operations with Optical Configurations

Once created, the configuration appears in the list and can be shared with other users by changing the *Private* option to *Shared* (see 2.3 User Rights). The following operations can be performed on the selected configuration:

- It can be deleted by pressing the *Remove* button. A confirmation dialog box appears.
- Its name can be changed by the *Rename* button. A button with the configuration name appears in the toolbar (if the *Show on toolbar* option is selected).
- The configuration settings may be transferred to another optical configuration by the **Copy to** button. Press the button and select the optical configuration to be overwritten with the current one.
- A copy of the configuration can be made via the *Duplicate* button.

- The configuration can be applied to by the *Set As Active* button.
- The settings of all optical configurations can be exported to an external XML file using the *Export* button.
- The previously exported optical configurations settings can be loaded from the XML file via the *Import* button.
- The list of optical configurations can be ordered manually using the arrow-up and arrow-down buttons.
- Each configuration can be arbitrarily modified within the right-side portion of the window.

The optical configuration buttons are available in the main toolbar when *Show on toolbar* option was selected during the setting process.

#### 4.1.4.5 Using Auto Exposure while Switching between Optical Configurations

You can use one optical configuration when working with live image and switch to another optical configuration just for the image acquisition, very often, this is performed automatically. However, if the optical configuration used for acquisition has the *Auto Exposure* camera mode turned on, it may produce over-saturated or under-saturated images.

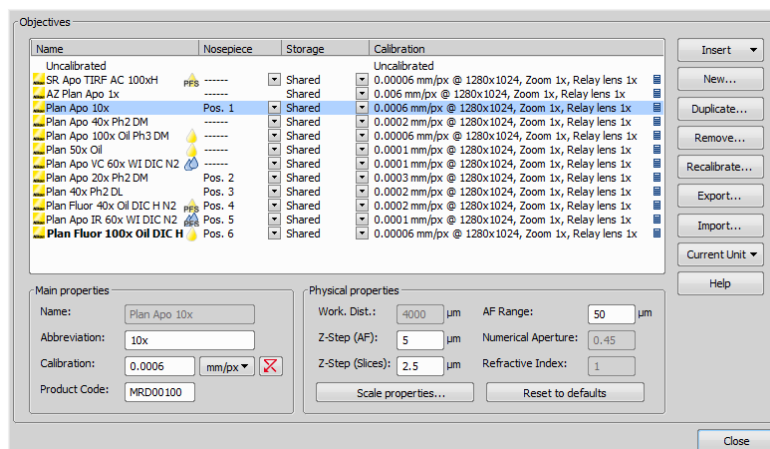
Since the Auto Exposure algorithm estimates the optimal exposure time by analyzing several last frames, it is not able to make the estimation correctly if the change of optical configurations would cause significant change in image brightness. To prevent this, we recommend to use *Manual Exposure* mode in the optical configuration used for capturing.


### 4.1.5 Objectives

If you want to perform measurements on captured images it is wise to calibrate all objectives used to capture images. Whenever an image is captured through a calibrated objective, the image inherits its calibration.

#### 4.1.5.1 Managing Objectives

Run the *Calibration > Objectives* command. The *Objectives* window appears:



A list of used objectives can be created within the Objectives window. For each objective, the objective name, position in the changer, storage status and the calibration are displayed. A *Calculator*  icon next to the calibration value indicates that the calibration has been calculated from objective properties. If the calculator icon is missing, the calibration has been performed manually. Use the buttons on the right side to manage the objectives:

**Insert** Click the button and select one of the objectives from the database (ordered by magnification). Custom objectives can be added to the database via an INI file.

**New** Press this button to create a custom objective. The new objective will be added to the list. Then, define the *Main properties* and the *Physical properties* of the objective.

**Duplicate** If zoom is used, the objective calibration must be re-calculated accordingly. Use this button to make a copy of the objective and define the zoom factor in the window which opens.

**Remove** This button deletes the selected objective.

**Recalibrate** Starts the calibration of the selected objective. See 4.1.5.3 Objective Calibration.

**Export** Enables you to export the complete list of objectives to an XML file. The standard Save As window appears.



**Import** This button enables you to import a complete list of objectives from an external XML file previously created by the *Export* button.

**Current Unit** This button invokes a pull-down menu where units for the whole application can be selected.

**Help** Displays a help page to the *Calibration > Objectives* command.


Edit the properties of the selected objective in the bottom part of the window. Press the *Close* button to finish this window.

#### 4.1.5.2 Assigning Objectives to Nosepiece Positions

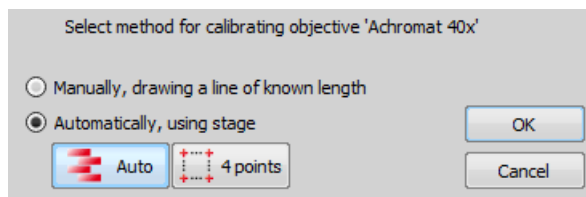1. Display the microscope control pad - *View > Acquisition Controls > Microscope Control Pad* 
2. Click the setup  button in the *nosepiece* section.
3. A window appears which enables you to select one of the available objectives to the corresponding position.
4. Each objective has its specifications displayed in the table. These specifications are not editable.

#### 4.1.5.3 Objective Calibration

After you press the *Recalibrate* button in the *Objectives* window or when creating a new custom objective, the following window appears:



Select one of the calibration methods:

- The *Manual* calibration lets you draw a distance into a picture and assign real length to it. The live image starts automatically and a window appears. Proceed to: 4.1.5.4 Manual Calibration
- If a motorized XY stage is available, the *Auto* and *4 points* buttons appear. Proceed to: 4.1.5.5 Objective Calibration Using XY stage.

Press **OK** to continue.

### **Superresolution Calibration**

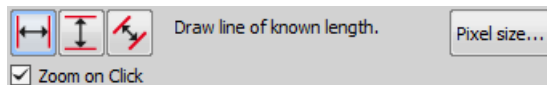
This type of calibration is available after successful manual or automatic calibration. It triples the camera's resolution and uses fine stage movements to improve the resulting objective calibration. Click the **Run SR Calibration** button and wait till the objective is calibrated.

#### 4.1.5.4 Manual Calibration

Manual calibration requires the user to specify a distance in the image. It can be either a measurable distance or the size of one pixel. The following procedure can be used for different purposes:


- Calibrating of an image (see 8.1.1 Calibration)
- (Re)calibrating of an objective (see 4.1.5.3 Objective Calibration)
- (Re)calibrating of a microscope zoom

1) The following window appears:






2)


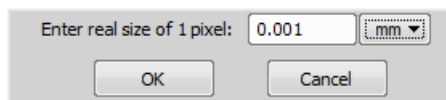
a) If you know the size of one pixel, click the **Pixel size...** button.




Enter the value and finish the calibration by the **OK** button.

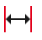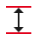b) If you do not know the pixel size, you will need to specify distance in the image. For this purpose, use an image of a calibration slide or a ruler captured with your current system configuration.

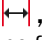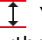
- 3) The distance is defined by placing lines to the image. If you are sure the calibration pattern is perpendicular to image edges, it is recommended to select either vertical  or horizontal  lines. Otherwise, select the parallel lines .
- 4) Click into the image to place the first line and then the second line to specify the distance:

,  You can modify the line position while holding the primary mouse button. After you release the button no further changes can be made. Place the second line in the intended position by another click.

 Draw the first line by clicking twice inside the image. The line can be moved and adjusted arbitrary by mouse. When satisfied, finish the first line by right-click. The second line can be placed by another click to the image, this time only to adjust the distance from the first line. The process is completed by right-click.

*Zoom on Click* zooms the image after the mouse button is clicked so that the position of the point can be set precisely.

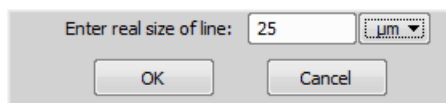- 5) Then the following dialog box appears:




- 6) Enter the distance between the two lines and select correct units.
- 7) Click **OK** to finish the calibration.

#### 4.1.5.5 Objective Calibration Using XY stage

This procedure requires a motorized stage. Select one of the following methods:


**Auto** The *Auto* method is fully automatic. Calibration is performed on a part of the live image marked by red square that is shown before the calibration is launched. You can also choose the channel to be used by auto-calibration.

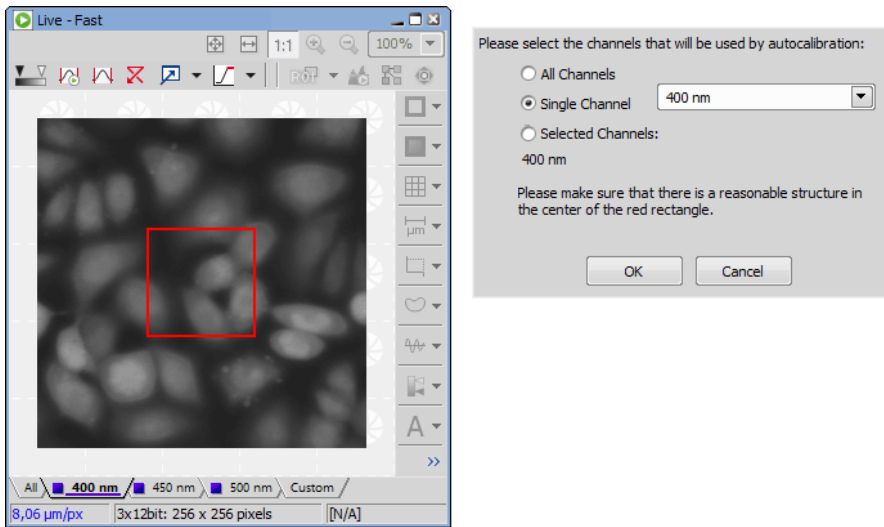Figure 4.6. Auto-calibration window



NIS-Elements moves the motorized stage, acquires two images, and calculates the calibration from the shift of the images. The auto-calibration performs an estimation of the real relay lens zoom factor. If the estimation does not match the value specified in the *Device Manager* window, a warning appears.

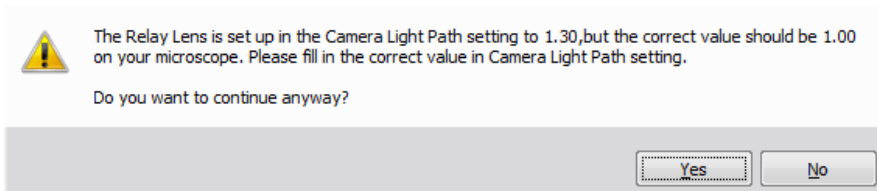Figure 4.7. Example of the warning message



It is recommended to cancel the auto calibration (click **No**), check the real zoom factor of your relay lens and set correct relay lens settings in the *Device Manager*.

#### Note

The success of this method depends on the texture of the specimen, its contrast, illumination, etc. If the combination of these factors is unsuitable, the auto calibration may fail. If it fails, try the following:

- Move the stage to another area of the specimen to get a better texture.
- Improve contrast by setting LUTs.
- Check focus, re-focus if needed.
- Turn on the shading correction (see 5.3 Shading Correction).


**4 points** If you select the *4 points* method, the system draws four points on the screen (subsequently) and asks user to move one significant part of the specimen to each position. After all four steps are completed, the calibration is calculated from the moves of the stage.

#### 4.1.5.6 Monitor Calibration

A real zoom factor can be displayed within the image window. In order for the value to be accurate, both the image calibration and the monitor calibration must be accurate. To calibrate the monitor:

1. Display a context menu over a calibrated image and select the *Show Physical Zoom* command. The computed value appears in the top left corner of the image. If the current image is not calibrated, the box will show *N/A*.
2. Right-click the box with the zoom value and select *Monitor Calibration*, the following window appears.

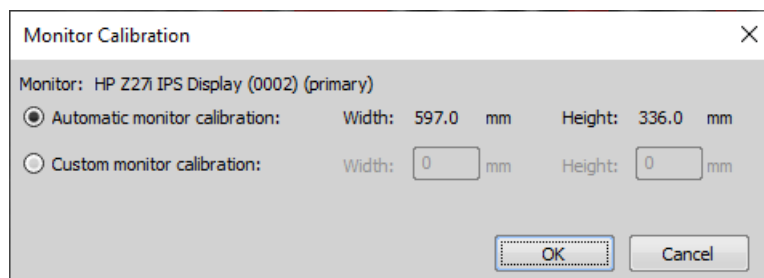Figure 4.8. Monitor Calibration Dialog



#### Caution

*The dialog displays calibration of the monitor on which it is displayed. If you have multiple monitors, move the dialog between them.*

3. Check the *Width* and *Height* values and whether they match the real size of your monitor.

#### Tip


*The detected monitor type is displayed in the window. If the information about the model is precise, it is likely that the automatic calibration is correct, otherwise, we recommend to switch to Custom monitor calibration.*

4. If you are not sure whether the detected calibration is correct, select *Custom monitor calibration*. Take a ruler or a measuring tape and measure width and height of your monitor. Enter these values and confirm the calibration by the **OK** button.

#### 4.1.6 Connecting a Device to NIS-Elements

Before you get to work with *NIS-Elements*, all hardware accessories should be connected properly to the system. In most cases, the following basic procedure is sufficient to connect a device successfully:

1. Install *NIS-Elements*, and select the appropriate device(s) during the installation.
2. Connect the device to the PC and switch the device ON.

3. Run *NIS-Elements*, and run the *Devices > Device Manager*  command to open the *Device Manager* window.
4. Select a microscope or create a new one.
5. Click the *Add* button next to the *Devices* caption to add the device to *NIS-Elements*.
6. Select the device from the list of installed devices and click Add Device.
7. Click on the device in the scheme and drag it to the highlighted hotspot so that the connection matches your physical microscope setup.
8. Create a hardware configuration with light paths suitable for the experiment.
9. Close the *Device Manager*.

If a peripheral device is connected to *NIS-Elements*, some extra commands and features become available. This concerns typically motorized XY stages and Z drives, but also other devices.

---

#### Note

*MULTIZOOM AZ100M, ECLIPSE LV series, ECLIPSE MA200, or ECLIPSE L200N/300N microscopes can not be connected from NIS-Elements while the setup tool of each microscope is active. To connect any of these microscopes to NIS-Elements, exit the setup tool and then run NIS-Elements.*

---

#### Renaming of devices

*Either logical or physical device can be renamed by the user. Just right-click the device name and select the *Rename...* command from the context menu. The user-defined names can be handy in two cases:*

- *You are using two logical devices having matching names, but you need to uniquely identify them (e.g. from macro).*
- *You are used to call a device with another name and would like to rename it in order not to be confused by the predefined name any more.*

### 4.1.7 What are “Logical Devices”?

*NIS-Elements* handles hardware accessories using the concept of logical devices. There are features of different hardware devices which equal and therefore can be controlled equally. Such features are called “logical devices”. A typical logical device is *Stage XY*. Different microscopes can be equipped with different XY motorized stages, although - regarding the user interface - they behave equally. One physical device (a piece of hardware) can contain one or more logical devices the list of which appear in the *Device Manager* after the connection is established.

#### **Available Logical Devices**





**Analyzer** The analyzer is a polarizing filter placed in the optical path between the specimen and the lamp. The logical device offers two states: ON (inserted) and OFF (extracted).

**Aperture** This logical device is used for controlling apertures in the light path. It is used in complex microscopes rather than as a standalone device. Two parameters can be typically set for aperture devices, the state (on/off) and aperture size.

**Condenser** A condenser is a two-lens combination located next to the illumination source in the optical path. Its purpose is to collect light and direct it to the specimen being examined. The corresponding logical device relates to a changer of different condensers.



**Filter** This logical device controls filter changer movements. There can be several filter changers connected to *NIS-Elements* at a time. Each filter changer needs to be set up - filter types shall be assigned to positions of the changer:

1. Display the filter changer control panel (*Devices > Filters and Shutters* or *Devices > Microscope Control Pad* ).
2. Click the settings button , a window appears.
3. Select one of the available positions which the filter will be assigned to.
4. Click the  button, a list of available filters appears.
5. Select the filter name from the list and confirm it by .
6. The filters can be moved within the already defined positions using the *Up/Down* arrow buttons.


---

Note

*When browsing the list of filters, details about the currently selected filter are displayed on the right side of the window.*

**Illuminator** This logical device is used for controlling the specimen illumination remotely. There is no standard dialog box for the illuminator control. Each device handles this logical device via a user interface specially designed for it - typically containing one button for switching it on/off and a slider for regulating intensity.

**Light Path** Some microscopes have more than one port where it is possible to attach a light source or a camera. This logical device can switch the illumination between these ports.

**Microscope** This logical device is used to group standalone logical devices used in certain microscopes. To control the logical devices of a microscope from one control panel, select the *Devices > Microscope Control Pad*  command.

**Nosepiece** This logical device serves for controlling microscope objective changers. There can be three nosepiece types attached to a microscope:

- *Manual* - it can not be controlled via the software.
- *Intelligent* - the current nosepiece position is read by the application, but can not be controlled.
- *Motorized* - such nosepieces can be fully controlled via the *Microscope Control Pad* or the *Nosepiece* control panel.

See 4.1.5.2 Assigning Objectives to Nosepiece Positions.

**ND Filter** A neutral-density filter is a light absorbing filter whose absorption spectrum is moderately flat. It is used to reduce the illumination intensity within the optical path. The logical device offers two states: ON (inserted) and OFF (extracted).

**PFS** Perfect Focus System - this logical device corresponds to the PFS physical device available with Nikon TE2000/TI microscopes.



**Shutter** This logical device can control shutters installed in your system. This device is handled via the *Devices > Filters and Shutters* control panel or straight from the microscope control pad. You can select the type and rename the shutter by running a contextual menu command either within the *Filters and Shutters* control panel, *Device Manager* window, or in the main toolbar.

---

Note

*There was a restricted number of shutters which could be connected to the system in previous releases of NIS-Elements and they were identified by Type (DIA, EPI, Aux1, etc...). The current version supports identification*

of shutters by custom Names, but the Type attribute can be still found in some windows (or used in some macro functions). This is to ensure backward compatibility.

**Zoom** This logical device is used for controlling the zoom factor. Run the *Devices > Zoom Configuration*   command to adjust the zoom settings.

**Stage XY** Stage XY enables the movement of specimen within X and Y axes. The system offers user to control stage movements.

**Stage Z** The Z Drive device enables movement throughout the Z axis direction.

### 4.1.8 XY Stages and Z Drives Tips


Please read the following tips for using motorized XY Stages and Z Drives.

#### **Motorized Stage Initialization**

A motorized stage being initialized could strike the objective. Before initializing the stage, make sure that the objective is away from the stage. If a motorized Z drive is available, use the *Devices > Objective Clearance* command to prevent collision.

#### **Setting Software Limits to Stage Movement**

Some microscopes enable you to reduce the range of movement of the motorized stage by setting limits within the configuration window.

- 1) Display the configuration dialog window (within the *Devices > Device Manager*  window).
- 2) Move the stage to the position where the limit shall be set.
- 3) Click the appropriate button.
- 4) Set all the limits by repeating this procedure.

#### Caution

*This procedure can not be used if you already set some limits and would like to broaden them at the same time. NIS-Elements will not allow you to move the stage to any out-of-limits position. What you need to do is to reset the limits within the configuration window beforehand.*

#### **Using two independent Z Drive devices**

Workstations can be equipped with two independent Z drive systems, one coarse (slow) and the other fine (fast). Typically, the first one is used for specimen manipulation in Z axis, when changing objectives, etc., the second one is used for auto-focusing. The following text explains how *NIS-Elements* handle the two Z drives.

**Absolute Z** The current positions of both Z drives (Z1, Z2) and the absolute Z position (sum of Z1 and Z2) are displayed in the main status bar. If there is not enough space to display all these values, only the *absolute Z* value is displayed (Z1 and Z2 still appear in a tool-tip). The absolute Z value appears in the Z Series Setup.

#### **Active Z**

The concept of *Active Z* enables user to select the preferred Z drive:

- The active Z drive is used when performing auto-focus.

- The *Devices > Enable Mouse Joystick Z in Live* command applies to the active Z. You can select the Active Z device in the *Devices > Mouse Joystick and Auto Focus Z* menu.


### Regular Z drive + Piezo Z drive

If you have a regular Z drive and a Piezo Z drive installed, some functionality is added:

- The *Move Piezo Z to* button is added to the *Z Series Setup* window when Piezo Z is selected.
- The *Devices > Keeps Z position and centers Piezo Z* and the *Devices > Move Piezo Z to Home Position* commands appear in the *Devices* menu

The button and the *Devices > Keeps Z position and centers Piezo Z* command provide corresponding functionality. They move the Piezo Z device to its home position, and compensate this shift by moving the second Z drive in order to maintain the absolute Z position.

## 4.1.9 Camera Settings

A digital camera mounted to a microscope port records images of the observed scene on a light-sensitive sensor, and transfers them to a computer. *NIS-Elements* supports various cameras differing from each other in resolution, frame-rate, sensor type, etc. Despite these differences, controlling different cameras is similar. The following features are included in the *View > Acquisition Controls > [Camera name] Settings*  window depending on a particular camera type.

### Note

*The complete list of cameras and devices supported by NIS-Elements is available in a separate document.*

### Reference of Common Camera Settings

**1 frame for Fast Timelapse** The camera head is equipped with 4 GB of memory available for fast capturing. Use this button to run acquisition with maximum available frame rate. However, the size of image sequence is limited to 4 GB.

### Note

*The actual frame rate depends on other settings (readout mode, sensor mode) and is limited by the read-out speed of the camera sCMOS chip. Maximum frame rate can be achieved with the combination of Rolling shutter and Overlap sensor mode.*

**AE Compensation** In automatic exposure modes, the compensation affects how optimum exposure settings (Exposure time and Gain) are calculated. The Compensation value is expressed in Exposure Values (EV). Setting the compensation to + 1.0 EV makes the image twice brighter (e.g. doubles the Exposure time or Gain).

**AE Lock** This option makes the automatic exposure mode to lock the current exposure settings (Exposure time and Gain).

**Analog Gain** Controls the Gain (Sensitivity) of the camera and strength of the camera analog signal before it is digitized. This setting affects image brightness.

**Auto White Balance** This button performs an automatic white balancing. It calculates the right values and then adjusts the red, green, and blue image components in order to get a neutral white color. The Auto White gives the best results on color neutral (gray) scenes.

**Averaging** Averaging is a commonly used technique of decreasing noise in the image. In this method 2, 4, 8 or 16 consecutive frames are averaged together.

---

**Binning** The binning mode provides considerably enhanced camera chip sensitivity by integrating more elements (pixels) together. E.g.: *binning* 4x4 integrates the signal from the area of 4x4 chip elements to one pixel of the resulting image. Smaller resolutions and faster frame rates are achieved using the binning modes.

**Capture Bit Depth Channel** Select per-channel bit depth used for capturing.

**Clear Cycles** This option enables you to decide, how many times the CCD chip should be cleared (reset). The set *Clear Mode* determines how often it is performed.

**Clear Mode** Sets how often the CCD chip will be reset. Please, see the camera user manual for further details. These options are available:

- Automatic
- Never
- Pre-exposure
- Pre-sequence
- Post-sequence
- Pre and Post-sequence
- Pre-Exposure and Post-Sequence

**Contrast** Affects dynamics of how the luminosity is rendered. There are several modes for different illumination (contrast) scene situation.

**Conversion Gain** This is a hardware pre-amplifier gain.

**Cooling** The Qi1/Ri1 camera heads are equipped with a cooling system that enables to cool down the CCD chip by  $-5^{\circ}$  or  $-10^{\circ}$  of the current room temperature.

---

Warning

*When cooling by  $-10^{\circ}$ , there might be a risk of water condensing inside the camera head (depending on humidity inside the room).*

**Desired Temperature** Sets the target temperature of the CCD chip.

**Detector Mode**

*Only in 12bit mode*

**Timed** A common mode for capturing images.

**Strobe** If the camera controller is equipped with *DS-RC* switch, this mode needs to be selected to use the switch for capturing images.

**Dynamic Range** Dynamic range of a camera sensor is defined by the largest possible signal that it can generate divided by the smallest possible signal. Higher bit dynamic ranges bring wider ranges of grayscale levels from each camera pixel.

**EM Gain Multiplier** EM gain allows weak signals to be multiplied before any readout noise is added by the output amplifier, hence rendering the read noise negligible. There is a complex relationship (approximated to an exponential) between the value set and the actual level of multiplication. Typically, optimal signal-to-noise ratio and dynamic range is achieved between x1 to x300 EM Gain. When the setting exceeds the value of 300, the field gets highlighted red in order to notify you.

**Exposure, Exposure Mode** Exposure mode determines how the Exposure time is calculated. Usually, the user can select between the automatic and the manual mode:

**Auto Exposure** Automatically calculates the exposure time in order to achieve optimal brightness of the scene. Image quality is the priority, therefore longer exposure times are preferred.

Note

*Depending on particular light conditions, the result of auto exposure may not be optimal for image acquisition.*

Warning

*Under- over-saturation may occur when switching between optical configurations. Please see 4.1.4.5 Using Auto Exposure while Switching between Optical Configurations.*

Note

*Available exposure time range depends on the selected Format and Readout Speed.*

**Manual Exposure** The user selects both *Exposure* time and *Gain* manually.

**Continuous AE** When selected, the automatic exposure time is calculated continuously. If brightness of the scene changes, the exposure time is adjusted so that the live image is not over- or under- saturated.

**Exposure Time** *Exposure time* is the time of charge accumulation in a camera chip between two adjacent frames. Prolonging the exposure time increases brightness of the image as well as its quality (there will be less noise).


**Fast (Focus), Quality (Capture)** The software enables to create two resolution presets - *Fast* used for live image and *Quality* used for capturing. The formats then differ in image size, and consequently in frame rate (number of frames per second - fps). The higher the resolution is, the lower frame rate can be achieved. Available resolutions depend on the camera type.

Note

*When switching between the formats while observing the live image, the size of the image on the screen is maintained - the zoom setting is changed instead. Only in some special cases, the behavior changes and the image size is changed instead of the zoom setting which is maintained.*

**FOV Size** FOV of a camera which has the sensor larger than the image from the microscope port (imaging circle) needs to be reduced. Otherwise, images captured with a too-large FOV Size would contain visible vignetting (will be darker at the periphery).

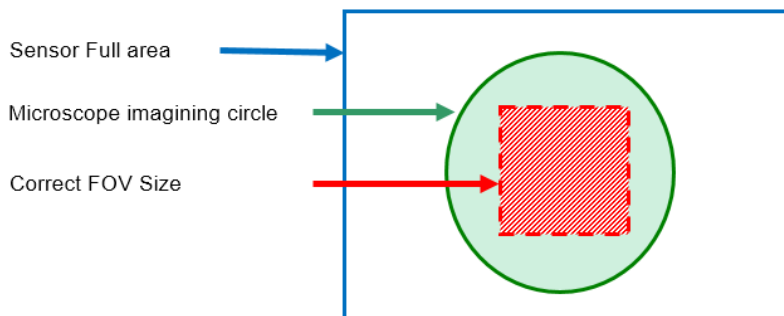Figure 4.9. Correct setting of FOV Size



---

**16 mm** Suits most conventional microscopes

**22 mm** Suits the *Nikon Ti2-E* microscope equipped with two filter-turrets.

**25 mm** Suits the *Nikon Ti2-E* microscope equipped with a single filter-turret.

**Gain** Controls the camera sensitivity. Increasing the gain increases brightness of the image, but decreases quality (the more random noise, the more streaky noise and color unevenness), and increases the frame rate indirectly by enabling shorter exposure times.

**Gamma** The gamma correction maps the intensity of live signal exponentially to the gamma parameter. For gamma < 1, dark portions of the image are enhanced whereas for gamma > 1, image parts of higher intensities are enhanced.

**High Quality Capture** When this option is checked, images are not processed by the camera control unit, but are sent to the computer in a RAW format. The computer then processes the images into the RGB format. This mode produces images with less noise and better color fidelity. Of course, transfer of the RAW data is slower than the normal mode which causes a lower maximal frame-rate.

**Hue** Hue shifts image colors across the rainbow.

**Internal Shutter** This option sets behavior of the camera internal shutter.

**Opened** Internal shutter is opened upon NIS-Elements and stays opened through the whole session. Only with one exception: if the *Readout Mode* is set to *Normal 80kHz at 16bit*, the shutter stays closed and opens only for exposures.

**Closed** Internal shutter is closed upon NIS-Elements startup and gets opened only for exposures.

**Automatic** Shutter is closed when the NIS-Elements is running. It opens only for acquisition.

**Isolated Crop Mode** This option turns on the *Isolated Crop Mode* described within the Andor documentation. In this special mode, the size of chip is reduced and very high frame-rates can be achieved.

**Keep Overlapped** If checked. The Clear Mode and Sensor Mode boxes are set to Automatic and become disabled. This mode produces maximal frame rate and decreases the list of exposure times (in the camera settings window). Only the times of 1 frame and longer become available.

**Limit Max. FPS to** Insert maximum frame rate (frames per second) you would like to use, this will prolong the total duration of the time-sequence in exchange for the reduced frame rate. This option may be useful especially when using the *1 frame for Fast Timelapse* button with a small region of interest. Then the frame rate could get unrealistically high.

---

Note

*The Limit MAX FPS is possible only when camera is in Normal mode (not overlapped), the maximum fps in normal mode is shown next to the edit box.*

**Live Acceleration** If the exposure time is too long, *Live Acceleration* helps you to make it shorter. The system automatically shortens the exposure time (so the frame rate rises), and the loss of intensity is compensated by increasing the gain (software multiplication). This procedure is not used upon *Capture*.

**Maintain Temperature on Camera Shutdown** If this option is checked, the camera maintains its temperature until the computer workstation shut down. The temperature will be maintained for example during NIS-Elements restart or when another camera is selected.

**Max gain** Maximal camera gain.

**Maximum Exposure** This is a safeguard of the time of the Auto Exposure. For quick exposure, it is convenient not to set this value too high.

**Metering Mode** If this option is available, auto exposure can be calculated with an emphasis on overexposed peaks (*Peak*) or average pixel intensity (*Average*).

**Mode** Defines how the Auto Exposure is performed. Average is suitable for bright field and Peak is better used for dark field.

**Multiplier** Controls on-chip multiplication gain. There is complex relationship (approximated to an exponential) between the multiplier value and the level of multiplication - in other words, the value 2 does not mean the signal is multiplied two times.

**Noise Reduction** If activated, it reduces noise generated by higher gain. NR works on interframe basis so you may notice some shadows on moving objects.

**Offset** Sets the brightness of the image. It is a constant additive (positive or negative) changing all pixel values of the image. With negative offset value the dark image areas become pure black. Considering fluorescence microscopy, an appropriate offset setting can create continuous black background and thus help to enhance (together with gain or illumination enhancement) the contrast.

**Overillumination Tolerance** Sets how many pixels should be white after the Auto exposure is performed. Use lower values (0.01%) for very bright (shining) objects (fluorescence). For common bright field, even 1% may be a good value.

**Readout Mode** Within this option you can select a combination of these parameters - readout speed and bit depth. The settings influences frame rate and image quality. Higher frequencies increase frame rate, lower frequencies offer better image quality.

**Readout Rate** Select the readout speed 100 or 280 MHz. If speed is not crucial, leave *100 MHz*.

**Readout Speed** The speed of digitization - influences frame rate and image quality. Higher frequencies increase frame rate, lower frequencies provide better image quality.

## Resolution

### Note

*If the resolution of 4076x3116 is used for capture and the scale is burnt to the image, the scale position is different from the Live image. This is a consequence. To keep the scale at the same position as in the Live image, use other resolutions for capture.*

### Note

*The High Quality Capture mode is also supported. If the Quality is 16bit or the resolution of pixel shift mode is selected, you cannot uncheck the High Quality Capture check box.*

**ROI** Select this option to activate the camera ROI defined via the *ROI Size* drop-down menu. When active, the current ROI size is displayed in brackets. If a binning is applied, this size which is the size of the ROI on the camera chip does not match the size of the resulting image. Also, if you define the ROI with the binning turned on, the pixel size displayed in the definition dialog does not match the value then displayed in the brackets because it includes the binning factor.

The user can choose a ROI size from the list or define a custom ROI via **Define ROI**. The ROI can also be saved/loaded to/from a `.camroi` file (*Save ROI/Load ROI*).

**Saturation** The amount of saturation determines how colors are rendered. More saturation produces richer colors. Less saturation makes the colors gray.

**Save Settings, Remove Settings, Load Settings** Save the current camera settings by clicking *Commands > Save Settings > New Settings...* and give it a name. Load the saved settings by selecting its name in the *Commands > Load Settings* list. To remove a setting, click on it inside the *Commands > Remove Settings* list.

**Scene Mode, Preset** There are several camera settings presets optimized for a specific usage. You can reset the settings using the *Reset* button.

**Neutral** Neutral preset.

---

**Industrial microscopy:**

**W (Water/IC)** Wafer IC-Chip.

**M (Metal)** Metal or ceramic.

**C** Circuit boards.

**FPD** Flat panel displays.

**Bioscience/Biological microscopy:**

**DF/FL, F** Dark-field fluorescence.

**DIC/PH** DIC or phase contrast.

**LED-Bright Field, B** LED illuminated bright field.

**HAL-Bright Field** Halogen illuminated bright field.

**HAL-HE, H** Halogen illuminated Hematoxylin and eosin stain.

**HAL-ELA, E** Halogen illuminated enzyme labeled antibody method.

**Asbestos** Asbestos.

**Sensor Mode** This option sets the method of reading the CCD chip.

**Automatic** This option makes the system to choose the best mode.

**Normal** A standard mode where the exposure phase is followed by the readout phase. Please, see the camera user manual for further details.

**Alternate Normal** Please, see the camera user manual for further details.

**Overlap** This mode enables the camera to achieve higher frame rates by reducing the exposure+readout time (the readout is performed simultaneously with the exposure of the next frame). Please, see the camera user manual for further details.

**Frame Transfer** This mode enables the camera to achieve higher frame rates by reducing the exposure+readout time (the readout is performed simultaneously with the exposure of the next frame). Please, see the camera user manual for further details.

**Separate Channel Settings** Having a color camera, influence of the settings described above on each color channel can be adjusted. In such case, three edit boxes with the default values set to 1.0 are displayed below the feature setting. These values determine the influence of the feature on each color channel.

**Sharpness** Some camera settings provide the sharpness control which affects how sharp edges in the image appear. Too much sharpness leads to over-saturated edges.

**Speed up Triggered Acquisition + Different Exposures** Connect the *Arm* and the *Ext. Trig* connectors of the camera and enable this option. When you run triggered acquisition with the *Use different exposures* option enabled, it will speed up the performance (the camera will not have to wait for software triggers but will be controlled by the *Arm* triggering signal).

**Spurious Noise Filter** This option reduces spurious noise. Try whether it has desirable influence on image quality.

**Target Maximum Intensity** Restricts the maximum of image intensity after the auto exposure was applied. The value represents a percentage of the whole camera dynamic range.

**Trigger Mode** Sets the exposure method.

**Internal** The exposure time and the beginning of each frame acquisition is controlled by the NIS-Elements AR settings.

**Strobe** In this mode, the beginning of each frame exposure is being controlled by external signal.



**Bulb** The exposure time and frames timing are controlled by external signal connected to the camera.

**Triggered by master camera** Having a dual camera system (or a special triggering device connected to the camera), you can use this mode. If the cameras *Interval* setting is set to *No delay* within an experiment or if the Live mode is run, the slave camera will be triggered only once so that both cameras will run simultaneously while single frames will not be synchronized. If the settings are different, the trigger will be engaged once and all the subsequent frames will be controlled by the NIS-Elements AR settings as in the *Internal* mode.

**Use Current ROI** The currently defined region of interest, set by the *Define ROI* command, can be switched on/off by checking this item.

**Using Probe** The probe  enables you to determine a small image area that serves as the data source area for LUTs, histogram etc. It also affects the AWB (auto white balance) and the AE (Auto Exposure) features of the cameras which support it. When the probe functionality is not supported, the AWB and AE algorithms are computed from the whole image.


#### Vertical Pixel Shift

**Shift Speed** Please, see the camera user manual for further details

**Vertical Clock Voltage Amplitude** Please, see the camera user manual for further details


**White Balance** There are Red, Green and Blue Gain properties that control how colors are rendered. It is used to eliminate a color cast from white areas.

### 4.1.10 Controlling Illumination Devices

The substance of an illumination device is the light source. The light emission is produced mainly by a laser, diode or a light bulb. *NIS-Elements* does not distinguish between wavelength modules of these light sources and all are controlled in a similar manner. Basically you select which wavelength(s) you want to use for illumination and set its light intensity. The selected wavelength modules are turned off until you activate the  shutter button usually placed in the bottom left corner of the control panel.

## 4.2 Visual Device Manager

### Key Features

- The full microscope configuration is displayed visually and controlled by the user using a mouse.
- Simplified work with multi-device hardware configurations. Double-clicking on a device in the scheme opens its control pad.
- Multiple hardware configurations (multiple independent microscope setups) can be set and running at the same time.
- Light path concept - optical paths are colored to match the defined wavelengths. The system computes the correct emission wavelength. Each light path can switch between the camera and the ocular. Switching between a single and a dual camera is safe.
- Simplified definition of Optical Configurations – only devices enabled in the current configuration are visible within the OC Panel (*View > Acquisition Controls > OC Panel* ).

---

## **Hierarchy of Configurations**

- *Hardware setup* - sets which microscope is used for the ongoing hardware configuration.
- *Hardware Configuration* - determines which devices are used together, other devices can be turned off.
- *Light Path* - groups devices (from the hardware configuration) used in one observation mode (EPI, DIA, TIRF, confocal, etc.).
- *Optical Configuration* - specifies the settings of each device in a light path. Each optical configuration is associated with one light path.

## **Modes**

**Hardware Configuration** By default, “Acquisition” hardware configuration is created. The user then adds the connected *Cameras* and *Devices* using the **+** buttons. Lightpath combinations are automatically calculated in the *Lightpath Selection* table. The user usually does not need all the devices connected so any camera/device can be turned off creating a new device combination. This new “hardware configuration” can be saved by clicking on the **+** *Create New Hardware Configuration* button and a new tab with a given name is created. The user can switch to a different configuration using the tab and clicking **Activate**, using the *Devices > Select HW Configuration...* command or on *NIS-Elements* startup. Devices which are disabled for the selected configuration are not loaded and can be switched off to make the startup faster.

To display the system overview with each start of *NIS-Elements*, check *Show Startup System Overview* in *Edit > Options > Visual Device Manager*.

**Light Path** The table inside each hardware configuration contains “Light Paths” which represent all available light paths (light tracks between the cameras and illumination devices present in the current configuration). Simply stated a light path tells which devices are used at the same time. Click on the name of a light path to highlight it in the image. Check *Edit* to select which light paths will be shown in the current hardware configuration.


Different light paths are usually saved under different hardware configurations to make the microscope operation most powerful and efficient.

### Note

---

*Hardware buttons of the Nikon Ti2-E microscope can be set to switch to a predefined light path. In the 4.2 *Visual Device Manager* select **Configure...** in the context menu over the microscope, switch to the **Programmable Buttons** tab, choose a button and click on its drop-down arrow, click **Select Lightpath** and set a light path from the list.*

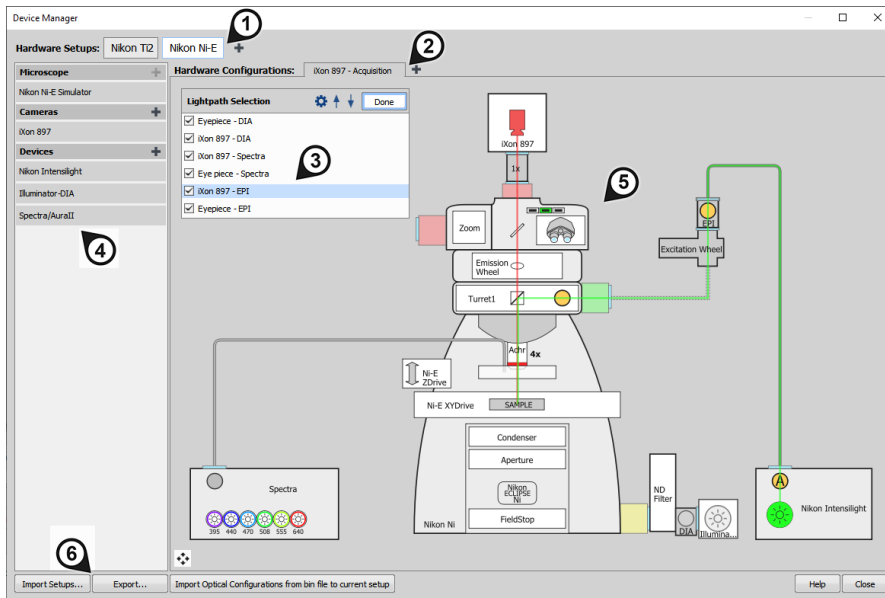Figure 4.10. Device Manager Overview



### **Device Manager window**

1. Hardware Setup tabs - select which microscope is used. More setups are added by the **+** button.
2. Hardware Configuration tabs - select a hardware configuration. More configurations are added by the **+** button.
3. Lightpath Selection table - displays all available light paths for the current hardware configuration and selects the active light path. Use **Edit** to select which light paths will be hidden/shown and change their order using the **↑ ↓** arrows. To add/remove the light path with all devices, click **Settings...** and check/uncheck *All Devices*.
4. Microscope/Cameras/Devices list - shows all connected microscopes, cameras and devices. To add more, use the **+** button.
5. Microscope scheme - shows a visual microscope scheme of the current setup. The system can be configured directly from this scheme using mouse dragging and context menus.
6. Device Manager Settings - **Export...** exports the whole device manager setup (hardware configurations, light paths, optical configurations, etc.) into a .bin file whereas **Import Setups...** can be used to load the previously exported setup. **Help** opens this help page.

### 4.2.1 Connecting Cameras

To correctly setup a camera using the *Visual Device Manager* (see 4.2 Visual Device Manager ), follow these steps:

- 1) Start *NIS-Elements*.

- 
- 2) Select a microscope - click on **Select Microscope** and select your microscope from the list. A new hardware configuration of the microscope is loaded.
  - 3) In the left panel, click on **Cameras +** .
  - 4) Select your camera from the list and click **Add Camera**.
  - 5) Make sure the camera (rectangle with the camera symbol and the camera name) is visually mounted to the proper port to match your physical setup. If not, drag and drop it to the correct place. Available positions are highlighted yellow.

#### Note

*Non-standard camera mounts can be replaced by “Coupler Universal” (found in Devices) which enables the user to connect any camera to any port. The system automatically detects non-standard camera connections. In such cases the system automatically prompts the user to Add Coupler. “Coupler Universal Switcher” enables to connect two devices to a single port.*

- 6) Reveal the context menu over the camera to:
  - disable the camera in the current hardware configuration,
  - remove the camera from the configuration,
  - change the camera to a different one,
  - select the camera mode,
  - rename the camera,
  - make a connection with a different camera,
  - open the camera light path settings.
  - hide the camera,
  - set the camera rules for active lightpath, or
  - open its configuration dialog (*Configure...*).

## 4.2.2 Connecting Devices

Connecting a device is similar to connecting a camera.

- 1) In the left panel, click on **Devices +** .
- 2) Select a device from the list and click **Add Device**.
- 3) Make sure the device is visually mounted to the proper port to match your physical setup. If not, drag and drop it to the correct place. Available positions are highlighted yellow.


## Note

Non-standard device mounts can be replaced by “Coupler Universal” (found in Devices) which enables the user to connect any device to any port. The system automatically detects non-standard device connections. In such cases the system automatically prompts the user to Add Coupler. “Coupler Universal Switcher” enables to connect two devices to a single port.



- 4) Reveal the context menu over the device to:
- disconnect the device,
  - disable the device in the current hardware configuration,
  - remove the device from the configuration,
  - show the connection dialog,
  - rename and change any parameters of the device, or
  - set the device rules for active lightpath, or
  - open its configuration dialog (*Configure...*).

### 4.2.3 Lightpath Scheme

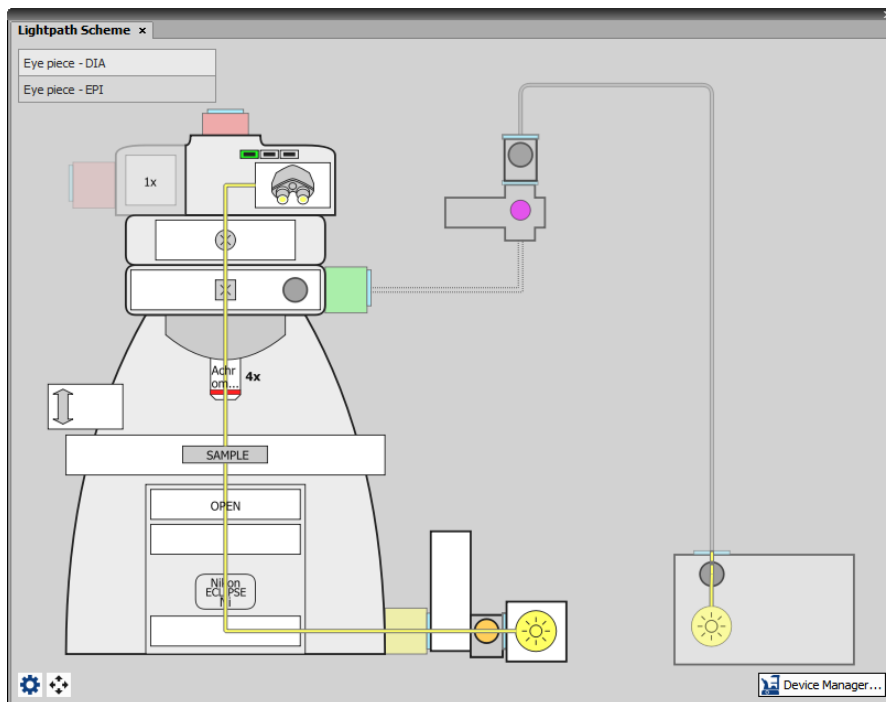Figure 4.11. Lightpath Scheme pad




---

This control pad invoked by the *Devices > Show Lightpath Scheme*  command shows the current microscope configuration with visible light paths which indicate the physical tracks of the excited and emitted light in real colors matching the defined wavelengths (*Real Color - Additive/Spectrum*), or in *Pseudo Colors*. Dashed paths indicate that the light does not pass through. A warning symbol (  ) points out what prevents light from passing through. The warning symbol next to a light path name indicates that device(s) in the particular light path are not set correctly. Such light paths are not saved. Thanks to this indication the user can quickly detect possible errors and setting mistakes (e.g. wrong filter cube assignment) and fix them directly in the pad.

Microscope buttons, shutters, lights, etc. can be controlled directly by clicking on them in the scheme. Double-clicking on a device opens its control pad. Devices not participating in the current light path are outlined in gray and can still be controlled.

#### Note

---

*Stimulation light paths are not shown in the Light Path Scheme. If you need to use the stimulation light path for acquisition and not only for stimulation, enable the acquisition light path under the [Edit](#) button in the 4.2 Visual Device Manager . If this light path is missing, make it visible by clicking the  Settings... button and checking it.*

### **Context menu over a device**

**Activate Lightpath with this device** Finds and activates a light path being the most similar to the currently selected light path and containing the device being clicked at the same time.

**Rules for Active Lightpath** Specifies further rules for the clicked device in the current light path:


*Set State/Position via OC* - controls the device state/position/settings using *Optical Configurations* when the current light path is selected.


*Keep Previous State/Position* - does not change the device state/position/settings when the current light path is selected.

*Active Device* - uses the device as an active device in the current light path and sets all other similar type devices to their default behaviour.

*Position* - always keeps the device in the position selected from the drop-down menu when the current light path is selected.

**Configure...** Opens the configuration dialog of the clicked device.

 **Lightpath Scheme Settings** Enables to change the color of the light paths (*Lightpath Colors*), turn on the table with listed light paths (*Show Lightpath List on Pad*), display labels on each device of the microscope ([Show Labels on Pad](#)) and turn on/off the warning message when light path changes and non-motorized devices positions do not match.

 **Fit to Screen** Adjusts the zoom so that the microscope scheme fits into the *Lightpath Scheme* pad. The scheme can be zoomed in/out using the mouse wheel.

 **Device Manager...** Opens the *Visual Device Manager* (see 4.2 Visual Device Manager ).

---

## 4.3 Legacy Cameras and Devices

Cameras and devices called “legacy” are no longer officially supported but should work with the current version of *NIS-Elements*. Drivers for these accessories are still available within the program setup, but are not tested any more.






## 5. Image Acquisition


### 5.1 Introduction to Image Acquisition


Having *NIS-Elements* installed and all hardware accessories set up, you can start capturing images. Let's begin with the simplest case.

#### How to capture a single image


- 1) Turn the connected camera and other devices ON and start *NIS-Elements*.
- 2) Run *Acquire > Camera Settings*  to display the *Camera Settings* control panel.
- 3) Switch camera to the *Live - Fast* mode (*Acquire > Live - Fast*)
- 4) Adjust resolution of the *Live - Fast* mode to get a continuous Live image. It is recommended to set a low resolution in order to achieve high frame rates. This is good when searching the specimen or when focusing manually.
- 5) Adjust *Exposure time* to get a nice image of the scene.
- 6) Focus on the scene.
- 7) Optionally, turn on the camera ROI. The *Define ROI* button appears on the toolbar if the camera ROI function is supported by your camera. See also 5.2 Camera ROI.
- 8) Capture the image by running the *Acquire > Capture* command
- 9) A new image is opened automatically and named "Captured".

#### Tools for Handling the Live Image

 **Live - Quality, `Ctrl` + `+`** | If a high frame rate is not crucial or if you would like to view the live image exactly as it will be captured, run the *Acquire > Live - Quality* command. It displays the live image in the *Quality (Capture)* resolution. This resolution is used whenever an image is captured.

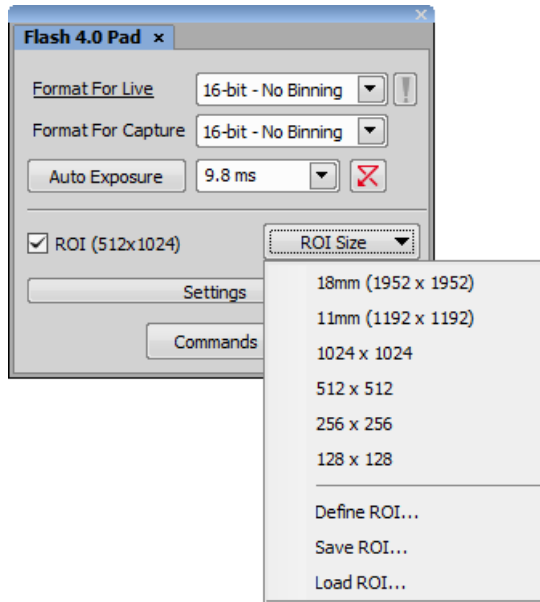
 **Capture, `Ctrl` + `-`** | When pressed, the camera exposure runs till the end of the current frame, and the next frame is captured and displayed on the screen (it is the first frame with the complete exposure after you have pressed the *Capture* button).

If you are using the *Live - Fast* mode and click *Acquire > Capture*, *NIS-Elements* will automatically switch to the quality mode to capture the image. When the image is captured, it is opened on the screen in a new image window.

 **Freeze, -** | The *Freeze* button interrupts the camera exposure, and displays the very last complete frame.

## 5.2 Camera ROI

You can specify a camera ROI on a live image by checking ROI in your camera pad. Please see 4.1.9 Camera Settings for more information about camera settings.



**ROI** Select this option to activate the camera ROI defined via the *ROI Size* drop-down menu. When active, the current ROI size is displayed in brackets. If a binning is applied, this size which is the size of the ROI on the camera chip does not match the size of the resulting image. Also, if you define the ROI with the binning turned on, the pixel size displayed in the definition dialog does not match the value then displayed in the brackets because it includes the binning factor.

The user can choose a ROI size from the list or define a custom ROI via **Define ROI**. The ROI can also be saved/loaded to/from a `.camroi` file (*Save ROI/Load ROI*).

## 5.3 Shading Correction


Shading correction is a method which can correct illumination inhomogeneities of the captured images. How does this work? First of all, a “shading image” must be obtained. It is an image which represents illumination intensities within the field of view.

Figure 5.2. Example of a shading image showing inhomogeneous illumination




As you can see in the image above, the illumination intensity is 100% in the center but gets darker near the edges. This can be corrected by applying the shading correction which will equalize intensities of the resulting image based on the captured “shading image”.

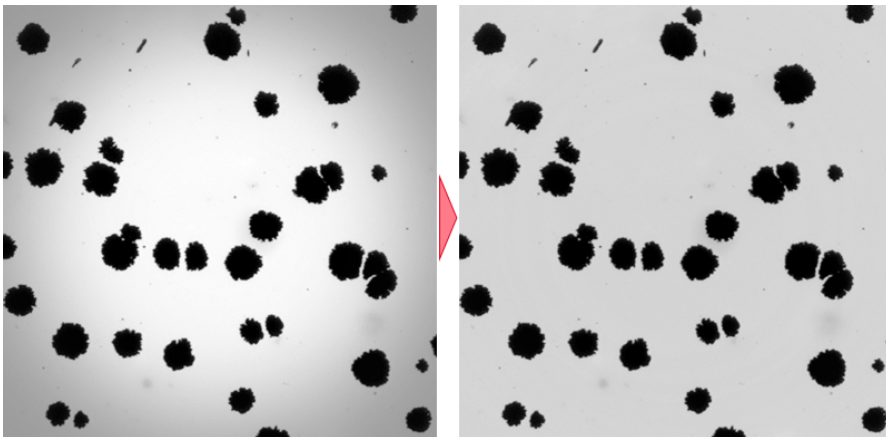Figure 5.3. Image before and after applying the shading correction



### 5.3.1 Simple Shading Correction

- 1) Run the *Acquire > Shading Correction Panel* command to display the panel.
- 2) Select the *Simple* mode in which one correction image is used for all optical configurations.
- 3) Click the  **Capture** button, a dialog appears.
- 4) Find an empty clean part of the glass and move slightly out of focus.

- 5) Select image modality (background brightness) in the dialog.
- 6) Capture multiple images (10 is reasonable) by pressing the **Next >>** button. It is optimal if you move the stage between the acquisitions. Images will be merged together to form the final correction image. The *Background Variability* parameter is taken into account..
- 7) Click the **Finish** button to save the correction image.
- 8) Start using the correction by clicking the  **Apply Shading Correction to All Acquisitions** button.

### 5.3.2 Shading Correction per Optical Configuration

- 1) Run the *Acquire > Shading Correction Panel* command to display the panel.
- 2) Select the *per Optical Configuration* mode in which one correction image is used for each optical configuration.
- 3) Click the **...** button in the *Capture Shading Image* section of the dialog and select the optical configuration for which you want to capture the correction image.
- 4) Click the **Capture** button and proceed in the same way described above (5.3.1 Simple Shading Correction)
- 5) Start using the correction by clicking the  **Apply Shading Correction to All Acquisitions** button.


### 5.3.3 Removing Dust from Camera Chip Using Shading Correction

If you have dust or dirt on the camera chip, the *Shading Correction* panel can be used to compensate it.

- 1) Run the *Acquire > Shading Correction Panel* command to display the panel.
- 2) Find an empty clean part of the glass and defocus the image
- 3) Run the *Acquire > Capture* command to capture the image. Make sure the captured image stays active.
- 4) Select the camera or optical configuration for which the correction image will be used In the *Shading Correction* panel.
- 5) Click the **Use Current Image as Shading Image** button at the bottom.

### 5.3.4 Shading Correction Panel

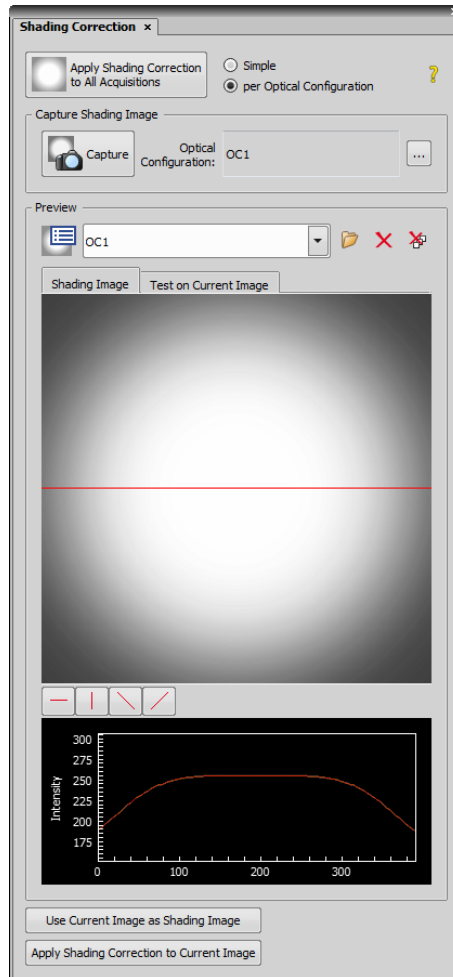Figure 5.4. Shading Correction panel




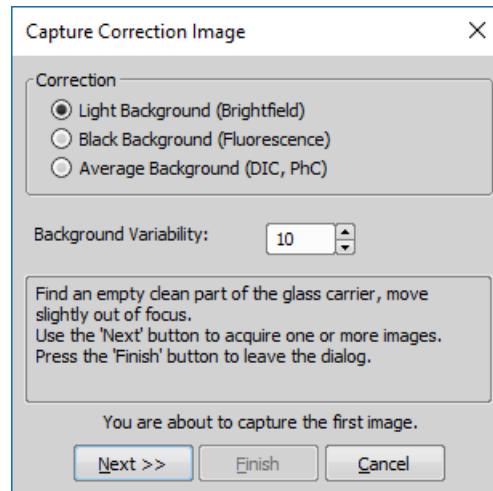
**Apply Shading Correction to All Acquisitions** This button turns the shading correction on/off on the current image (*Simple*) or on the current optical configuration (*per Optical Configuration*).

**Simple** shading correction is applied globally. If multiple cameras are used, select the one used for capturing the shading image in the *Camera* drop-down menu.

**per Optical Configuration** applies the shading correction to the currently selected optical configuration. Use the  button to select an optical configuration and use  to capture the shading image.

**Help** Opens this help page.




 **Capture** Opens the *Capture Shading Image* dialog. The information box gives advice on how to properly capture a shading image. Follow the instruction and click **Next >>** to capture the image. Repeat the capture multiple times (10 times optimally) to capture an accurate shading image, then click **Finish**.




**Correction** Select the type of background to be corrected. It is important for the algorithm to know whether the background is bright, dark, or neutral.

**Background Variability** This value determines the level of detail of the correction algorithm (smoothness of the curve). Homogeneous shading such as in our example above has low variability, so the default value (10) should work fine. Higher values are suitable for more variable shading images. The range is from 5 to 50.

**Preview** The drop-down contains a list of captured correction images. In the *Simple* mode, the correction image has the name of the camera (there can be two items if you have two cameras). In the *per Optical configuration* mode, each correction image carries the name of the optical configuration used during acquisition. Once you select a correction image, its preview appears in the *Shading Image* tab. You can:

- Open it to the main window by the  button.
- Delete it from the list by the  button.
- Examine its intensity profile at the bottom of the panel. Below the preview area there is a set of buttons for changing the rotation of the profile line providing data to the intensity graph below. The profile line can be moved manually using the mouse.
- Test the correction algorithm on the current image by switching to the *Test On Current Image* tab and clicking the  **Test** button. The profile graph changes accordingly.

 This button deletes all shading correction images for both modes.

**Use Current Image as Shading Image** If the user has a previously saved shading image, it can be loaded into this dialog and used as a shading image.

**Apply Shading Correction to Current Image** Applies the currently selected shading image to the currently opened image.

## 5.4 About ND Acquisition

### 5.4.1 Introduction to ND Acquisition

*NIS-Elements* being a multi-purpose imaging system can be used as a handy tool to study objects, live organisms, processes, etc. The universal ND2 (N-dimensional) file format is what makes this possible. One *ND2* file can contain multiple images organized according to what type of acquisition they came from. There are the following acquisition types:

#### Note

*Each acquisition type gives a name to the corresponding dimension of the resulting ND2 file. For example, we can speak about two-dimensional TZ ND2 file, which means the file contains the Time-lapse dimension and the Z-series dimension.*

**Time-lapse - T** A sequence of images can be captured over a period of time and create a time-lapse image. See 5.5 Time-lapse Acquisition.

**Multi-point - XY** Images from different coordinates of the XY stage can be acquired. See 5.6 Multi-point Acquisition.

**Z-series - Z** Several techniques utilize Z-stacks of images. Such Z-stack can be for example converted to a 3D model of the specimen. See 5.7 Z-series Acquisition.

- Single-dimensional documents can be created automatically (using motorized accessories) or manually using commands from the *Acquire* menu.

### 5.4.2 Common ND Experiment Options

The following options can be applied to all types of ND experiments:

**Experiment** You can enter a custom name of the experiment. Also name of the phase is editable.

**Path** Browse to a folder where your ND2 images will be stored.

**Filename** Specifies the name of the file to be saved.

**Custom Metadata** Check this option to add custom metadata to the image(s) being captured. Select the metadata set from the drop-down menu and the metadata items present in the set are automatically listed below.

For information about creating and editing custom metadata please see 5.4.2.1 Custom Metadata.

**Autofocus** Automatic focusing can be used during the experiment. You can select the auto-focus method that best meets your needs. The *Define* button shows a window where you can define parameters of the selected focusing method.

**Closing Shutter** The active shutter can be closed between single image acquisitions. Just select the *Close active Shutter when idle / Close active shutter during filter change* option in the experiment window.

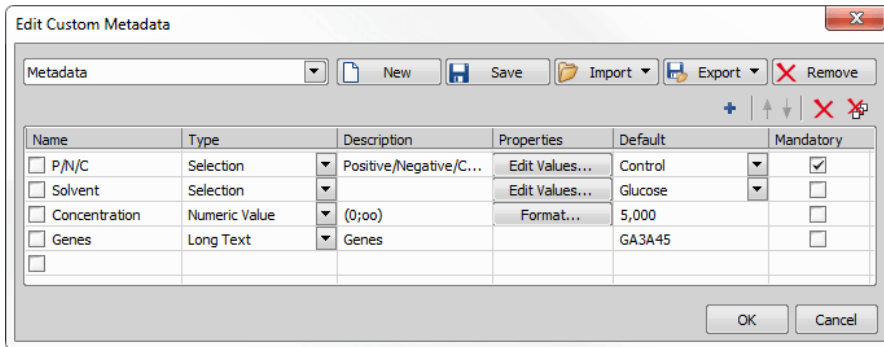
**Running Macro Commands** You can define a command (or a macro) to be run in various stages of the experiment. Select the timing in the *Advanced* section of the experiment window and enter the command to be run.

**Load** Loads a configuration XML file. The file with the current configuration can be created using the *Save* button.

**Save** Saves the current configuration of this dialog to an XML file. A standard *Save As* window appears.

### 5.4.2.1 Custom Metadata

To create or edit existing metadata, check the *Custom Metadata* checkbox in your ND experiment and click **Edit**. The following dialog window appears.



To create a new custom metadata set - click on **New** and give it a name. Click **+** Add New and *Name* your first metadata item, choose its *Type*, enter a text *Description*, edit its *Properties*, choose the *Default* value and select whether it is *Mandatory* or not. Add as many metadata as you like and confirm their definition by clicking **OK**.

To edit an existing metadata set, simply change parameters of the items in the table, add more items ( **+** ), change their order ( **↑ ↓** ), or remove one ( **×** ) or all ( **✖** ).

You can save the whole metadata set as a .xml file by clicking **Export**. To import your saved metadata set, use **Import**. To delete the current metadata set, click **Remove**.

Note

The created metadata can be changed after acquisition in the image properties dialog (File > Image Properties).

### 5.4.3 ND Sequence Options

The following options can be applied to time, multipoint and multichannel sequences in ND experiments:

**+** **Incubation** Adds a new incubation phase into the sequence (Time-lapse acquisition only, (requires: Stage Incubator)). For more information please see 5.5.1 Special Options.

**+** **Add** Adds a new step into the sequence.

**☑** **Select All** Selects all steps in the sequence.

**☐** **Clear Selection** Clears the step selection.


**↑** **Move Up**, **↓** **Move Down** These buttons move the current step up/down in the sequence.

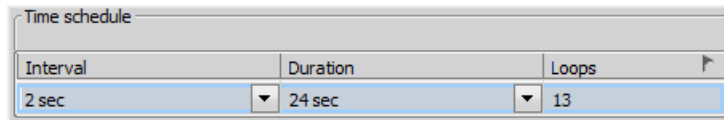
**×** **Remove Current** This button removes the selected step from the sequence.

**✖** **Remove All** This button removes all steps from the sequence.



## 5.5 Time-lapse Acquisition

Detailed studying of long-lasting processes is enabled by the time lapse acquisition mode of *NIS-Elements*. The achievable experiment duration is limited by the hardware abilities of your PC only. Invoke the *Acquire > Capture Timelapse*  > *Capture Automatically* command to set the experiment up:



The *Time schedule* table enables you to define one time phase where duration, interval between single images, and number of images of the phase can be adjusted. The *Interval*, *Duration*, and *Loops* settings are bound together, so you just need to set two of these parameters. The remaining parameter is calculated automatically.

The flag mark indicates the Duration/Loops priority. For example: if the camera exposure time exceeds the defined *Interval* between frames, than the experiment settings are not achievable - only the marked column will be pushed to be correct. You can set the priority column by clicking on its caption.

### 5.5.1 Special Options

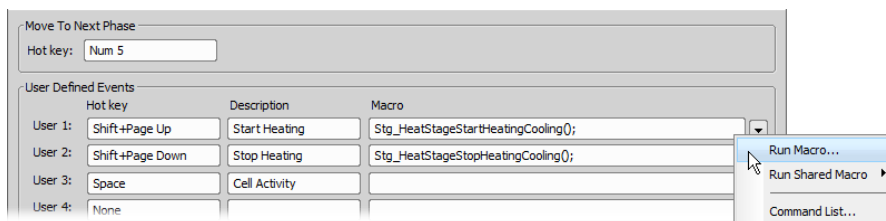
The following options can be used with this type of ND acquisition. Options common to all types are described here.

**Switch Transmitted Illuminator off When Idle** If this option is selected, the lamp is turned off whenever there is a long-enough gap between two acquisitions.

#### Warning

*It takes a little time for some lamps to reach the set intensity. In such case, we recommend to use the `Wait()` ; macro function before the capture is performed.*

**Events** Press the **Events...** button to display the following window:



Define user hot keys which run user events during time-lapse acquisition. Simply click inside the *Hot key* field of the User events window and press any key combination you want to assign to the event. Then you can enter the event description. A macro command can be specified which will be executed when the event occurs (is triggered by the user). When the event occurs, the specified command is run and a marker is placed to the time line of the ND2 file.

### 5.5.1.1 Advanced options

These options are shown after clicking **Advanced >>**. Following features are available.

**Auto Focus** Automatic focusing can be used during the experiment. You can select the auto-focus method that best meets your needs. The *Define* button shows a window where you can define parameters of the selected focusing method.

#### Note

*It can be determined whether the focusing shall be performed at the beginning of the whole experiment or at the beginning of each phase. If the auto-focus is set to be performed Before each Phase and the ND Experiment contains just a single time phase, the automatic focus will not be performed. To run auto focus, use the At the Beginning option.*

**Execute before/after each Time** This field enables to run a command which is executed before or after each loop. Click on the right drop-down button to enter a macro from a file (*Run macro...*), a macro command from the list (*Command List...*) or a macro command from the list performed on every N-th loop (*Command List Every N...*).

#### Note

*Image processing macro commands are not executed during the acquisition.*


**Execute before/after Time phase** This field enables to run a command which is executed before or after each time phase. Click on the right drop-down button to enter a macro from a file (*Run macro...*) or a macro command from the list (*Command List...*). *Advanced* for combo box specifies to which phases the macro commands are assigned. To assign them to all phases present in the current settings, click **Apply to All**.

#### Note

*Image processing macro commands are not executed during the acquisition.*

## 5.6 Multi-point Acquisition

(requires: Stage)

This window serves for defining XY(Z) points to be scanned during the multipoint capture experiment. This feature is available when a motorized stage XY(Z) is present in the system. Check the *Include Z* box to display the Z column. The list of defined points can be saved (and loaded later) to an XML file by the *Save (Load)* button. Run the *Acquire > Capture Multipoint*  > *Capture Automatically* command.

### 5.6.1 Point by Point (Manual) Multi-Point

1. Move the stage to the first position.
2. Press the **+** *Add New* button. A new line containing current coordinates of the stage appears in the list.
3. Move to the next position and repeat the steps until you have all the intended points defined.

## 5.6.2 Well Plate (Rectangular) Multi-Point

You can insert a pattern designed to cover a wellplate. Of course it can be used to create any rectangular pattern.

- 1) Click the  button. The following window appears
- 2) Select *Manual* if you know the distances between wells. Otherwise select the *Interactive* and continue.

**Manual** This method will create a pattern which will scan *Rows* x *Columns* fields. The *Distance X* and *Y* parameters specify distances between two fields. The scanning will begin at the current position.

**Interactive** The interactive method lets you specify the number of fields to be scanned and starts a wizard where you will specify the top-left and the bottom-right corner of the well plate. The distances between wells are calculated automatically.

- 3) Click .

## 5.6.3 Large Image (Covering) Multi-Point

A Multi-point which will cover certain area can be created:

- 1) Click the  button and select the *Large Image* tab
- 2) Set the large image settings:

**Scan Area** Select whether to define the *Scan Area* by number of fields or the actual field size.

**Camera** Select the camera to be used. This field is enabled only if you have a dual-camera setup.

**Objective** Select the objective for which the multi-point will be calculated. The resulting number of fields / area size changes depending on the selected objective magnification.

**Overlap** Overlap of neighboring images in %.

- 3) Click .

## 5.6.4 Random Multi-Point

Specify the area where a random set of points will be generated.

- 1) Click the  button and select the *Random* tab.
- 2) Select shape of the area to be filled with random points

**Rectangle** The area will be defined by a rectangle. Define X and Y coordinates of left top corner of the area, width and height of the area.

---

**Radius** The area will be defined by a circle. Define X and Y coordinates of center of the circle.

- 3) Specify *Number of points* in the multi-point set.
- 4) Click .

## 5.6.5 To Change a Single Z Coordinate

1. Make sure the *Move stage to selected point* is pressed. This button ensures that the motorized stage moves to the coordinates of the currently selected point.
2. Click inside the line you would like to change.
3. Move the Z drive to the new position.
4. Click the <- button.

### Note

---

*The XY coordinates of one point cannot be adjusted (unless you delete it and add a new point).*

## 5.6.6 Special Options

**Autofocus** None, Steps in Range, Steps in Continuous.

**Custom** Displays a tool for creating predefined multi-point patterns.

See 5.6.2 Well Plate (Rectangular) Multi-Point, 5.6.3 Large Image (Covering) Multi-Point, 5.6.4 Random Multi-Point.

**Define Stimulation Laser Powers** If checked, *Laser Power Settings* dialog window appears when defining stimulation ROIs. This window enables to define power of each laser line on each stimulation group.

**Execute Command after Capture** You can either Run Macro, or select a command from the Command list. The command is executed after capturing.

**Execute Command before Capture** You can either Run Macro, or select a command from the Command list. The command is executed before capturing.

**Include Z** If selected, also the Z coordinate will be taken into account during XY(Z) acquisition.

### Note

---

*This option is disabled if the Relative XY option is used.*

**Leave PFS offset ON between points** Check this item to keep PFS on while moving between the points.

**Move Stage to Selected Point** This button ensures that the motorized stage moves to the coordinates of the currently selected point.

### Offset All

The  button can shift the XY coordinates of all points in the same way:

1. Make sure the *Move stage to selected point* is pressed. This button ensures that the motorized stage moves to the coordinates of the currently selected point.

2. Select one point of the list.
3. Move the XY(Z) stage to a new position (define the offset).
4. Press the *Offset All* button.
5. The coordinates of all points change - the same shift which you made with the stage is applied to them.

**Optimize** If the *Optimize* button is pressed, the system will re-order the defined points in order to minimize the XY stage trajectory.


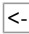

**Point name** Displays the name of the point. You can change the name (default names are #1, #2 #3, etc.).

**Redefine reference Z after Auto Focus** Check this option to redefine reference Z position after performing autofocus or PFS (if AF or PFS is turned on).

**Relative XY** Check the *Relative XY* item to consider all coordinates as relative with respect to the current stage position. Any of the points may be used as the reference point, just right-click it and select *Set this point as a reference position*.


**Stim. Device** Defines which connected device will be used for stimulation.

**Use Focus Surface** Uses Focus Surface for capturing.

**X, Y, Z** Displays X, Y and Z coordinates of the point. The arrow buttons   assign the current XYZ position to the point. The  button shifts the XYZ coordinates of all points by the same offset which is defined as a difference between the current stage coordinates and coordinates of the selected point.



## 5.7 Z-series Acquisition

(requires: Z Drive)

Automatic capturing of images from different focal planes using a motorized Z drive can be performed. Run the *Acquire > Capture Z-Series*  *> Capture Automatically* command to display the setup window. There are three different approaches of how to set the experiment:

### Defined by top bottom



*Required values: top position, bottom position, step size or number of steps.*

- 1) Click on the  *Defined by top bottom* button, the cube turns blue.
- 2) Run the  Live camera signal.

Note

*If using two Z devices, choose one Z device from the pull-down menu.*

- 3) Define the Z-range of interest:

- Move the Z drive to the top position and press  *Top*.
- Move the Z drive to the bottom position and press  *Bottom*.


---

Note

Both values can be adjusted in the Bottom and Top edit boxes below.

- 4) Enter the Step size in  $\mu\text{m}$  (distance between adjacent steps) or specify the number of Steps to be captured.



Note

This mode assigns the Home position automatically to the middle step (e.g. the third step out of the whole 5 steps).  **Reset** discards the Top, Home, and Bottom position settings.

Depending on the direction of Z-acquisition, the last Z position (top or bottom) may slightly differ from the user settings. However, you can select which position will be preserved exactly. Right-click on one of the buttons (Top or Bottom) and select Keep exact Top Position / Keep exact Bottom Position. This setting is indicated by underlining/overlining the button text.




### **Symmetric mode defined by range**

Required value: range.

- 1) Click on the  Symmetric mode defined by range button.
- 2) Run the  Live camera signal.



Note

If using two Z devices, choose one Z device from the pull-down menu.

- 3) Move the Z drive to the absolute home position (center of your Z series) and click  Home. If you do not know the exact center position yet, you can then switch to  Relative and the home position will be taken from the current Z position prior to the actual acquisition.
- 4) Enter the Range to be captured and then adjust the Step size in  $\mu\text{m}$  (distance between adjacent steps, for the recommended value click  Use suggested step size) or enter the number of Steps to be captured within this range.



### **Asymmetric mode defined by range**

Required values: above position, below position.

- 1) Click on the  Asymmetric mode defined by range button.
- 2) Run the  Live camera signal.

Note


If using two Z devices, choose one Z device from the pull-down menu.


- 3) Move the Z drive to the absolute home position (center of your Z series) and click  Home. If you do not know the exact center position yet, you can then switch to  Relative and the home position will be taken from the current Z position prior to the actual acquisition.

- 4) Specify the range from the home position by entering the *Below* and *Above* values and adjust the *Step* size in  $\mu\text{m}$  (distance between adjacent steps, for the recommended value click  $\leftarrow$  *Use suggested step size*) or enter the number of *Steps* to be captured within this range.

### **Special Options**

**Piezo Z** If a Piezo Z drive device is connected, the *Piezo* button appears in the panel. Use the pull-down menu to select the action which will be performed upon pressing the button:

 **Keeps Z position and centers Piezo Z** This option moves the Piezo Z drive to the home position, but compensates the movement by the second Z drive in order to keep the original absolute Z position (sum of Z1 and Z2).

 **Move Piezo Z to Home position** Moves the Piezo Z drive to the home position regardless of the absolute Z drive position.

**Direction** You can define the Z-stack scanning direction - either from *Top to Bottom*, or vice versa (*Bottom to Top*).

**Two Z devices** Two Z devices (Main Z + Piezo Z) can be used for Z-Series acquisition. Z device combination is used automatically if AF/Z-Stack range exceeds the range of Piezo Z. Combining two Z devices speeds up the Z stack acquisition, however the user can still decide whether or not to use the Z device combination. Only the Main Z device is used if it is selected in the *Z device* combo box. Only the Piezo Z device is used if it is selected in the *Z device* combo box and the defined *Range* is within the device range. Combination of both Main Z and Piezo Z is used if Piezo Z is selected in the *Z device* combo box and the defined *Range* is bigger than the device range.







## 6. Displaying Images


### 6.1 Open, Save, Close Images

#### 6.1.1 Opening Image Files

##### 6.1.1.1 Different Ways to Open an Image File



*NIS-Elements* offers several ways to open an image file, using either:

**File > Open**  **command** To invoke the Open dialog box where you can select the file to be opened, run the *File > Open*  command. This command is also called when you click the *Open* button on the main toolbar.

**Organizer** An image can be opened by double clicking its filename within *Organizer*. Run the *View > Organizer Layout*  command **F10** to switch to the Organizer. See 6.6 Organizer.

**Recent Files List** You can quickly access the last opened images using the *File > Recent Files* menu.

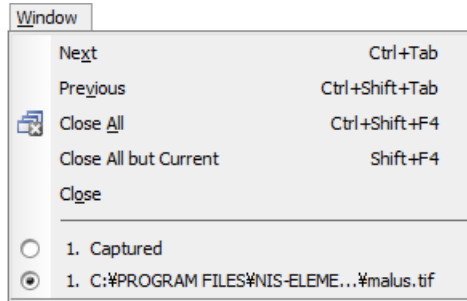
**Open next/previous/first/last Commands** These commands enable you to continuously open the subsequent images from a particular directory or a database table. The *File > Open/Save Next > Open Previous*, *File > Open/Save Next > Open Next*, *File > Open/Save Next > Open First*, *File > Open/Save Next > Open Last* commands may be used.

**Auto Capture Folder** The *Auto Capture Folder* is a control panel that can be displayed by calling the *View > Acquisition Controls > Auto Capture Folder*  command. It displays images within a selected folder. To change the folder click the  button in the top left corner of the control panel and browse for another folder. Any image of this folder can be opened by a double-click.

**Any File Manager** During installation, *NIS-Elements* creates file associations to files that are considered its native format for storing images (JP2, ND2). The JP2 and ND2 files can be then opened to *NIS-Elements* by double clicking their names within any file manager.

##### 6.1.1.2 Switching Between Loaded Images

Commands for managing the opened images are grouped in the *Window* menu. The recently opened files are listed in its bottom part. The currently displayed image is indicated by the selected radio button. To change the current image, select it from the list or use the *Next* or the *Previous* commands (represented by **Ctrl + Tab** and **Ctrl + Shift + Tab** shortcuts).




### 6.1.1.3 Automatic Image Sequence Recognition

When you select to open an image that is a part of image sequence such as 001.JP2, 002.JP2..., it is recognized automatically and you are offered to convert it to an ND2 file.


## 6.1.2 Saving Image Files

### 6.1.2.1 Basic Workflows

*NIS-Elements* offers several ways to save an image file, using either:

**The File > Save  command** Saves changes in the current image. If the current image is new (was never saved before), the *File > Save As* command is called instead.

**The File > Save As command** Saves the current image and lets the user specify the destination folder, format and file name.

**The File > Open/Save Next > Save Next  command** This command saves the current image (live or static) automatically according to the settings defined within the general options window.


**During image-capturing experiments** Images and image sequences can be saved automatically during experiments such as ND Acquisition (5.4 About ND Acquisition).

### 6.1.2.2 Saving Images with UAC

Because of the security enhancement with the UAC (user account control) of *Windows Vista* and *Windows 7*, it is not possible to save images in those folders that require Windows administrator user rights. Those folders are:

- C:\Windows and subsequent folders
- C:\Users and subsequent folders excluding C:\Users\[login-user-name]
- C:\Program Files and subsequent folders excluding C:\Program Files\NIS-Elements\Images

### 6.1.3 Closing Images

- The currently displayed image can be quickly closed by pressing the cross button in the top-right corner of the image window.
- The image can be also closed by invoking the *File > Close* command.
- If you want to close all images, use the *Window > Close All*  command.
- If you try to close an image that has been changed, *NIS-Elements* will display a confirmation dialog box, offering to save the changes.
- Use the *Window > Close All but Current* command to close all opened documents but keep the current one opened.

### 6.1.4 Supported Image Formats

*NIS-Elements* supports a number of file formats for opening. In addition, *NIS-Elements* uses its own image file format (ND2) to fulfil specific application requirements. The following list contains file formats which can be displayed in the software, however, if you do some changes to the files which you want to save, we recommend to use the native *NIS-Elements* format ND2 for that purpose - use the *File > Save As* command.

**AVI, MP4** *NIS-Elements* imports this movie format as a time-lapse sequence of images. Sound is not supported.

**JP2 - JPEG2000 Format** An advanced format with optional compression rates. Image calibration, text descriptions, and other meta-data can be saved together with the image in this format.

**ND2 - N-Dimensional Image** This is the special format for storing sequences of images acquired during ND experiments. It contains various information about the hardware settings and the experiment conditions and settings. It also maintains all image layers of course.

**NEF, NRW** These RAW images produced by Nikon photo cameras can be displayed in *NIS-Elements* keeping the original bit depth (10 or 12 bits per channel).

**JPG, JFF, JTF - Joint Photo Expert Group Format** Standard JPEG files (JPEG File Interchange Format, Progressive JPEG, JPEG Tagged Interchange Format) used in many image processing applications.

**TIFF, TIF - Tagged Image File Format** This format can save the same amount of meta-data as JPEG2000. TIFF files are larger than JPEG2000 files but are loaded faster. TIFF files have several ways to store image data, therefore there are many versions of TIFF. *NIS-Elements* supports the most common TIFF modalities. TIFF image format in *NIS-Elements* supports also floating point images.

**OME.TIF, OME.TIFF** OME metadata (OME-XML) are included for increased compatibility between applications. This format can handle multi-dimensional images (Z-stack, Timelapse...). Multipoint (XY) images are split and saved as separate files automatically.

**GIF - CompuServe Graphic Interchange Format** This is a file format commonly used on the Internet. It uses a lossless compression and stores images in 8-bit color scheme. GIF supports single-color transparency and animation. GIF does not support layers or alpha channels.

**PNG - Portable Network Graphics Format** This is a replacement for the GIF format. It is a full-featured (non-LZW) compressed format intended for a widespread use without any legal restraints. *NIS-Elements* does not support the interlaced version of this format.

**PSD - Adobe Photoshop** *NIS-Elements* can display Photoshop images. Specialized features of the format such as vector or textual layers will be rasterized.

**BMP - Windows Bitmap** This is the standard Windows file format. This format does not include additional image description information such as author, sample, subject or calibration.

**LIM - LIM Format** Developed for the needs of laboratory image analysis software. Nowadays, all its features (and more) are provided by the JPEG 2000 format.

**LSM - Zeiss LSM Image** *NIS-Elements* can display LSM images. Specialized features of the format may be omitted.

**ICS/IDS image sequence** ICS/IDS image sequences are generated by some microscopes and consist of two files: the ICS file with information about the sequence; the IDS file containing the image data. The ICS file must be stored in the same directory together with the IDS file.

---

Caution

*ND Images containing multi-point XY dimension cannot be saved to a file in the ICS/IDS format. See also 5.4 About ND Acquisition.*

**NDPI, VMS - NanoZoomer files** We can open these files produced by Hamamatsu devices and save the image information to an ND2 file.

**ShuttlePix files** Nikon ShuttlePix Digital Microscopes produce standard TIFF and JPEG files and save the calibration info to a separate file. *NIS-Elements* recognizes the calibration file and loads it along with the image data.

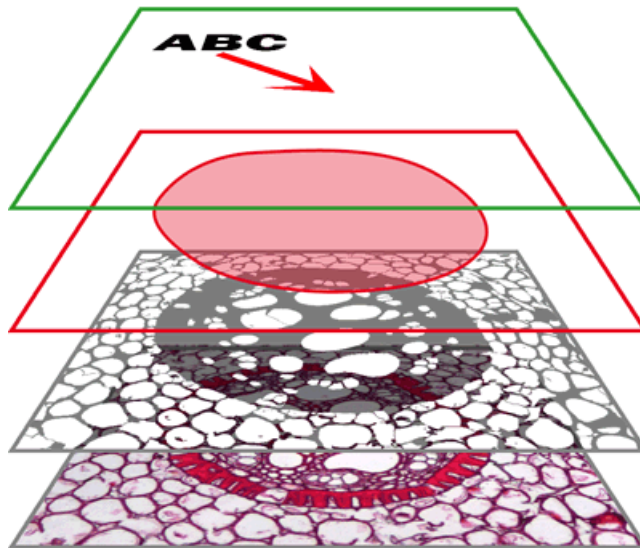
## 6.1.5 Floating Point Images

In *NIS-Elements*, a “floating point image” (FPI) is any image which has its pixel intensities represented by “floating point numbers” instead of integers. The format used is called “single-precision floating-point format” and it extends the available intensity range to 32 bits per channel (intensity can range from  $-3.402823 \times 10^{38}$  to  $3.402823 \times 10^{38}$ ). In short, floating point images allow to store high-precision quantitative information and can be saved in TIFF or ND2 file formats.

## 6.2 Image Layers

### 6.2.1 Introduction to Image Layers

The following image layers can be saved within a *NIS-Elements* image.



**Annotation layer** In this layer, vector objects are stored. The results of manual and automatic measurements, text labels and other annotations can be included.

**Binary layer(s)** The binary layer is usually the result of segmentation. By segmentation, you distinguish objects of interest from background. A picture of night sky can make a nice example - shining stars may be “thresholded” and compose a binary layer. Then, by means of automatic measurement, you can count them. More than one binary layer can be placed over one image. See 7.3 Binary Layers. (*requires: MEAS (Automatic Measurement)*)

**ROI layer** ROIs (regions of interest) is a strong tool for distinguishing objects from background similar to the binary layer. The advantage over the binary layer is that ROIs are vector objects and therefore provide different way of manipulation. See 7.4.1 Introduction to ROIs.

**Color layer** The color layer contains the image data captured by a camera. It can handle images with the depth of up to 16 bits per color component. The dimensions of this layer determines the dimensions of the other layers.

#### Note

*When saving an image, only some file formats are capable of saving all the layers. The other image formats will save the content of the color layer only. See 6.1.4 Supported Image Formats.*

## 6.2.2 Image Types

### RGB Images

Images acquired by a color camera typically consist of three components that represent red, green and blue channel intensities. You can display a single color channel using the tabs located in the bottom-left corner of the image window. Or, an arbitrary combination of them can be selected while holding the `Ctrl` key down.




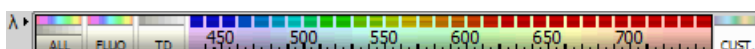
## Multi-channel Images

These documents usually arise from fluorescence microscopy. Instead of 3 color components (RGB), multichannel images can be composed of arbitrary number of user-definable color channels.

### Note

*If an image contains more than 8 components, the tabs in the bottom left corner of the image are replaced by the wavelength dimension, similarly to how other dimension loops of ND2 files are displayed.*

Click the  *Treat As Spectral* button in the top image toolbar to use the current image as a spectral image revealing more channel switching features in the bottom control bar.



**ALL** Selects the full spectrum of the image including spectral (FLUO) and transmitted detector (TD) channels.


**FLUO** Selects only spectral channels. This button is useful for quick switching between FLUO and TD.

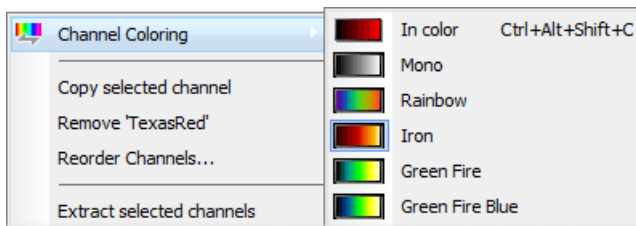
**TD** Displays only the transmitted detector (TD) image.

**CUST.** This button displays custom channels. To set a channel (multiple channels) as custom, make a channel selection in the control bar, right-click over the selection and choose *Set selected channels as Custom*.

Please see 6.3.1 Navigation in ND2 Files, 5.4 About ND Acquisition.

## 6.2.3 Channel Coloring

Single channels and monochromatic images can be visualized using a predefined color scale. Right-click the “channel tab” in the bottom left corner of the image window and select a color scale from the  *Channel Coloring...* sub-menu.



### Brightfield and Fluorescence channel image

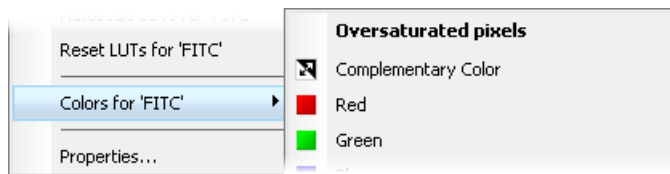
If your ND2 image contains only one brightfield channel with its *Modality* set to *Brightfield* in *Channel Properties*, the behaviour of viewing colors of these channels slightly changes. The brightfield channel is set to *Mono* color by default but can be switched to any other pseudocolor. The fluorescence channels are displayed in their native fluorescence color by default and can be switched to any other pseudocolor. Once the brightfield and fluorescence channels are displayed together in an overlay, the pseudocolor of the fluorescence channel is ignored and is shown in its true color.

## 6.2.4 Assigning Colors to Channels

Despite the color in which a channel is displayed, other per-channel colors can be assigned to it. This regards colors for:

- indication of oversaturated pixels
- indication of undersaturated pixels
- indication of binary pixels

In the context menu over a channel, select *Colors for* and choose a color for oversaturated pixels, undersaturated pixels (these colors are used globally for all images in *NIS-Elements* until next change) and for binary pixels. Selected color for binary pixels is linked to the channel name and is shown in all images having the same channel name.




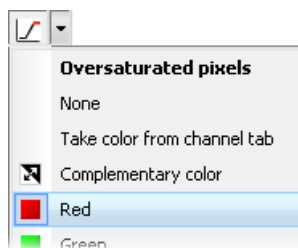
Color of oversaturated and undersaturated pixels may also be selected from the pull down menu next to the *Pixel Saturation Indication* (Ctrl+Shift+S)  button shown in the image toolbar.

Figure 6.7. Pixel Saturation Indication sub-menu



**None** The pixel saturation will not be indicated.

**Take color from channel tab** Color which has been previously assigned to oversaturated and undersaturated pixels in the menu over the channel tab will be used. See 6.2.4 Assigning Colors to Channels.

**Complementary color** Colors complementary to the display color of each channel will be used in order to ensure good visibility of the highlighted pixels.

## 6.2.5 Copying Channels by Drag and Drop

Drag one of the channel tabs and drop it to the left portion of the image window in order to create a new image.

Note


- Even the All and RGB tab can be copied.
- When extracting a channel from Live image, it does not freeze the camera signal.

## 6.3 ND Dimensions

### 6.3.1 Navigation in ND2 Files

#### 6.3.1.1 Control Bar

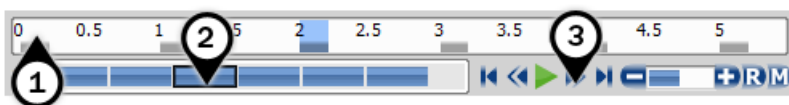Figure 6.8. A Time-lapse image.



When an ND2 file is opened, its structure is pictured at the bottom of the image window.

1. There is a time line with all captured images indicated by gray markers. The blue-highlighted marker indicates the currently observed image frame. You can browse the ND2 file by clicking inside the time line.
2. Below the time line, loops of each dimension are indicated by blue rectangles. Loops can be selected by mouse while holding `Ctrl` or `Shift` keys and are highlighted green.

Note

*ND2 files may contain a very large number of frames. It may be useful to omit some of the frames from the selection. See Selection Context Menu.*

3. Each dimension can be browsed frame-by-frame or you can use the playback controls on the right side.

In specific cases the ND2 file control bar looks differently.

If the number of channels exceeds 10, the image is considered to be a spectral image and a *lambda* dimension is displayed instead of the standard channel tabs.

Figure 6.9. T/Z/multi-channel image containing 11 channels and the time line hidden












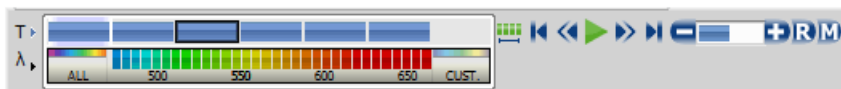



Figure 6.10. HDR multi-exposure image



Figure 6.11. Multi-channel image containing 11 channels



## Playback Controls


-  **Select All Frames** Selects all frames of the dimension and highlights them green.
-  **Play Sequence** Plays all images of the dimension at a defined speed. If a selection is applied, only the selected images are included in the playback.
-  **Stop Playing** Stops playing the sequence at the last displayed frame.
-  **Previous Position**,  **Next Position** Displays the previous/next frame of the dimension.
-  **Decrease Playing Speed**,  **Increase Playing Speed** Changes the playback speed by one step down/up.
-  **Real Time Playing Speed On** Sets the playing speed to real-time (as the image was captured).
-  **Maximum Playing Speed for Every Frame On** Sets the playing speed to maximum while the display of every frame is guaranteed (when the speed is set to maximum by the + button, some frames are usually omitted when playing the sequence depending on your graphics card).
-  **Home Position** Displays the frame of the Z dimension that was set as “home” during the acquisition.

## Selection Context Menu

Right-click a loop (blue rectangle) to display the selection context menu. There are the following commands:

Note

*The placeholder X is the name of the current (clicked) dimension, i.e.: T, Z, M, ...*

- Select All Frames in X** Selects all frames within the current dimension. The *Select Every* option is reset to *Every* automatically.
- Select Frames in X** Opens a dialog where the user can define a range of frames to be selected.
- Invert Selection** Inverts the current selection while preserving the *Select Every* option.
- Select Every** Every Nth frame of the selection is used. First make a selection of frames either by mouse (while holding down `Shift` or `Ctrl`) or select all frames using the *Select All Frames* command or the  button. Then choose the *Select Every* option. Selection is changed to every Nth frame of your previous frame selection.

Note

*Once this option is turned on, it is applied all over again even after making a different frame selection. To turn this function off, switch it to *Every* or *select all frames*.*

**Keep Selected Frames in X, Delete Selected Frames in X** When one of these commands is called, a new image file is created containing only the selected/deselected frames.

**Select Frame (N)** Selects the clicked frame and de-selects all other frames of the current dimension.

**Delete Frame (N)** Deletes the clicked frame of the current dimension and also the custom selection of frames/loops of the other dimensions.

Note




*If you want to keep all frames/loops of other dimensions, make sure you applied the Select All Frames command to them.*

**Tips**



- Detailed info about dimensions will be displayed after you click the leftmost button of the control bar (T>, Z>, ...).
- Place the cursor over one of the dimensions. A tool-tip which displays statistics of the dimension appears.

**Playing options**

Right-click the speed bar to display the following context menu and select the playing mode.

 **Back and Forth**,  **Forward**,  **Backward** Define the direction how the image sequence will be played.

 **Repeat** Sets the repetitive infinite playback.

 **Stop on Events** When you select this option and then press the  *Play Sequence* button, a dialog window appears once the first user event is reached. The dialog window displays information about the reached user event. Use the *Continue* button to continue playing to the next event, or press the *Stop* button to stop the playback at the current frame. Within the window, you can also select the *Do not ask again in this session (Always stop)* option. Playing will then stop automatically on every user event but the window will not be shown.


**Maximum Speed** This command corresponds to the  **Maximum Playing Speed for Every Frame On** button described above.

**Real Time** This command corresponds to the  **Real-Time Playing Speed On** button described above.

### 6.3.2 ND Views

There are several views which can display the ND2 files in various ways. Some views are available for some dimensions only. If a view is available for two or three dimensions of the ND2 file, a pull-down menu appears in the top image toolbar. There you can select the dimension to be displayed.

Note

*When you switch the view, a new image window opens by default. This behavior can be changed within the **Edit > Options**  command window so that only one view of an image will be opened at a time.*


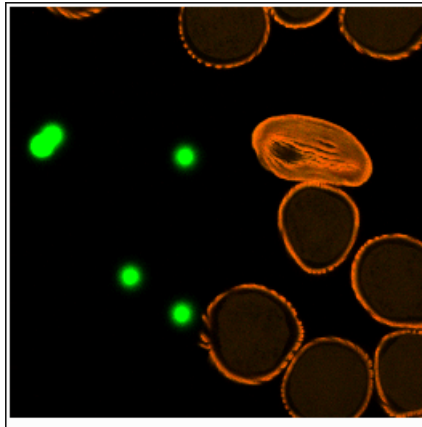
 **Main View** When you open an ND2 file, it opens in this view.

Figure 6.12. Main View




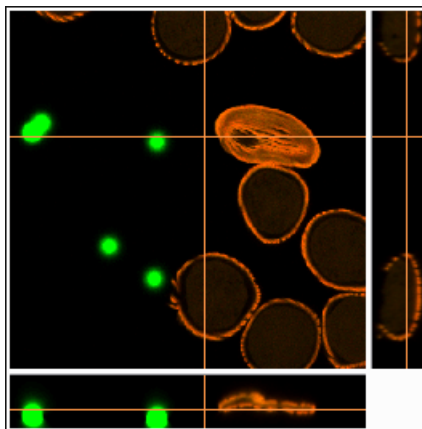

 **Slices View** (requires: EDF Module (Extended Depth of Focus)) This view displays orthogonal XY, XZ, and YZ projections of the image sequence. (Requires Z or T dimension)

Figure 6.13. Slices View



### 6.3.2.1 ND2 Information

Information about current frame of an ND2 document or live image can be displayed using the *Show ND Information* command from the image contextual menu. It displays information about current frame of the ND2 document or live image. All possible information (metadata) are displayed by default.

You can edit which information is displayed in the *Label Properties* dialog window. Right click the ND2 information field and select *ND Info Properties...* Select displayed features from the *Available ND information* area and use the  button to insert them into the *ND information text scheme* area which defines the visible information. Use

*Format options* tools to specify the data format (e.g. numeric precision, date format, etc.) and adjust the text appearance in the *Text Box* tab. *Always show Acq. Time* forbids removing the Acquisition Time information from the ND information text box.

To view the physical zoom of your image (how many times the image is magnified on your monitor compared to the real dimensions of the calibrated sample), click on *Show Physical Zoom* in the context menu over your image. The magnification is placed as an overlay information in the corner of your image.

## 6.4 Large Images

### 6.4.1 Opening Files in Progressive Mode

If a single image (or one ND2 frame) is too big and it can not be loaded to RAM in one piece - you will be asked whether to open it in a *Progressive Mode*. This means that only a thumbnail of the image will be loaded to the RAM. When zooming-in to view image details, the image data of the particular portion of the image will be loaded to RAM progressively (piece by piece).

Many image processing functions and commands are disabled in this mode. This problem usually occurs with *Large Images* created by stitching several frames together. In case you need to process such an image, a smooth solution is to split it into tiles and create a multi-point ND2 file out of it. Please see 6.4.2 Splitting Large Images.

### 6.4.2 Splitting Large Images

You can split an existing large image to tiles easily.

- 1) Open the large image to be split.
- 2) Run the *Image > Split Image* command. A dialog window appears.
- 3) Set split options:

**Split to separate files - output format** Select the output format (ND2 or TIFF). A separate file will be saved for each tile. Define *Output folder* and *Prefix* where the files will be saved.

**Create multipoint ND Document** Only one multi-point ND2 file will be created and opened after the split.

**Set measurement frame after splitting** The measurement frame will be turned ON after the split in order to exclude overlapping areas. This is to prevent distortion of data in case of measurement.

**Tile Size, Number of tiles** Define width and height of a single tile in pixels or set the number of columns and rows to which to split.

**Overlap** Set overlap value of neighboring tiles either in % or the current calibration units (presumably  $\mu\text{m}$ ). Overlap up to 50% of the tile size is allowed, higher values are reduced to the maximum.

**Fill background** Select a color to fill empty spaces which will appear the total area of tiles exceeds the image. Use **Optimize** to prevent the empty spaces.

- 4) If you do not require the tiles to have exact size, click **Optimize**. The system will arrange/resize the tiles automatically so that the whole image is covered precisely.
- 5) Click **OK**.

### 6.4.3 Locating XY Positions Between Images

If two images have matching XY coordinates you can locate the exact position in one image inside the other. This would be the case especially after using the *Image > Split Image* command. Right click the location within the first (multi-point) image and select *Find this Point in Paired Document* from the context menu.

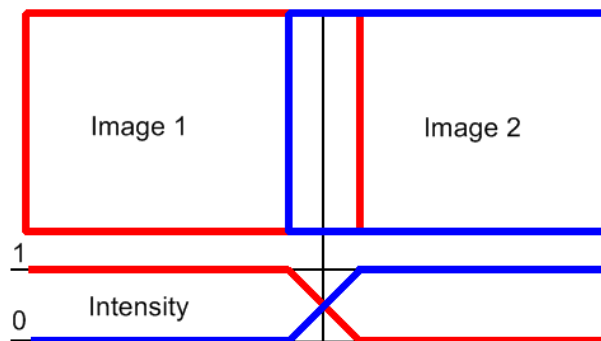
If you do this right after splitting the large image (the two images are “paired”), the system displays it and highlights the XY position by a flashing cross. If it is not clear to the system which images shall be paired (e.g. three or more images with matching coordinates are open), the user will be prompted to select the second image to the pair.

### 6.4.4 Methods Used for Stitching Large Images

Some *NIS-Elements* functions for creating large images by stitching multiple image frames enables the user to select a stitching method. Select the one you serves better your purpose:

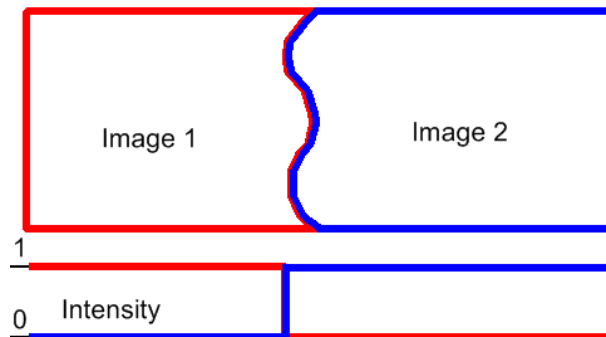
**Blending** Overlapping image parts will be blended.

Figure 6.14. Stitching via "Blending" (pixel intensity is changed in the overlapped region)





**Optimal Path** A contour where two overlapping images are least different will be computed. The images will be stitched copying this contour.

Figure 6.15. Stitching via "Optimal path" (pixel intensity is preserved)



## 6.5 LUTs - Non-destructive Image Enhancement

### 6.5.1 Introduction to Look-Up Tables (LUTs)

LUTs is a useful tool for image color and brightness modifications. You can use LUTs to enhance images for observation purposes so the color modifications will be non-destructive to the image data. LUTs settings are saved along with the image file. If required, the LUTs settings can be applied to the image data by pressing the  button. Use the  Show LUTs window button in the image toolbar to display the LUTs window.

The following actions are available within the LUTs window:

- Adjusting the *Gamma* parameter. The gamma curve is drawn in gray over each graph. You can drag the middle point and move it up or down, or enter an exact value to the G: field at the top of the graph area.
- Adjusting the input intensity range. The input intensity range can be restricted by moving the black and white triangular sliders to the center. All pixels with values outside of the modified input range will be set to maximum/minimum values and the remaining pixels will be approximated in order to fill the output (full) range.

#### Note

*This is the way how images can be equalized. For example, if the image is very dark and most of the histogram is located to the left while the right part of it is flat, you may move the white slider(s) to the position where the histogram line starts to rise. This will incredibly brighten the image and reveal details which were previously hidden in a dark.*

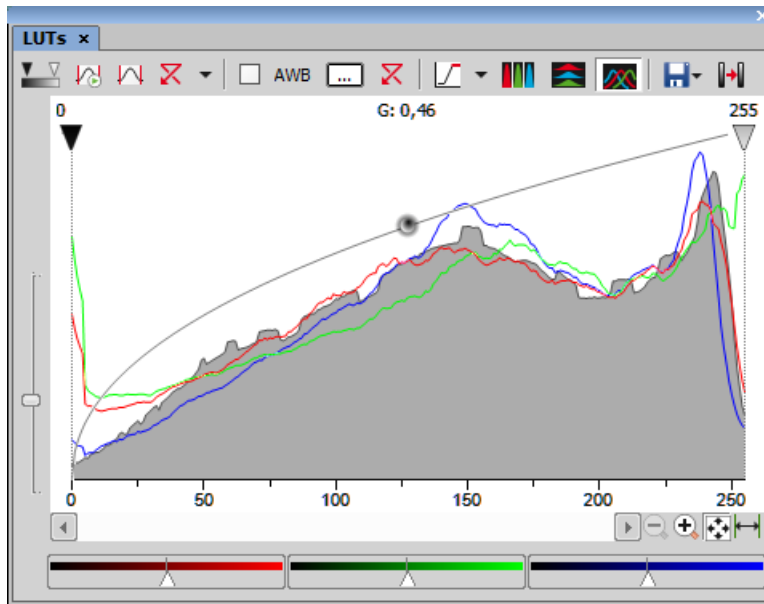
- You can adjust the histograms height non-proportionally by moving the slider on the left side of the window up and down.

#### Tips

- Right-click the graph area and (de)select *Draw trend style*. When ON, the LUTs curves will be smoothed to display the data trends rather than represent the actual image data values.
- The position of the black, white and Gamma sliders can be reset by a double click.

Different controls will be available when applying LUTs to monochromatic, RGB, multi-channel or spectral image. When LUTs are active, the LUTs button in the top-left corner of the image window is highlighted red.


## 6.5.2 LUTs on RGB Images




There are 3 separate windows for each of the RGB channels. All channels are controlled at the same time by default. To control each channel separately, press and hold the **Shift** key while you drag the sliders or adjust Gamma.

### LUTs Tools

 **Enable/Disable LUTs** This button applies LUTs to the current image.

 **Keep Auto Scale** Press this button to run the auto scale procedure permanently (on the live image). When you turn this button OFF, the settings remain as if the *Auto Scale* button was pressed only once.

 **Auto Scale** This button adjusts the white slider position of all channels automatically with the purpose to enhance the image reasonably. If you have selected the *Use Black Level* option from the *Settings* pull-down menu, the black slider will be affected too.

 **Reset All Components** Discard all LUTs settings and turn LUTs OFF by pressing this button.

**Settings pull-down menu** Display the pull down menu and select one of the following commands:


- *Brightfield Opacity* - You can change opacity of a brightfield channel. Define the opacity in the *Brightfield Opacity* dialog window that appears after you run this command.
- *Use Black Level* - Check this item to ensure that the black slider will be affected by the auto scale functions.
- *Settings* - opens the *Auto Scale Settings* window. See 6.5.2.1 Auto Scale Settings.


**Keep Auto White Balance Check Box** Check this option to switch the Auto White Balance functionality permanently (on the live image).


**AWB** Press this button to perform the Auto White Balance operation once.


**Auto White Balance Color (...)** This button Opens the *AWB Color* window and enables you to select a color shade which the system shall eliminate (make white). See 6.5.2.2 AWB.


 **Reset AWB** If *Keep Auto White Balance* or *Auto White Balance* function is applied, this button turns it off.

 **Color Oversaturation** Switch this button ON and the system highlights all pixels with values reaching maximal values. Select colors to highlight the pixels with in the pull-down menu. See also 6.2.3 Channel Coloring.


 These buttons arrange histograms of single channels next to each other horizontally or vertically.

 This button arranges histograms of single channels overlapped. Use sliders at the bottom of the graph window to amplify or reduce the green/red/blue component display in the view.

 **Save/Load LUTs** This pull-down menu enables user to handle LUTs settings in various ways. They can be saved to a \*.lut file and loaded later. Or, LUTs settings saved along with an image can be loaded directly from this image by the *Reuse LUTs from File* command. Or you can copy/paste the settings within the documents opened in *NIS-Elements*.

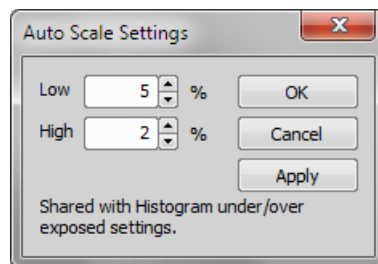
**Modify Image Through LUTs** Press this  button to apply the LUTs settings to the image data - the original image will be overwritten. Until you press this button, no changes are made to the image data.

 **Reset Zoom** Zooms the histogram to fit the preview window.

 **Auto Range** Zooms the histogram so that the “high” and “low” limits are distinguishable. For example, if a small intensity range is defined on a 16bit image the low and high lines are displayed as one-pixel line. Pressing this button will stretch the histogram in order to display the lines separately.

### 6.5.2.1 Auto Scale Settings

Press the arrow next to the *Reset LUTs* button and a pull-down menu appears. Invoke the *Settings* command.




The *Low* and *High* fields determine how many of all pixels of the picture are left outside the sliders when *Auto Scale* is applied (0-10%).



### 6.5.2.2 AWB

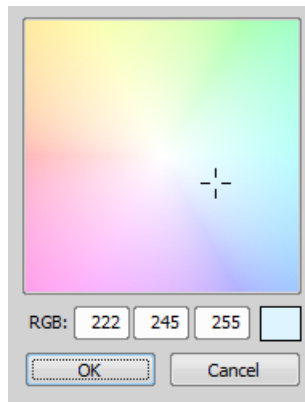White color in some images may suffer from having a color tint. The AWB (Auto White Balance) mode adjusts the image to get pure white instead of this tint. Similarly to LUTs auto scale, the AWB function can be used once, or permanently on the live image. If you know the tint your “white” color has, you can select this color by the color picker that appears after pressing the ... button:




Choose the color by mouse. You can also enter the RGB values to the fields below. A preview of the chosen color appears in the rightmost rectangle.

### 6.5.3 LUTs on Monochromatic Images

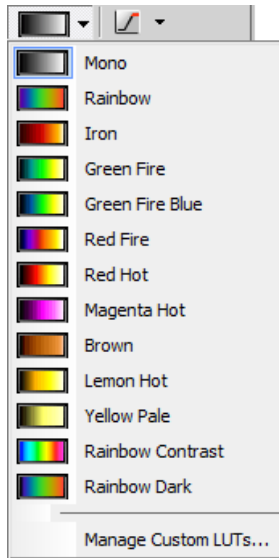All features mentioned above are also valid for monochromatic images, except that the AWB function is not available. There are some other features added when in “mono” mode.


**Gradient** Mapping monochromatic images to pseudo-color gradients is often used to enhance details in the image which would otherwise not be obvious. This button indicates the currently selected gradient. When pressed, a pull-down menu appears where you can select the color scheme you would like to use. Try a few of the schemes to see which one suits best.



**Manage Custom LUTs** The *Manage Custom LUTs* command in the pull down menu enables to define a custom LUTs gradient. Use the tools available in the dialog window to define your custom gradients. They then appear in the pull down menu ready to be applied. Please see *View > Image > LUTs > Create Custom LUTs* for details about the custom LUTs definition.

## 6.5.4 LUTs on Multichannel Images

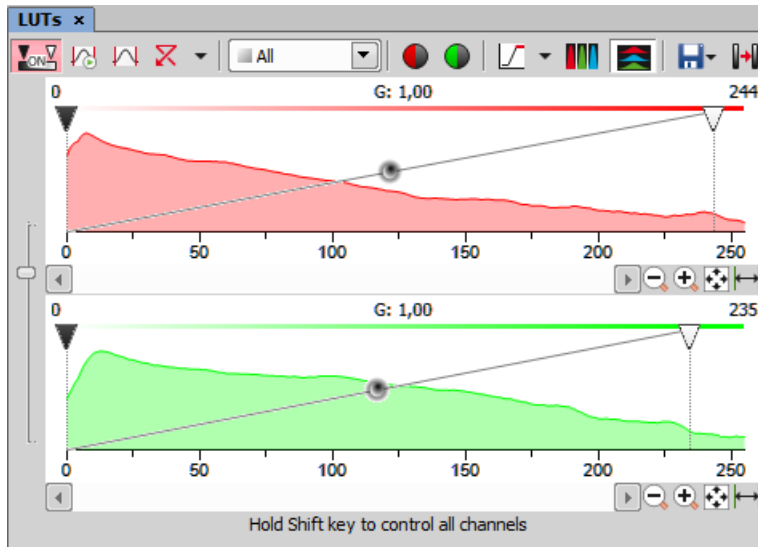Figure 6.20. LUTs window on two-channel image



Most of features mentioned above are also valid for multi-channel images. There are some other features added when in “multi-channel” mode. Each channel is controlled separately by default. To control all channels at the same time, press and hold the **Shift** key while you drag the sliders or adjust Gamma.

When the image contains up to three channels, all channels can be visible at the same time and you can handle them similarly to the RGB mode. When there are more than three channels, only one channel is displayed at a time. A pull-down menu in the toolbar enables you to select the channel to be displayed.

● Next to the channel selection pull-down menu, there is this *Auto Scale* button. By pressing it, the system will automatically adjust only the current channel settings.

## 6.6 Organizer

### 6.6.1 About Organizer



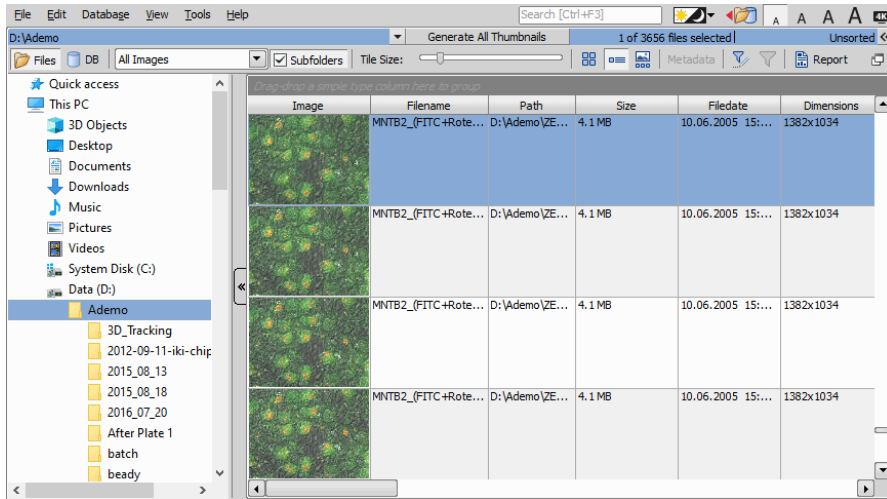
Apart from the main application layout used for capturing and image analysis, *NIS-Elements* provides an extra layout called *Organizer*. *Organizer* makes the work with image files and databases as easy as possible. To activate it, run the **View > Organizer Layout**  command or click the  button located in the top right corner of the application window. The screen opens divided into two identical panes. To switch from one pane to the other, use the **F6** key. To copy files between the panes, simply drag the images from one side to the other side.

Figure 6.21. Organizer layout



- To navigate to a folder containing your image files, use the tree view on the left or enter a path into the first blue edit box. The drop-down menu next to it (▼) reveals a list of recently used folders.  
`Generate All Thumbnails` calculates image thumbnails for all images present in the current folder.
- The next pull-down menu sets the file type. Only images with the selected type will be displayed or *All Images* can be shown.
- If the *Subfolders* check box is selected, all images from the included sub-folders are displayed.

## 6.6.2 Image Filter

You can define a filter which enables you to display only images which fit the defined conditions. One or two conditions may be applied.

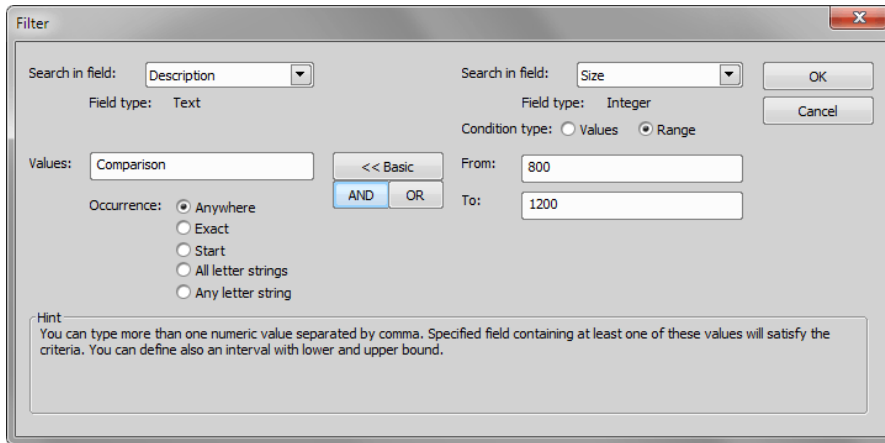
- Pressing this button invokes the filter setup window.
- This button activates the filter.

### **Basic Mode**

This mode enables you to filter files according to one condition.

### **Advanced Mode**

This mode enables you to define two conditions with a relationship between them. Either select *OR* to display files matching at least one condition, or select *AND* to display files which match both conditions.




1. *Search in field* - select the field where *NIS-Elements* should search for a given expression.
2. If the selected field is of a numerical type (e.g. *Size*, *Calibration*, *File date*, etc.) you can specify, whether you want to find the exact *value* or a value in a given *range*. This is selected by the *Condition type* radio button.
3. If the field type is *Text*, the *Occurrence* setting determines the way of evaluating the expression:
  - *Anywhere* - If the given sequence of characters is found anywhere in the field, the system will evaluate it as a match. For example: you have entered “set” to the *Values* field. The filter will select records with the following field values: “set”, “reset”, “settings”, “preset”, etc.
  - *Exact* - If the given sequence of characters exactly equals to the content of the field, it is evaluated as a match. Fields containing the “set” value only will match.
  - *Start* - If the entered string is found at the beginning of a field, it passes the condition. For example: fields containing “set”, “settings”, “setup” will match.
  - *All letter strings* - It is possible to search for more expressions. These should be entered separated by commas. If you want to enter an expression with a space, insert it quoted. If this option is selected, only records where all of those expressions appear (anywhere) will match.
  - *Any letter string* - This option is for entering multiple expressions as above, but this time every field with an occurrence of at least one from the given expressions is matched.

### 6.6.3 Operations with Images within Organizer

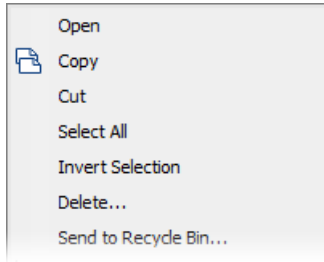
- To open an image double click its thumbnail. *NIS-Elements* will close *Organizer* and display it in the main window.
- To select multiple images, either click on the first and the last image holding the **Shift** key (continuous group selection) or click individual images with the **Ctrl** key down.
- You can copy one or more images from one folder to another by “drag and drop”.
- To delete selected images press the **Delete** key.


All these operations and some other can be invoked also from the context menu, which appears after you right-click on the image thumbnail:





### Thumbnail Displaying Options

You can adjust the way images are shown in the organizer.

 **Show thumbnails** Click this button to display image thumbnails and use the *Tile Size* slider to adjust their size.


 **Show image details** This mode displays images below each other with all available information aside.

 **Show film strip** Images are shown as a film strip below the currently selected image which is displayed big.

 **Metadata** Displays a metadata summary list related to the selected image. This mode is disabled for the image details view.
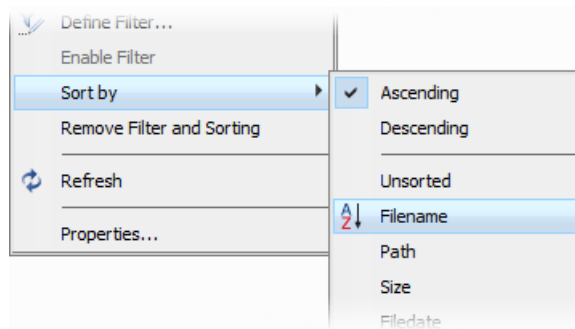
### Sorting of Images



To order displayed images, right click anywhere in the pane - a context menu will appear. Move to the *Sort by* sub-menu which offers several ordering criteria. If the ordering is turned ON already, the icon is displayed on the left side of the applied criterion.



### Grouping of Images



To arrange the view of images efficiently, you can use the capability of grouping of images. Drag the column caption to the grouping bar right above the table. All files with matching field values of the selected column will be grouped together. This can be undone by dragging the column caption back to the others. See Figure 6.21, “Organizer layout” (the Calibration column is grouped).


Note

If you have the Database module installed, the Organizer becomes switchable between two modes: the  Files view and the  Database view.


## 6.6.4 Resizing the Organizer Panes

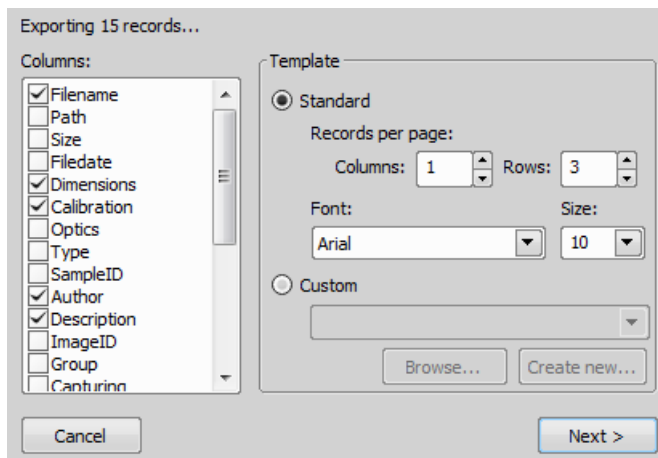
The pane size is adjustable. To resize it, you can either:

- 1. Place the mouse cursor over the dividing line in the middle.
- 2. The cursor becomes an arrow with two tips.
- 3. Drag it left or right to the new position.
-  Resizes the panes to achieve the same size for both of them.
-  Resizes the pane to its maximal/minimal size (one pane is then displayed on the whole screen).

To change the size of the whole organizer window and restore it down, click  *Restore Organizer* ( **F10** ).

## 6.6.5 Reporting

Images together with their metadata can be exported into a report. Selecting image(s) and clicking on  **Report** opens the *Create Report from File Listing* dialog.



Here you can select which columns will be transferred to the report and which template is used. Click **Next >** to generate the template and then adjust it according to your needs. For more information about reports, please see 9. Creating Reports.




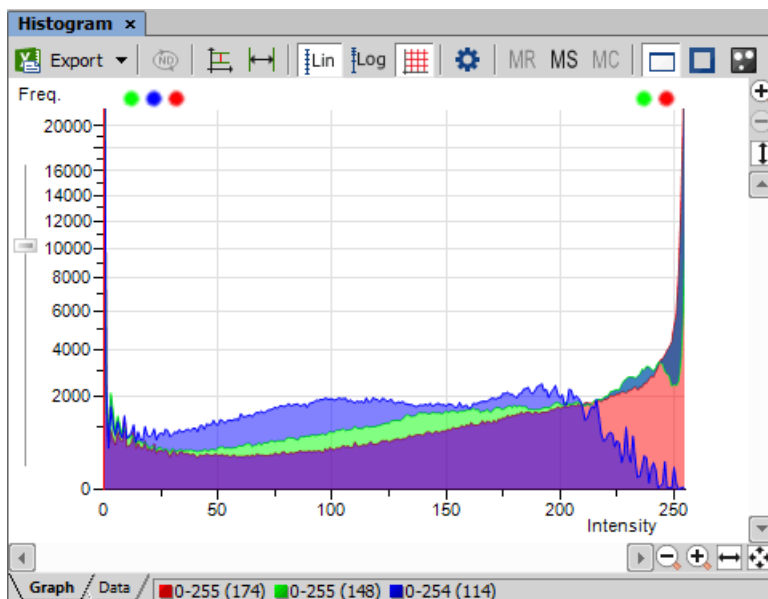


## 7. Image Analysis

### 7.1 Preparation

#### 7.1.1 Histogram

A histogram is a graphical representation of pixel intensity distribution in the image. It is a basic tool which is used during all phases of an experiment to assess the image information. It is especially useful to make sure there are not too many over- or under-saturated (overrun) pixels. Run the *View > Visualization Controls > Histogram*  command to display the histogram control panel:





##### 7.1.1.1 Source Data

Source data of the histogram can be viewed if you swap to the *Data* tab in the bottom-left corner of the window. The area from where the histogram is calculated can be restricted according to the selected histogram mode:

- The data are obtained from the whole image.
- The data are obtained from within the probe.
- The data are obtained from within the current ROI.
- The data are obtained from the image parts under the binary layer.

### 7.1.1.2 Export

 The source data or the histogram image can be exported to an external file. Display the *Export* pull-down menu and select a suitable destination. Click the button to perform the export. The *Export ND histogram*  button enables to export histogram data of all frames of the current ND2 file. The same destination which is selected in the *Export* pull-down menu is applied.


Note


*Please, see the 8.1.7 Exporting Results chapter for further details.*


### 7.1.1.3 Histogram Scaling

The histogram can be zoomed in and out using the zoom buttons on sides of the window. There are also other options to adjust the graph appearance:


**Left slider** There is a slider along the Frequency axis. Drag it in order to stretch the view of either lower or upper values of the axis.

 **Auto Scale Vertical** Zooms the graph of each channel separately to fit the available area. When this function is ON, the histogram is not proportional.

 **Auto Scale Horizontal** Zooms the graph so that the marginal zero frequencies, if there are some, are excluded from display.

 **Graph Linear** Displays the linear scale on the Y axis.

 **Graph Logarithmic** Displays logarithmic scale on the Y axis.

 **Show Grid** Turns on/off the grid in the background.

### 7.1.1.4 Overrun Indication

If certain amount of pixels with maximum/minimum intensity values (black/white pixels) are spotted in the image, color dots appears above the graph. The color indicates which channel is affected. When you place the mouse cursor over the dots, a tool tip message appears with details about the percentage of under/overexposed pixels. The percentage limit setting is loaded from the LUTs settings (the color dots do not appear unless the set percentage of black/white pixels is exceeded).

Note

*If the overrun concerns more than three channels (considering multi-channel images), only one white dot is displayed instead of many color dots.*

### 7.1.1.5 Graph Memorizing

You can save the current graph to memory and display it later for comparison with another graph.

MR Displays memorized graph.

MS Memorizes the current graph. Only one graph can be memorized at a time. If any graph has already been memorized before, it will be overwritten by the current graph.

MC Clears the memory.


### 7.1.1.6 Histogram Options

#### **Drawing Style**

There are two ways the histogram can be drawn:

- *Raw Data* - tries to draw the source data precisely to the graph
- *Trend Style* - interpolates the data so that the histogram lines appear smoother.

#### **Histogram Options Window**

The graph appearance can be modified. Press the   button - a window appears where the following settings can be adjusted:

**Colors** Graph background and axes colors can be selected.

**Pen Width** Set the width of the histogram(s) line to 1, 2, or 3 pixels.

**Fill Graph Area** The area below the histogram line can be filled with the channel color.

**Graph Area Opacity** Select the opacity of the Graph Area color in %.

**Vertical/Horizontal AutoScale, Show Grid** These options equals the corresponding buttons of the histogram toolbar.

**Interpolation method** Select the way of drawing the graph line. The *Linear* (smooth) and *Quick* (precise) options are available.

**Show grid** Shows the grid of the graph.

**AntiAlias** Smooths the edges of the graph line.

**Horizontal axis always visible** If checked, the axis does not leave the graph area while zooming in the graph.

## 7.1.2 Image processing

### 7.1.2.1 Processing On: Intensity/RGB/Channels

If you work with RGB images, you can decide whether to perform the processing on single RGB channels of the image, or on the intensity component. The RGB/Intensity setting is global so it is applied to all the other processing commands automatically. When processing a multichannel image, this option is disabled - the operation is applied to channels automatically.

Note

*In case the command does not open a window where it would be possible to make the choice, the setting from the last performed processing command (the global setting) is applied.*

## 7.2 Image Segmentation

### 7.2.1 Thresholding

(requires: MEAS (Automatic Measurement))

#### 7.2.1.1 A Basic Way to Threshold an Image

Specifying correct threshold limits is a crucial procedure of the automated image analysis. The point is to determine which pixels will and which will not be included in the binary layer and thereby distinguish objects to be analyzed from background. Thresholding can be performed in the following modes:


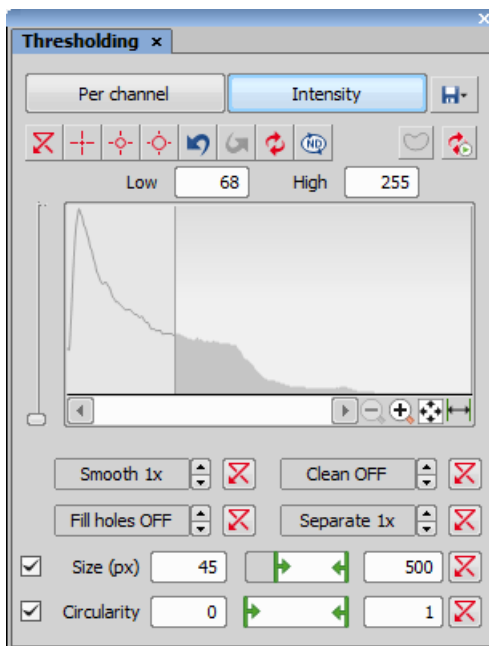

- 1) Open or capture the image to be thresholded.
- 2) Run the *View > Analysis Controls > Thresholding*  command to display the *Thresholding* panel:

Figure 7.2. Thresholding panel



- 3) In the top toolbar, select the color mode which will be used for thresholding. According to the type of image opened, the following modes could be available: *RGB*, *HSI*, *Intensity* and *MCH*. See 7.2.1.2 Color Modes.

- 4) Select one of the pixel-picking tools (e.g. ) and start clicking on objects in the image you want to detect. Color values of each of the clicked pixel(s) are used to adjust the low and the high threshold values so that the clicked pixel would stay inside the thresholded interval.  
Apart from this “clicking” method, the thresholds can be adjusted directly within the histogram by dragging the vertical lines which highlight the thresholded interval.
- 5) All pixels having their numeric values inside the thresholded interval become a part of the resulting binary layer.
- 6) Very often, also other binary objects appear in the image which you did not intend to detect, for example dust particles with intensity similar to the specimen, etc. To eliminate these mis-detected object, use one or more tools available at the bottom of the panel. You can either apply some morphology functions to the binary layer or apply size or circularity restrictions to detected objects. See 7.2.1.3 Tips and Tricks.

### 7.2.1.2 Color Modes

**RGB** The *RGB* thresholding mode is available only if an RGB image is opened. There are two ways how to define the threshold limits, by choosing reference points within the image, or by defining the limit values for each color channel.

**HSI** The *HSI* thresholding mode is available only if an RGB image is opened. The thresholding procedure works the same way as in the RGB mode except that the pixel values are displayed in the HSI (Hue, Saturation, Intensity) color space and that the *Saturation* and *Intensity* channels can be switched OFF. Therefore you can threshold over the *H*, *HS*, or *HI* channels only. Switch the channel off by deselecting the check box in the top-left corner of the channel histogram.

**Per channel** In this mode, the channels are handled separately, a binary layer is created for each channel and also, the binary processing functions (*Smooth*, ...) can be adjusted differently for each channel. Switch between the channels by selecting their names in the bottom pull-down menu.

This mode is dedicated to perform threshold on multichannel images, but can be applied to RGB images as well. It is not available when working with a monochromatic (single channel) image.

**Intensity** The thresholding procedure works the same way as in the RGB mode except that it is performed on the intensity pixel values.

### 7.2.1.3 Tips and Tricks

#### **Threshold Adjustments**

Thresholding parameters can be adjusted very precisely in the following way:

- Threshold limits of each channel can be adjusted by rewriting the values in the top-left/top-right corner of each channel histogram.
- Threshold range of each channel (the colored stripe) can be shifted by mouse. Place the cursor in the middle of the range (a circle appears) and drag it left or right.
- Threshold limits of each channel can be adjusted by mouse. Place the cursor on the edge of the thresholding range and drag it left or right.

---

The two last behaviors can be further modified by pressing additional keys:

- Move the threshold using the right mouse button - the binary layer in the image is not updated continuously, but is updated only once after the mouse button is released. Such operation saves some computing and therefore is quicker.
- Hold **Shift** while moving the threshold limits - the threshold adjustment is performed on all channels together (RGB images only).
- Hold **Ctrl** while moving the threshold limits - both the low and the high threshold limits move in opposite directions.

### **Binary Operations**

Four basic operations can be performed on the binary layer before it is displayed on screen. To turn the operation on, click on the up arrow button in order to define the number of its iterations:

**Clean** Removes small objects from binary image.

**Smooth** Smooths the binary image contours.

**Fill Holes** Fills holes within binary objects.

**Separate** Separates objects.

### **Restrictions**


**Size** Define size range using min/max value in the edit boxes or using the slider.

**Circularity** Define circularity range using min/max value in the edit boxes or using the slider.

### **Thresholding Large Images**

If thresholding images larger than 5000 x 5000 pixels, a new check box called *Preview on selected area only* is displayed. This feature enables to display the threshold preview just on a limited area to speed up the thresholding adjustment.

### **Changing Appearance of the Binary Layer**

To change the color or transparency of the layer, right-click the  *Overlay* button in the right image toolbar. A contextual menu appears, where you can select one of the predefined transparency levels, or run the *Colorize Binary Objects* command. This command displays the binary objects in several different colors. The algorithm ensures that two neighboring objects are never colored by similar colors.


See 7.3.4 Mathematical Morphology Basics.

## 7.3 Binary Layers

### 7.3.1 Introduction to Binary Layers

#### 7.3.1.1 How Do I Get a Binary Layer?


Commands for creating and editing binary layers are placed in the Binary menu. Basic methods of binary layer creation are:

**Thresholding** This is the most common way of creating a binary layer for image analysis purposes. Use the *View > Analysis Controls > Thresholding*  command to specify threshold limits, a binary layer will be created automatically. See 7.2.1 Thresholding.

**Drawing Manually in Binary Editor** A binary layer can be edited or even created manually using the built-in binary layer editor. To enter it, click *Binary > Binary Editor*. See 7.3.3 Binary Editor.


#### 7.3.1.2 How Do I Display the Binary Layer?

There are the following buttons at the bottom of the right image toolbar.

 **View Binary** This button displays the binary layer only of the image.

 **View Overlay** Binary layers can be displayed together with the color layer using an overlay mode.

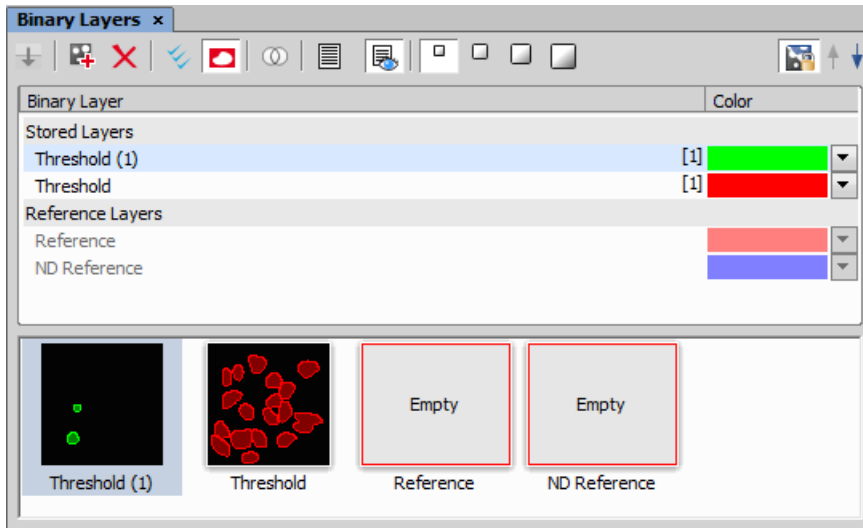
 **View Color** This button switches to view the color image (only).


To manage display of multiple binary layers, use the *View > Analysis Controls > Binary Layers*  control panel. See 7.3.2 Managing Binary Layers.

## 7.3.2 Managing Binary Layers

### 7.3.2.1 Binary Layers Control Panel

Run the *View > Analysis Controls > Binary Layers*  command to display the following panel.




Each time you create a new binary layer, it is automatically placed into the *Working Layers* area of the *Layer List*. Working Layers may be overwritten by the application each time a new binary process is performed. Therefore once you are satisfied with the layer content it is highly recommended to move it into the *Stored Layers* section using the  *Store Current Working Layers*.

### 7.3.2.2 Operations with Binary Layers

**Change color** Color of a binary layer can be changed using the color pull down menu. It offers a list of predefined colors. Besides the list of predefined colors it is possible to create a custom color (*More...*) or to colorize objects in a single layer using multiple colors (*By Object/By 3D Object*).



**Change order** Use the arrow buttons (   ) to change the order of the layers in the list.


**Combine Layers using Binary Operations** If you drag-drop a binary layer name over another binary layer name, the *Binary operations* window appears. The same window can be displayed by the  *Binary Operations Dialog* button.

#### Note

*If multiple layers are selected in the layer list and drag-dropped over another layer, the Binary Operation dialog calculates with multiple (more than two) layers. Available functions are dependent on the number of selected layers.*

See *Binary > Binary Operations*.

**Create 3D Binary Objects** To create 3D objects in the selected binary layer, click  *Connect objects to 3D*. If the binary layers are not defined for all frames of the ND document having more than one frame (e.g. only the current frame contains binary information), you can use  *Fill missing binary frames* to copy the binary from the current frame to all frames where the binary layer is missing.


**Duplicate** To duplicate one or more binary layers, select them by mouse and drag-drop them to the empty area below the layer list, or click the  button.




---

#### Note


If you drag and drop a binary layer to some already existing reference layer, content of the dragged layer is copied to it.

**Make Selection** Binary layers can be selected by mouse (pick more layers holding **Ctrl** or select a range holding **Shift**) or you can select all layers by clicking  *Select All Layers*.

**Remove** Select the binary layers to be removed and click the  button.

**Rename** Double click a layer name to rename it. Only name of the *Reference Layers* can not be edited.

### 7.3.2.3 Attaching Binary Layers to Channels

- 1) If a binary layer is created by thresholding in the **Per channel** mode, such binary layer is automatically attached to the channel on which it has been created. Other binary layers can be attached to a channel manually:
  1. Right-click a binary layer in the *Binary Layers* panel.
  2. Select the *Attach to Component* sub-menu and select to which channel/component, the layer should be attached.
- 2) Visibility of binary layers attached to channels can be synchronized with channels. Turn ON the  *Synchronize binaries with channels* button. Now if a single channel is selected for display, only binary layers attached to it are displayed and vice versa.
- 3) To make an attached binary layer global again, right click the binary layer name or thumbnail and select *Detach from Component*.

### 7.3.3 Binary Editor

(requires: MEAS (Automatic Measurement))

The binary layer, as a result of thresholding, can be modified by hand using the binary layer editor. It is a built-in application providing various drawing tools and morphology commands. Go for the *Binary > Binary Editor* command or press *Tab*. New buttons appear on the toolbars:

#### 7.3.3.1 Drawing tools

The binary image can be modified using various drawing tools. Although the way of use of some tools differs, there are some general principles:

- Make sure you are in the right drawing mode (drawing background **BG** /foreground **FG** )
- Drawing of any object which has not been completed yet can be canceled by pressing *Esc*.
- The polygon-like shapes are drawn by clicks of the primary mouse button. The right button finishes the shape.

- The automatic drawing tools (threshold, auto detect) have a changeable parameter. It can be modified by **+** and **-** keys or by mouse wheel.
- The scene can be magnified by the UP/DOWN arrows when mouse wheel serves another purposes.
- You can drag a magnified image by right mouse button.
- A line width can be set in the top toolbar.
- Hints are displayed in the second top toolbar.



## Binary Layer Color and Transparency

When in overlay mode:



- The **Insert** key switches between predefined overlay colors.
- **Ctrl + Up/Down** increases/decreases the binary layer transparency.

## Erasing Single objects

Single binary objects can be erased in the following way:

- Run *View > Analysis Controls > Binary Toolbar* .
- Select the *Delete Object*  tool from the toolbar.
- Click inside the objects to be erased.

## Multiple Binary Layers

An arbitrary number of binary layers can be created within one image. Click this *Create New Binary Layer*  button in the image toolbar to add a new binary layer. The binary layer that you are currently editing can be selected in the nearby pull-down menu. Binary layers can be managed from the *View > Analysis Controls > Binary Layers*  control panel.

## 7.3.4 Mathematical Morphology Basics










(requires: MEAS (Automatic Measurement))

The binary image as a result of thresholding often needs to be modified before a measurement is performed. Edges of the objects can be smoothed, holes in the objects filled, etc. by using the mathematical morphology commands.

Note

*“Image Analysis and Mathematical Morphology” by J. Serra (Academic Press, London, 1982) was used as a reference publication for the following overview.*

The basic processes of mathematical morphology are: erosion, dilation, open, close and homotopic transformations.

-  **Erosion** After performing erosion, the objects shrink. Marginal pixels of the objects are subtracted. If an object or a narrow shape is thinner than the border to be subtracted, they disappear from the image.
-  **Dilation** After performing dilation, the objects enlarge. Pixels are added around the objects. If the distance between two objects is shorter than twice the thickness of the border to be added, these objects will be merged. If a hole is smaller than twice the thickness of the border, it disappears from the image.
-  **Open** Open is erosion followed by dilation so the size of the objects is not significantly affected. Contours are smoothed, small objects are suppressed and gently connected, particles are disconnected.
-  **Close** Closing is a dilation followed by erosion so the size of objects is not significantly affected. Contours are smoothed, small holes and small depressions are suppressed. Very close objects may be connected together.
-  **Clean** This transformation is also called geodesic opening. The image is eroded first so small objects disappear. Then, the remaining eroded objects are reconstructed to their original size and shape. The advantage of this algorithm is that small objects disappear but the rest of the image is not affected.
-  **Fill Holes** Fills the holes inside objects. This transformation is handy when objects have a rich inner structure with intensities typical for background. After applying the *Fill Holes* transformation, objects become homogeneous.
-  **Contour** This transformation converts objects into their contours.
-  **Smooth** Smooth affects rough edges of the objects makes them smooth.
-  **Morpho Separate Objects** This transformation detects standalone objects that are connected together and isolates them.

#### 7.3.4.1 Homotopic Transformations

Homotopic transformations preserve topological relations between objects and holes inside them. Using a homotopic transformation, an object with 5 holes will be transformed to another object with 5 holes. Two objects without any holes, even if they are very near each other, will become again two objects without holes, but likely with a different shape and size. Opening, Closing, Erosion and Dilation are not homotopic transformations. Typical homotopic transformations in *NIS-Elements* are: *Skeletonize*, *Homotopic Marking* and *Thickening*.

#### 7.3.4.2 Connectivity


Applying the above mentioned transformations has some limitations due to digital images rasterization.

When speaking about binary image processing, a binary image is a set of pixels where values 1 represent objects and values 0 represent background. In the square grid of the image, two possible connectivities can be used for processing - the 8-connectivity or the 4-connectivity. Look at the picture below. If the 8-connectivity is used, the two pixels represent one object. If the 4-connectivity is applied, there are two objects in the picture. *NIS-Elements* works with the 8-connectivity model, so all pixels neighboring by the corner belong to one object.



### 7.3.4.3 Structuring Element = Kernel = Matrix

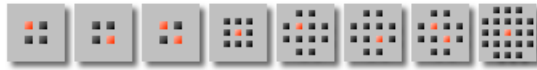When applying Erosion, Dilation, Opening or Closing, one of the parameters which determines the transformation result is the selection of kernel (structuring element, matrix) type. There are the following kernels used in *NIS-Elements*:



The bright pixel in the center or near the center of the kernel represents its “midpoint”.

#### Example 7.1. Erosion

Let's assume 1 and 0 represent object(1) and background(0) pixels of the binary layer. You can imagine the erosion as the following algorithm:

Move the midpoint of the kernel to every point of the image. Each time, look at the neighboring pixels of the kernel and make the following decision:

- If there are object(1) pixels in all the positions of the kernel, set the midpoint to object(1).
- If there is at least one background(0) pixel in the kernel, set the midpoint to background(0).

#### Example 7.2. Dilation

You can imagine the dilation as the following algorithm:

Move the midpoint of the kernel to every point of the image. Each time, look at the neighboring pixels of the kernel and make the following decision:

- If there is at least one object(1) pixel in any position of the kernel, set the midpoint to object(1).
- If there are background(0) pixels in all the positions of the kernel, set the midpoint to background(0).

#### Example 7.3. Open and Close

Open is performed by eroding the image and then applying a dilation to the eroded image. On the contrary, Closing is performed as a dilation followed by erosion.


### Repetition issues

If the midpoint is not in the center, applying erosions or dilations by odd number of iterations causes image to shift by 1 pixel. Normally, the total image shift would be determined by the number of iterations (in pixels). *NIS-Elements* eliminates this shift: it changes the position of the midpoint 1 pixel down-rightwards within the kernel for even operations. For opening and closing it is possible to eliminate this shift totally. However, if you run the erode or dilate processes again and again using the kernel with even dimensions and odd number of iterations, then the shift becomes significant.

### 7.3.4.4 Mathematical Morphology Examples


Please see the following examples of some of the binary functions applied to an image. In the following sequence of images, the functions were applied subsequently:

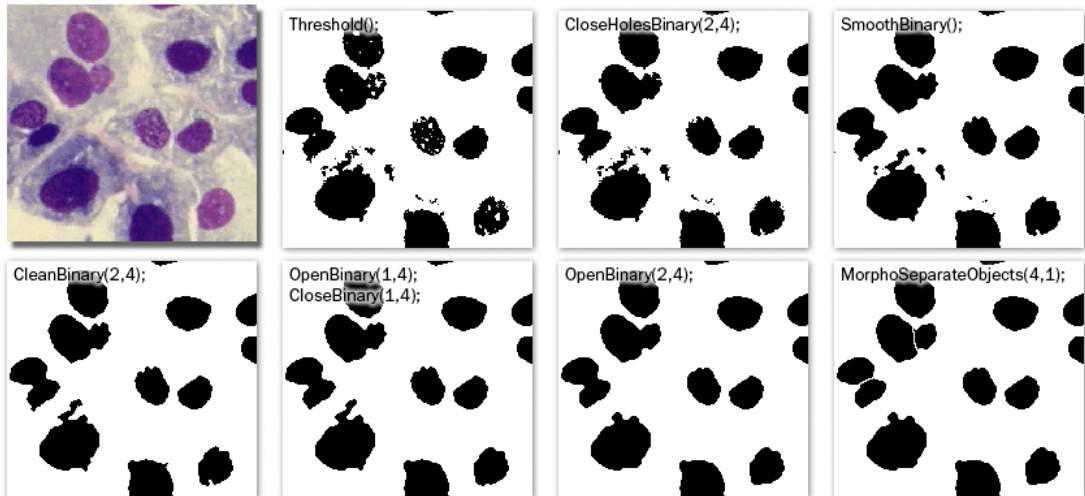Example 7.4. Functions applied subsequently



Example 7.5. Single function applied to the original

The following images were created by calling each of the functions separately:

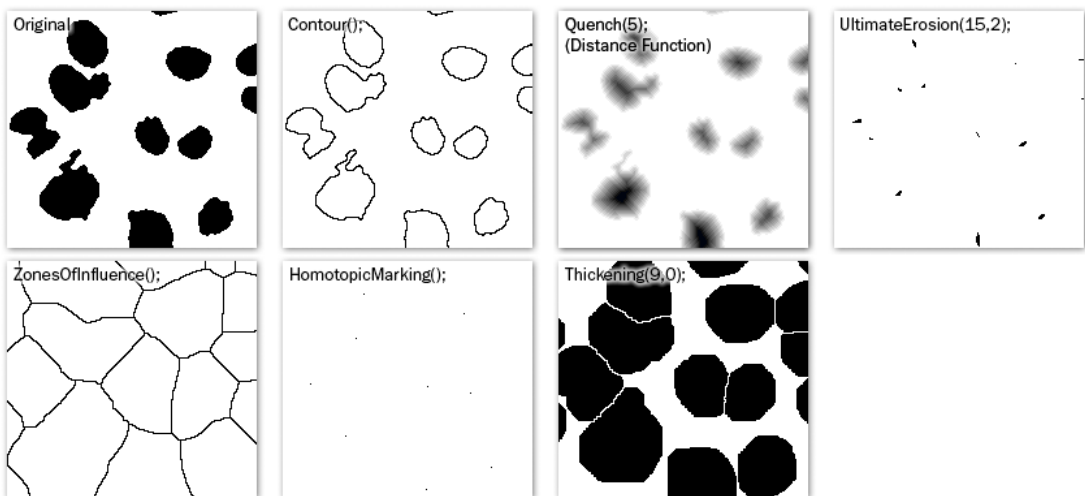

## 7.4 Regions of Interest - ROIs

### 7.4.1 Introduction to ROIs

In laboratory imaging experiments, users are often interested in just a part of the image. To define such an interesting part of the image we use one or more Regions of Interest (ROIs). These regions are used later in analysis and measurements.

#### Note

*Do not confuse image ROIs with Camera ROI which is used to reduce the active area to just a part of the camera sensor (CCD or CMOS). Such reduction usually results in higher frame rate. See 5.2 Camera ROI.*

#### 7.4.1.1 Types of ROI

Individual ROIs (ROI objects) are 2D objects. They can have different shape: rectangular, elliptical polygonal or Bezier. One shape can be created with one or more tools. For instance polygon ROI can be defined using both "Draw Polygonal ROI" and "Autodetect ROI".

In case of an ND file which has time axis (e.g. time-lapse) individual ROI objects can be defined as "Global" or as "Changing over time". Global objects do not change any of their characteristics over time. Objects that are defined as changing over time can change their position and shape in time (however objects cannot change their shape - e.g. from ellipse to polygon). This feature is useful for track moving or warping objects.

In case of Multi Point ND2 file there is one more option: ROI can be defined per Multi Point (which is the default). It means that ROIs from one point are completely independent of ROIs defined on the other point (even if they have the same number). This applies also to ROIs defined as "Changing over Time" (in case the image has time axis too). On the other hand if the ROI is changed to a "Global" it will be shared among all points of the Multi Point set.

Depending on the application purposes ROI objects can have different usage. NIS can handle "Background ROIs", "Reference ROIs", "Stimulation ROIs" and "Standard ROIs".

#### 7.4.1.2 Interaction with ROIs

ROIs can be manipulated anytime directly in the image window. Many of the following techniques work in ROI editors too. Objects cannot be moved or resized when they are Locked (some analysis tools or users may lock them to prevent accidental modification). They must be unlocked (from the ROIs' context menu) before any manipulation. Note that the mouse cursor changes over the ROI to indicate which action will be started after mouse button is pressed (it changes when modification keys like Alt and Ctrl are pressed).

**Selection** ROI objects can be selected by clicking on an unselected object (other previously selected are unselected). Clicking on a ROI object with Ctrl key pressed results in altering the selection state. More than one object can be selected in this way. Double-clicking on any object results in only that object being selected. All objects can be selected by pressing Ctrl+A or by choosing "Select All ROIs" from the menu over any ROI. Alternatively, it is also possible to select a set of objects by dragging a rectangle around them with Ctrl key being pressed.

**Moving** Objects are moved by simply dragging them (cursor indicates that). When more objects are selected, only the one below the cursor position is moved by default. This can be changed by pressing the **Alt** key and move all selected objects. When object is changing over time, any moving results in a new key-frame. Individual dots

representing the key-frames can be moved too. It is possible to duplicate existing objects by Shift dragging. It works on selected objects. In case of a key-framed object, object is duplicated with all its' key-frames unless Alt is pressed.


**Resizing** Resizing a ROI is done by dragging its' contour (the cursor changes accordingly). Note that the hot-zone (the area where mouse cursor changes into different tool) for resizing is outside of the object contour while the hot-zone for moving is inside. Resizing always affects only one object (selection is not taken into account). The behavior slightly varies depending on the object shape. By default resize does not hold aspect of the object - width and height can be changed independently. By holding Shift key aspect is maintained (useful for circles) and the center is not moved. Rectangle can be also rotated by using Alt key while dragging.

Exact values can be entered in "ROI Properties" dialog which is accessible from the ROI context menu.

### 7.4.1.3 Simple ROI Editor

When creating more than one ROI, the Simple ROI editor is the tool of choice. With this tool it is possible to make all possible shapes with different methods. All interaction techniques work here as well.

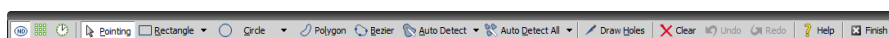Figure 7.8. Simple ROI Editor toolbar



Note that many actions have a hot-key (P for Pointing tool, R for Rectangle, E for ellipse, L for Polygon, B for Bezier, A for Auto detection, H for drawing holes and others) to speed up the process of interaction.

### 7.4.1.4 Auto-Detect Tool

Auto-detect tool is available from ROI menu, Simple ROI editor and some other places. It produces a Polygon ROI. Using the tool is two step process: The first click on an image structure results in preliminary object contour, that is drawn on the screen. After that, user can further tweak the object contour and then finish whole auto-detect by confirming with right mouse button. User can always cancel auto-detect by pressing the **Esc** key.

The very first click is important, because these pixels under the cursor determine intensities from which the algorithm guesses the whole object. As a general rule it is better to choose brightest pixels in the bright objects (fluorescence) and dimmest in the dark object. It is also good to click near the center of the object.

After the click the algorithm presents its guess of the object contour and goes into the interactive phase. In this phase it is possible to:

**Redefine the first point** by clicking again in the image

**Cancel auto-detect** by pressing **Esc** key

**Grow or shrink the object area** (changes the intensity range of pixels making the object) by scrolling mouse wheel or pressing PageUp/PageDown (for finer steps use Ctrl + mouse wheel)

**Erode or Dilate** morphologically the object by 'E'/'D' or Down/Up keys

**Open or Close** morphologically the object by 'O'/'C' keys

**Morpho-separate** portions of the object by 'S' or 'P' key (if it is not possible to separate the key does nothing)

---

#### 7.4.1.5 Using ROIs for analysis

ROIs are used mainly in Automated Measurement to measure ROIs features or number of binary objects inside each ROI, Object Count to restrict binary objects to areas of interest only, ROI statistics to measure intensities interactively. In all these cases ROI selection is taken into account as well as intersection options in Measurement Options.



## 8. Measurement

### 8.1 Basic Workflows

#### 8.1.1 Calibration

##### 8.1.1.1 Introduction

Calibration of the image is of crucial importance to measurement. On calibrated images, realistic measurements can be performed and objects of different images could be compared to each other. A correct calibration has to be made before you begin to measure.

There are two ways how to obtain a calibrated image:

- To capture images with a calibrated system (objective). Please see the 4.1.5.3 Objective Calibration chapter which explains how to calibrate the *NIS-Elements* system.
- To calibrate an uncalibrated image manually:

##### 8.1.1.2 Calibrating an (Uncalibrated) Image

There are two ways to calibrate an image

**Manually** After you run the *Calibration > Recalibrate Document* command, you can specify a distance in the image or set the current pixel size. Please see 4.1.5.4 Manual Calibration

**Using Objective** When you have an uncalibrated image, you can use an already calibrated objective to calibrate it. Right-click the *Uncalibrated* field on the image status bar and choose one of the objectives from the *Calibrate using Objective* sub-menu. Of course, the same objective which was used to capture the image shall be selected.

##### 8.1.1.3 Units

*NIS-Elements* supports the following units: pixels, nanometers, micrometers, millimeters, centimeters, decimeters, meters, kilometers, inches, mils

If the image is uncalibrated, pixels are the only units available. In case of a calibrated image, it is possible to select other units which are then used to display all values (e.g. measured length/area). There are two ways of how to select the desired units:









- Right click the image status bar where the calibration is displayed. Select the units and precision from the context menu.
- Or, click the *Current unit* button located in the *Calibration > Objectives* command window and select the units.

## 8.1.2 Rough Measurement

Quick and approximate measurements can be performed using the *graticule* layer. They behave like adjustable floating rulers. User can simply align a graticule with the measured object and read the distance value (e.g. the diameter). To activate the graticule, press the *Graticules* button.

### 1) Select Graticule Type

The type of the graticule ruler is indicated by a picture on the *Graticules* button placed on the right image toolbar. To change the graticule type, click on the graticules button with the right mouse button and select the appropriate item from the context menu:

-  Rectangular Grid
-  Circle
-  Simple Circle
-  Cross
-  Industrial Cross
-  Simple Cross
-  Vertical Ruler
-  Horizontal Ruler

### 2) Define Graticule Properties


Right click the *Graticules* button and select the *Graticules properties* command from the pull-down menu. A window appears where display parameters (shape, color, line width, density of lines) of all graticules can be adjusted. Please see *View > Layers Properties > Graticules Properties*.

#### Graticules Density

The density value - the closest distance between two line intersections of the graticule - can be set, or you can let *NIS-Elements* adjust it automatically according to the current zoom factor. The units selection depends on the image calibration (calibrated/uncalibrated).

#### Note

---

*Concerning the Cross  graticule. The distance defined by Density is divided into smaller sections automatically. However, for odd density values greater than 10, the markings remain hidden because it is not possible to display them accurately.*

### 3) Measure the Image


The graticule measurement provides the following options in the graticule menu:

- A graticule can be moved around by clicking on it and dragging.
- A binary layer can be created from the current graticule using the *Copy Graticules to Binary* command.
- A *Graticule Mask* can be created from the current binary layer using the *Copy / Add Binary to Graticule Mask* command.
- A *Graticule Mask* can be stored/loaded to/from an external file via the *Save Graticule Mask As* and *Open Graticule Mask* commands.

- Position of the graticule can be reset by the *Move Graticules to Center* command.
- If *Graticule Mask* is set as a graticule type, its transparency can be adjusted using the *Mask Transparency*. In the *Graticules properties*, the opacity can be precisely set, mask can be rotated (*Angle*) and *Calibration* changed.
- A new image containing graticules can be created by the *Edit > Create Full View Snapshot (8bit RGB)* command.

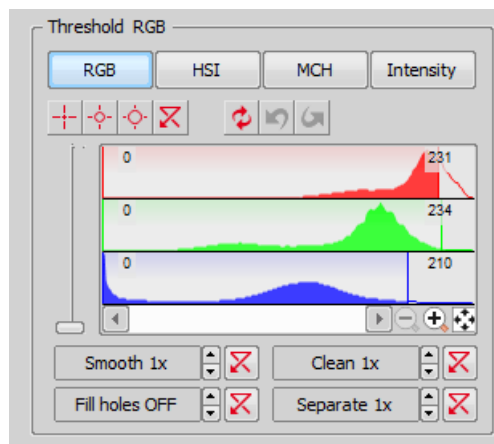
### 8.1.3 Object Count

(requires: MEAS (Automatic Measurement))




The *Object Count* tool is designed for automated object detection and counting. It enables users to threshold the image, automatically measure the binary objects, and export the measured data to a file in a straightforward way. Object count can be performed even on Live image. Run the *View > Analysis Controls > Object Count*  command to display the control panel.

The recommended workflow through the dialog is to move from left to right, starting with thresholding, then using restrictions and finally reviewing the result and export out of *NIS-Elements* to a text file or an MS Excel spreadsheet.

#### 8.1.3.1 Thresholding the Image



The threshold limits shall be defined by picking object-typical colors from the image. Select one of the following tools:


-  1 point threshold tool
-  3 points circle threshold tool
-  6 points circle threshold tool

Click inside the image to define typical object areas. The system will detect similar parts of the image and highlight it by color. The threshold limits are indicated in the histogram and can be further modified by dragging the limit lines by mouse.


Please, refer to the 7.2.1 Thresholding chapter for further details on thresholding.

### 8.1.3.2 Applying Restrictions

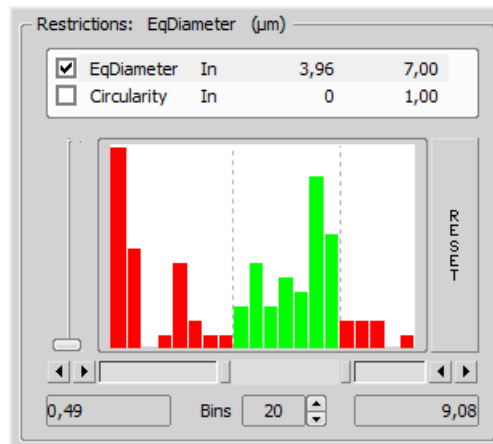The number of objects that are included in the results table can be restricted by defining limits. Objects which do not fit these limits will be left out of the result table.



- Right click to the restrictions field to select one or more of the available measurement features.
- Select the restriction feature you would like to define.
- Move sliders below the histogram to set the limits. The limit values are indicated next to the feature name, and can be modified by double clicking the indicated value.
- Decide whether the defined interval will be excluded or included - this is done by setting the *Inside / Outside* option next to the feature name.
- The nearby check box indicates whether the restriction is currently being applied or not. If applied, the histogram below is color, otherwise it is gray.

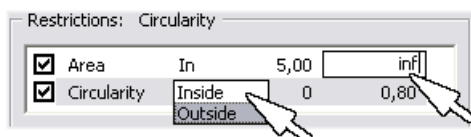

Note

*Infinitude can be defined by entering "oo" or "inf".*


**Reset** The *Reset* button on the right side discards restriction settings of all features.

**Bins** This setting determines the number of columns in the histogram.

### **Restrictions from the Image**

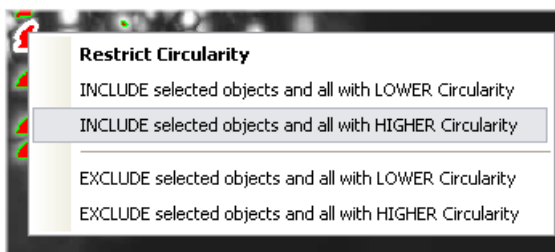A reference object can be picked from the image and used to set the limit values.

- Select the restriction feature you would like to set by mouse. Let's assume you have selected *Circularity*.
- Right click the thresholded object which is going to be used as source. A context menu appears.
- Select whether to *INCLUDE* or *EXCLUDE* the selected object and all others with higher/lower circularity.
- The restriction will be applied and indicated by colors within the image.



### 8.1.3.3 Results

The following results appear in the results table:

**Total Area** The area of the ROI. It is shown in square pixels or square *calibration units*.

**Measured Area** The area of the binaries inside the ROI.

**Number of objects** Number of objects after applying restrictions.

**Table of results** Measurement results for each object are displayed in this table. Select the columns to display/hide from the context menu which appears when you right-click one of the column captions.

**#** Number of selected objects is displayed in the left part of the toolbar. One or more objects can be selected by mouse while holding the Ctr1| or the Shift| key.

Result

Total Area: 0,007 mm<sup>2</sup> **42**

Measured Area: 23,57 %

0 Export  Use Standard EqDia 0,001 mm

ROI ID	Binary ID	#	Area	Perimeter	Circularity	MeanIntensity
Whole Image	Cell Count 3 (In...	1	0,00002	0,02	0,38	84,91
Whole Image	Threshold	1	0,00002	0,02	0,52	178,82
Whole Image	Cell Count 3 (In...	2	0	0,01	0,34	75,04
Whole Image	Cell Count 3 (In...	3	0	0,010	0,69	90,45
Whole Image	Cell Count 3 (In...	4	0,00001	0,01	0,87	87,74
Whole Image	Threshold	5	0	0,01	0,44	178,47
Whole Image	Threshold	6	0	0,01	0,49	186,11
Whole Image	Threshold	7	0	0,009	0,76	183,43
Whole Image	Threshold	8	0	0,02	0,36	118,03

### Toolbar

There is a toolbar above the results table which contains the following buttons:

**Delete selected objects** Removes the selected objects. The objects can be selected by mouse while holding Ctrl.

**Invert selection** Deselect objects selected by mouse and vice versa.

**Generate binary** Erases binary objects which do not meet the restrictions settings.

**Show Object Catalog** Opens the *Object Catalog* control panel with data source automatically set to “Object Count”.

**Export** The measured data can be exported to clipboard or a file using the *Export* button. Please refer to the 8.1.7 Exporting Results chapter for further details.

**Save/Load Object Count configuration** Invokes a pull-down menu which enables you to load/save the current object count settings from/to an external file (\*.counting).

**Use Standard EqDia** Check this option to enable counting of objects using standard *EqDiameter*. Value of the standard *EqDiameter* parameter can be edited in the field nearby.

Note



*Please be aware that the number of objects calculated by this function is only an approximation and not an accurate count of objects.*

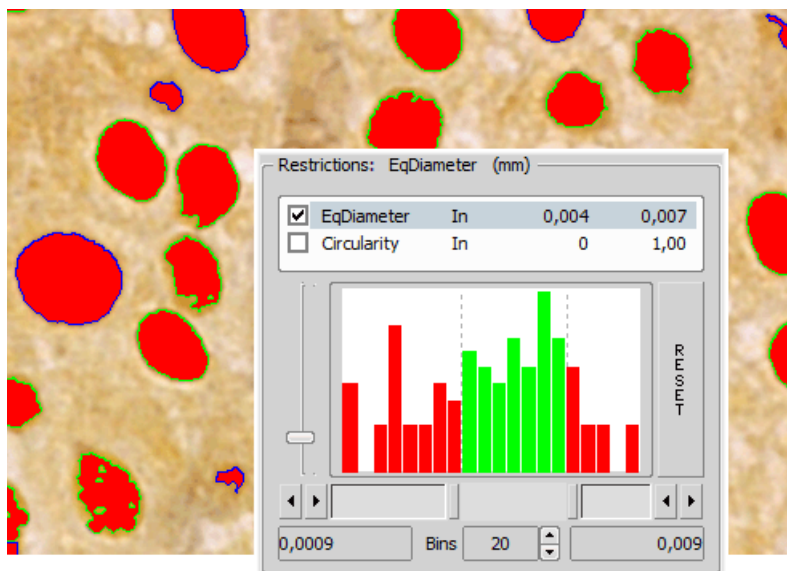
**Keep updating count** Refreshes the measured results each time the binary layer changes (e.g. when the threshold is modified).

### 8.1.3.4 Object Count Procedure Example

1. Open the *Object Count* control panel by the *View > Analysis Controls > Object Count* command.
2. Threshold the image. Use the thresholding picking tools and click on the image to sample areas of interest. These pixels selected will determine what intensity/ part of histogram is considered accepted. The threshold


will detect several specs or smaller objects such as noise or background. If desired, the *Clean* and *Smooth* filters will exclude smaller objects.


3. You can also measure using the defined region of interest (the  button) or measurement frame ( the  button). Check the options are properly set in the *Measure > Options* dialog window.
4. Apply restrictions. Select *Area* from the list of available restriction features. Use the sliders in the histogram to specify the lower limit and the upper limit. The lower and upper limit will update in the display.



The green area of the histogram represents the accepted range areas. Any object with an area in the green range will be considered an object. Alternatively, any object with an area in the red ranges will not be considered an object.

The outlines of objects on the image differ according to whether the object is restricted or accepted. In this case, accepted objects have a green outline and restricted objects have a blue outline.

To ensure that that count is updating with any change, click the  icon:


5. Remove any unwanted object from the object count by selecting the **Delete Selected Objects**  icon from the Result section. The object will lose its thresholded overlay and will be excluded from the measurement.

#### Warning

*You cannot Undo this action. To restore an object repeat the original threshold procedure or load a saved Object Count configuration.*

6. View the result data and export them to a file or spreadsheet application. See 8.1.7 Exporting Results for further information.

## 8.1.4 Measurement Options

Run the *Edit > Options*  command and select the *Measurement* tab.

**Rules for Excluding Objects** This section regards automatic measurement. Select the options how to treat the objects touching the area borders when you are performing the automatic measurement:


**When using measurement frame** Pick up the option how to treat the object touching the measurement frame. Run the automatic measurement by the *Measure > Use Measurement Frame* command.

**When using ROI** Pick up the option how to treat the object touching the ROI frame. Run the automatic measurement by the *View > Layers > Measurement ROI* command.

**Without Measurement Frame and ROI** Pick up the option how to treat the object touching the image border. Turn off the measurement frame and ROI and run the automatic measurement.


### Note

*Rules for excluding objects are valid for all features of objects measurements. For field, frame and ROI measurement, these rules are taken into account only for the number of objects and do not affect the field measurement results. Field measurement is always applied inside the defined area without exceptions.*

**Object Colors** After the automatic measurement is performed, all measured objects are highlighted by color borders. Here you can select colors for the objects excluded from measurement according to the ROI/measurement frame settings. The *Out of limits* color will be used to highlight objects which do not fit the applied restrictions (*View > Analysis Controls > Automated Measurement* ).

### Manual Measurement

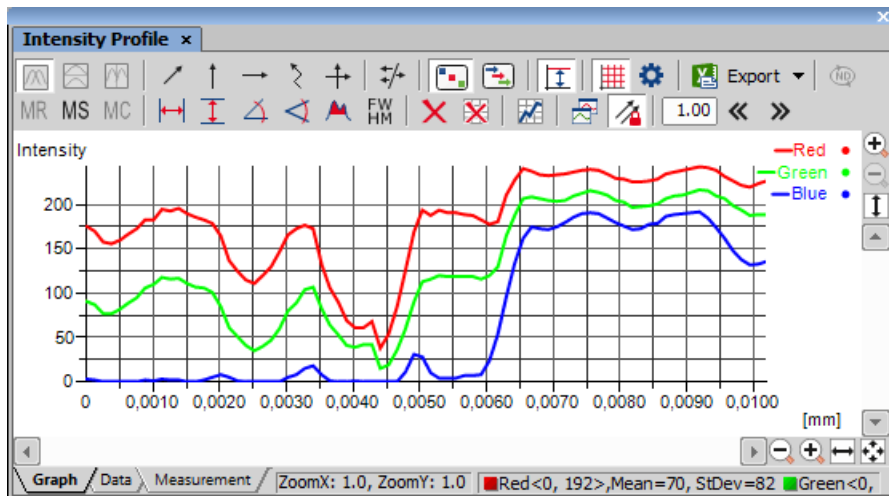
**Automatically zoom in Length manual measurements** This option zooms in while placing the measurement points into the image (click and hold the primary mouse button). Only the indicated measurement tools are affected.

**Copy manual measurement objects from live image to captured** *NIS-Elements* enables to measure manually on live camera signal ( *View > Analysis Controls > Annotations and Measurements*  ). After the *Acquire > Capture* command is used, the measurement objects in the image are copied (or not) to the captured image according to this setting.



**Label** The manual measurement objects can be labeled. Select what information will be attached to every manual measurement object in an image.





## 8.1.5 Measurement on Graph





The following interactive measurements can be performed within graphs.

 **Vertical**,  **Horizontal** Pick this tool to measure distances within the graph.

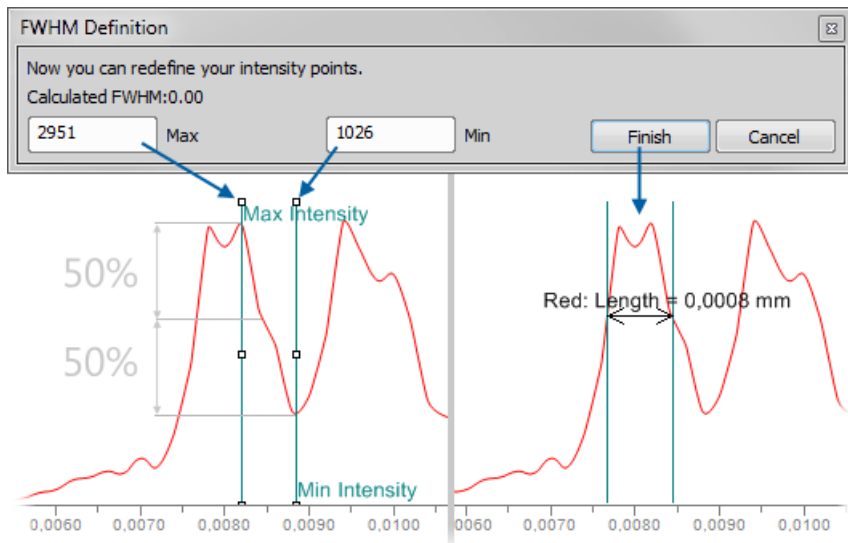
 **Angle** This tool enables selecting an angle and its both line lengths. The resulting number represents a ratio between the projection of the y arrow on y axis and projection of the x arrow on x axis. The resulting number is a dimensionless value. This tool does not allow adjusting the x line length.

 **Free Angle** This tool measures the same features as the previous angle tool but allows to adjust all of the angle's anchor points.

 **Area of Graph** This tool enables you to draw a rectangle in the graph. Area will be measured of intersection of the rectangle with the area below the graph curve.

 **FWHM** This tool measures the *Full Width at Half Maximum* value on the given graph range.

1. Click to the graph to define the maximum, then the minimum.
2. Half of this interval defines the value in which the graph width will be measured.



Note

If the minimum is defined in a different part of graph (e.g.: different peak), the width is measured on the peak where the maximum value was defined.


- ✗ Clear Measurement Objects** Deletes all the measured objects in the graph.
- ✗ Reset Data** Clears the table with the measurement results.


### 8.1.6 Pixel Classifier

A typical task for Pixel Classifier may be counting the area of a cross-section of a blood vessel. Since there are at least three different types of tissue whose area should be calculated while background should not be counted at all.


1. Eliminate the background. You can:
  - Define a region of interest. See 7.4.1 Introduction to ROIs for more information.
  - Use the method of accepted region. Run the *Edit > Region > Region Settings* command first. Choose color of the background in the *Region Settings* dialog that appears. You can fill in the value 0, which is easy to threshold. Next run the *Edit > Region > Accepted Region* command. Draw the boundary around the area to keep. Finish drawing with a right click. The image is cropped according to the boundary. The area left outside the border is deleted and filled with the set background color.

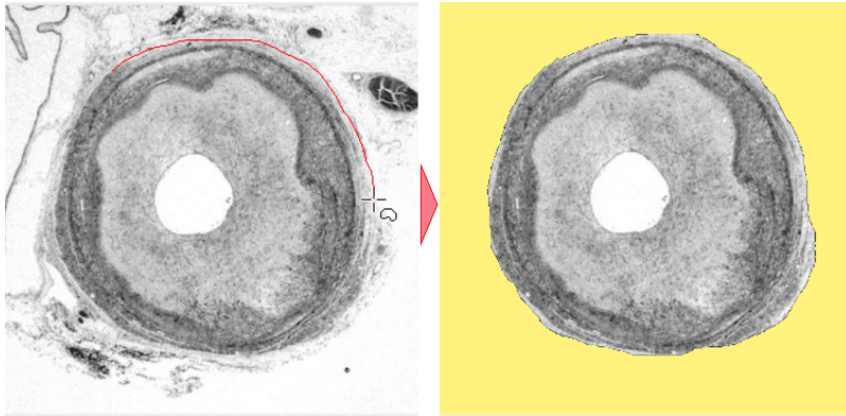Figure 8.9. Accepted Region Example



2. Run *View > Analysis Controls > Pixel Classifier*  to display the classifier panel. Press the *Define* button and edit all phases you want to detect separately. Use the **+** and **X** buttons and commands in the contextual menu to *Add, Rename, Reset* or *Delete* any class.

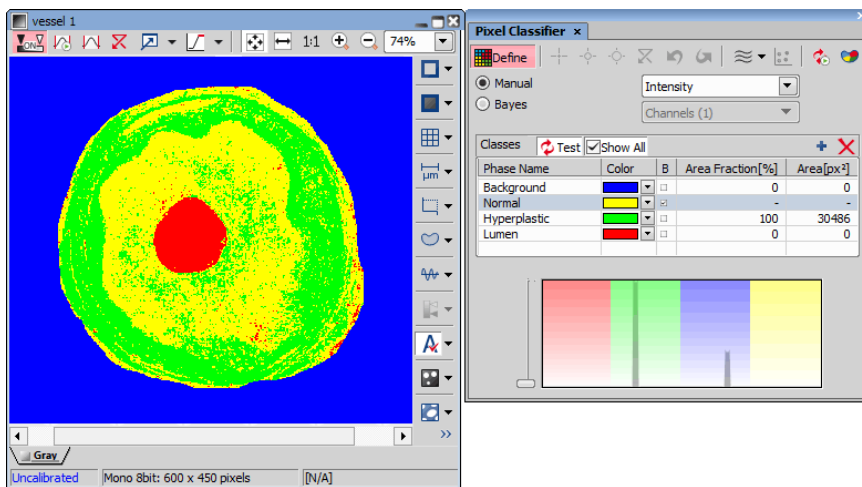If you choose the *Manual* method of classifying, you can manually choose the intensity ranges of all phases.




Alternatively, you can sample areas and train the other algorithms to find phase boundaries.

3. Once all phases have been defined properly, the settings can be reused to classify another images.

## 8.1.7 Exporting Results

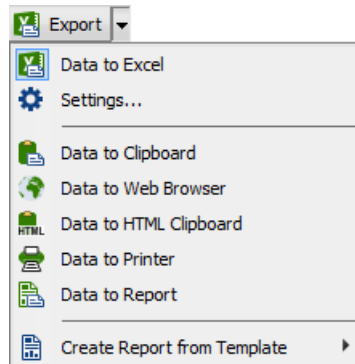Some image analysis data and measurement results can be transferred out of *NIS-Elements* so they could be used further. There is a standard *Export* pull-down menu in some of the control panels (intensity profile, histogram, measurement, etc...) which enables the export.



The set of commands of the menu differs according to the type of data (tables or graphs) which can be exported. Click the arrow button to display the pull-down menu. When you select the destination, the pull down menu hides and the icon of the selected export type appears on the *Export* button. The export action is performed after the button is pressed once more.

### 8.1.7.1 Export Destinations

It is possible to export tabular data or a graph:

**To Excel** (*requires: MS Excel*) Tabular data can be exported to MS Excel. A new XLS sheet opens and the table is copied to it automatically. There may be also the *Export All To Excel* option, which copies the data table, measurement table, and the graph image into Excel at once.

**To File** Tabular data can be exported to an MS Excel sheet (xls) or a text file (tab delimited txt).

**To Raster File** Graphs can be saved as a \*.bmp image file. A save as dialog appears each time so you can name the file properly.

**To Clipboard** Data tables and graphs can be exported (copied) to MS Windows clipboard. Then the data or the image can be inserted into any appropriate application (text editor, spreadsheet processor, graphics editor) typically by the *Paste* command.

**To Web Browser, To HTML Clipboard** Some tabular data can be exported as a HTML table. It can be displayed in a default Internet browser (the *Web Browser* option), or the HTML code can be copied to Windows clipboard - ready to be inserted to any HTML/text editor.


**To Printer** Tabular data can be printed directly. This option opens a standard *Print* dialog window where it is possible to select a printer and print the data.

**To Report** Tabular data and graphs can be exported to *NIS-Elements* report. If a report is already opened for editing, the data/graph will be appended to it. Otherwise, a new report will be created and the exported data inserted to it.

**To Report Template** The *Create Report from Template* option Inserts the result data into an existing report template. Choose an existing report template from the list in the pull-down menu or click on [Browse for Template...](#) to open

a report template file (.rtt) from your hard drive. The report opens and dynamic data are inserted automatically. Then you can adjust the report, save it, print it, or export it to PDF using the menu commands.

### 8.1.7.2 Data Export Options

Run the *Options* command and switch to the *Data export* tab or press the *Settings*  button within the export pull-down menu. There are two tabs available:

#### **Global Settings**

**Export into Microsoft Excel application** Select this option if you want to export the data to MS Excel application. You can choose if the data will be exported into currently opened file or a new file will be created. Or you can set to export the data to a specific file. If this file does not exist, it will be created. Choose the path and name of the file to be created/opened.

Set the *Start at column* values (column and row indexes). These values mark the first cell in the MS Excel application which will hold the exported data.

**Export text files into folder** Select this item if you prefer to export the data to a text file. Press the ... button and specify the target folder and the file name.

**Data delimiter** Pick up the type (tabulator, space, semicolon) of data delimiter which will be used when exporting data to MS Excel or clipboard.

**Append current data to previous exported data** If you check this item the existing data will not be overwritten by the new data.


**Export also column headers if not appending** Include the titles of columns in the files which are being created or overwritten.

**Activate target application after export** If you check this item the exported file is opened in the default application (MS Excel or the default text editor) automatically.

**Insert empty lines** Check this item to allow insertion of empty lines in the exported data.

#### **Data export**

You can modify the amount of data included in each export type. The control panels which contain the Export to file functionality are listed in the top pull-down menu. Select the control panel of which the export shall be modified.

**Export Name** A control panel such as *View > Analysis Controls > Annotations and Measurements*  stores information about the measurement in multiple tables according to the measurement type (area, length, etc.). In this column, you can select the table you would like to modify.

**Sheet name** In this column you can define an arbitrary name which will be used to name a sheet or a txt file to which the data will be exported.


**Options for:** This section of the window enables you to select the items to be included/excluded from the selected table during export.

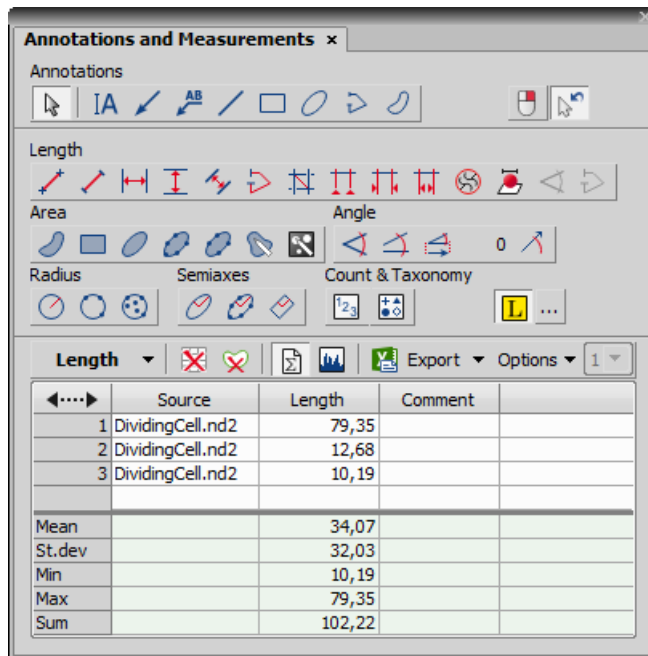
## 8.2 Manual Measurement


### 8.2.1 Basic Workflow

Length, area, angles, taxonomy, counts, circle radius, and ellipse semiaxes can be measured manually over an image. There are several tools for measuring each feature. The results are being recorded to a simple result table, which can be exported to a file or clipboard. The data can be presented as a graph, too.


#### **Example - Measurement of a Grain Thickness**

- 1) Run the *View > Analysis Controls > Annotations and Measurements*  command. The manual measurement control panel appears.



- 2) Select the *Vertical parallel lines*  tool.
- 3) Place the first line on the top edge of the crystal by clicking into the image. The position of the line can be adjusted while you hold the primary mouse button down. After you release it, the line is positioned.

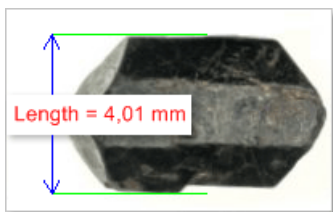- 4) Repeat this to place the second line on the bottom edge of the crystal.




- 5) When finished, an arrow is drawn between the lines, and the result of the measurement is attached. A record with the measurement type and the measured value is added to the results table.
- 6) After finishing the measurement (e.g. of multiple grains or other images), select where to export results in the *Export* pull-down menu.
- 7) Export the results using the *Export* button.


## 8.2.2 Working with the Measurement and Annotation Objects

There are two types of vector objects which can be placed over the image to a separate layer:

**Annotations** Vector objects used to emphasize or label interesting areas in the image, such as an arrow, a text box, etc.

**Measurement objects** After a measurement object is placed to the image the measured value is appended to the table of results. See 8.2.1 Basic Workflow.

**Displaying the vector layer** Use the  *Show Annotations (A)* icon on the right image tool bar to display all vector objects in the layer.


**Select or Clear Objects** Objects of the annotation layer can be selected by mouse. Multiple selection is enabled when the C**trl** or S**hift** keys are pressed at the same time. The selected objects can be deleted by the D**el** key. Or, right-click the  *Show Annotations (A)* button to display a context menu containing these commands:

**Clear All Objects, Clear Annotation Objects, Clear Measurement Objects** Deletes objects of the selected type from the layer.

**Select All Objects, Select Annotation Objects, Select Measurement Objects** Selects objects of the selected type in the layer.


**Visibility of objects** If an ND document contains a time dimension, you can set visibility of any annotation or measurement object throughout the time. Use the secondary mouse click over the object and select *Object Visibility...* The following dialog appears.

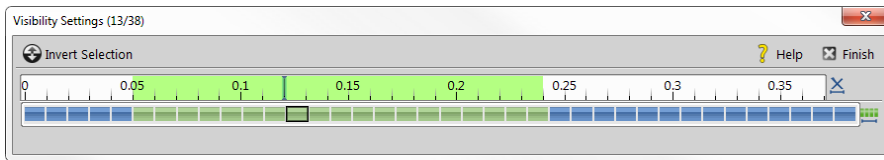Figure 8.14. Visibility Settings






Use **Shift** and **Ctrl** keys and primary mouse click to select frames over which the object will be visible. Once you define the selection, click **Finish**.




## 8.2.3 Measurement Tools


Run the *View > Analysis Controls > Annotations and Measurements*  command to display the following measurement tools:

### 8.2.3.1 Angle



 **Measure Free Angle** Angle between the two connected lines is measured. Click the primary mouse button and drag to draw the first line and release the button. Then use the secondary mouse button click to confirm the first line and repeat the procedure for the second line.



 **Measure 2 Lines** Angle between the two separated lines is measured. Click the primary mouse button and drag to draw the first line and release the button. Then use the secondary mouse button click to confirm the first line. Click the primary mouse button and drag to draw the second line. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).

 **Measure Via Reference Angle** Angle between two lines (one is a reference line) is measured. First use the  *Define Reference Angle* tool to define the reference line. Then select  *Measure Via Reference Angle*, draw the second line using the primary mouse button and confirm the object by the secondary mouse click.



 **Define Reference Angle** Draw a line in the image (using **Draw**) and confirm it using the secondary mouse button or enter a precise value in degrees into the edit box and click **OK**.



### 8.2.3.2 Area



 **Polygon** Polygon area is measured. Draw a polygon by clicking the primary mouse button inside the image. The secondary mouse button or a double-click will place the last node of the polygon and close the object. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).


 **Rectangle** Rectangle area is measured. Draw a rectangle holding down the primary mouse button. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).




 **Ellipse** Ellipse area is measured. Draw a circle holding down the primary mouse button. Click on the edge of the circle and drag to form an ellipse. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).



 **5 Point Ellipse** 5 point ellipse area is measured. Define five different points into the image using the primary mouse button. The system automatically interpolates these points and finishes the ellipse. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).



 **4 Point Ellipse** 4 point ellipse area is measured. Define four different points into the image using the primary mouse button. The system automatically interpolates these points and finishes the ellipse. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).



 **Auto Detect** Area of the detected object is measured. The system automatically detects a homogeneous area around the clicked point. Tolerance (size of the area) can be adjusted by rolling the mouse wheel or by using the Page Up/Page Down keys. Confirm the object by the secondary mouse click.

 **Select Binary** Area of objects detected on the binary layer is measured. The system automatically detects a homogeneous area on the binary layer around the clicked point.

### 8.2.3.3 Length

 **2 Points** Distance between two points is measured. Use the primary mouse button to define the starting and ending point. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).


 **Simple Line** Line length is measured. Click and hold the primary mouse button over the starting position and drag to the ending position and release the button. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).

 **Horizontal, Vertical** Horizontal/vertical distance between two lines is measured. Click the primary mouse button over the starting point and move to the ending point and click again. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).


#### Note



*You can modify default length of the measurement lines. Click the measurement line and resize it by mouse. The new length will be remembered and used for the next measurement.*

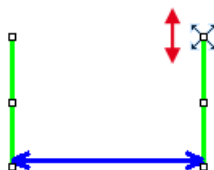

Figure 8.15. Resizing of the measurement line







Alternatively, you can hold the mouse button when placing the first measurement line and draw the line long as you like. The default length will not be affected by this action.



 **Parallel Lines** Distance between two parallel lines is measured. Click the primary mouse button over the first point of the line, move to the second point of the line, adjust the position of both points and confirm it using the secondary mouse click. Then move to the position of the second line and click the primary mouse button, adjust its position and confirm the whole measurement by clicking the secondary mouse button.


 **Polyline** Polyline length is measured. A polyline consists of several line segments. Use the primary mouse button to draw the line segments. Each mouse-click creates a new node. Complete the total length definition by the clicking the secondary mouse button. Confirm the object again by the secondary mouse click if  *Confirm with R-Click* is turned on.

 **Crosses** Distance between two crosses is measured. Click the primary mouse button to define two points between which the distance is measured. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).

 **Auto Default** Distance between two points estimated by the computer is measured. Click and hold the primary mouse button, drag and release the button to draw the line. The system automatically detects the sharpest changes of the pixel intensities and adjusts the point position. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).

 **Auto Outer** Distance between two points estimated by the computer is measured. Click and hold the primary mouse button, drag and release the button to draw the line. The system detects several sharp changes of the pixel intensities and selects the first and the last one as the measurement points. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).

 **Auto Inner** Distance between two points estimated by the computer is measured. Click and hold the primary mouse button, drag and release the button to draw the line. The system detects several sharp changes of pixel intensities and selects the second and the next to the last one as the measurement points. Confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).



 **2 points on XY stage** Measures the distance between two points on a live image which do not fit the camera FOV. Click the primary mouse button to mark the first point, move the stage to the second point on the sample and click in the image to mark the second point.


### 8.2.3.4 3D

 **Length 3D** (requires: EDF Module (Extended Depth of Focus))







This tool enables to perform interactive 3D measurement on ND2 files which include the Z dimension.

#### **Measurement on Slices View**






1. Open an ND2 file containing a Z dimension.
2. Run the `View > Image > ND View > Slices View`  command to display the image in *Slices View*.
3. Select the  *Length 3D* tool. The cursor changes.

4. Select a Z plane (use the mouse wheel) on which the first point lies. Click the primary mouse button to mark that point (crosshair is shown). Confirm its position by the secondary mouse click. The color of the crosshair changes.
5. Select a Z plane on which the second point of your measured line lies. Place the second crosshair as in the previous step and confirm it using the secondary mouse click.
6. Measured value is written into the results table of the *View > Analysis Controls > Annotations and Measurements*  window.





### **Measurement on a Focused Image, Surface View or Volume View**



1. Open an ND2 file containing a Z dimension and run the *Applications > EDF > Create Focused Image*  command.
2. You may switch to  *Show EDF 3D Surface View*,  *Volume View* or stay in the  *Show EDF Focused Image* view which appears automatically. The 3D Measurement behaves the same way in all of the views.
3. Invoke the  *Length 3D* measurement.
4. Define both points of the measured line by the primary mouse click. Measurement is performed automatically and the length is displayed over the line and in the results table of the *View > Analysis Controls > Annotations and Measurements*  window.

#### 8.2.3.5 Radius


-  **Circle** Radius of the circle is measured. Click and hold the primary mouse button and drag to create a circle and release the button. Adjust the circle size and position and confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).
-  **3 Pts. Circle** Radius of the circle is measured. Define three different points into the image using the primary mouse button. Adjust the circle size and position and confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).
-  **N Pts. Circle** Radius of the circle is measured. Define as many points as you like into the image using the primary mouse button and confirm the object by the secondary mouse click. The circle is created by interpolating the defined points. Adjust the circle size and position if needed.


#### 8.2.3.6 Semiaxes


-  **Ellipse** Ellipse semiaxes are measured. Click and hold the primary mouse button and drag to create a circle and release the button. Click on the edge of the circle and drag to form an ellipse. Adjust the ellipse size and position if needed and confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).
-  **5 Pts. Ellipse** 5 point ellipse semiaxes are measured. Define five different points into the image using the primary mouse button. The system automatically interpolates these points and finishes the ellipse. Adjust the ellipse size and position if needed and confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).

 **Free Rectangle** Rectangle semiaxes are measured. Click and hold the primary mouse button and drag to create a rectangle. Adjust its size, position or rotation (using the green anchor point) and confirm the object by the secondary mouse click (only if  *Confirm with R-Click* is turned on).

### 8.2.3.7 Count and Taxonomy

 **Count** Click inside the image to insert counted points. The current count is shown next to each point and its position and number is recorded. Finish counting by the secondary mouse click.

 **Taxonomy** This tool enables to manually count objects of up to 70 classes.

1. Click the  icon.
2. Use the secondary mouse click over the results table and select *Taxonomy Options*.
3. In the *Taxonomy Options* window, define the number of classes (from 2 to 70) and the way of inserting markers:
  - On left button down** Inserts a marker to the current cursor position after a mouse click. If this option is selected, the function keys enable to switch between the classes. In this mode, the **Fn** keys works for switching between classes.
  - On Fn key at cursor position** After the **Fn** key is pressed, the count is performed and the corresponding marker is placed to the current cursor position.
  - On Fn key at screen center** After the **Fn** key is pressed, the count is performed and the corresponding marker is placed to the center of the image view.
  - On Fn key at cursor position, no marker inserted** After the **Fn** key is pressed, the count is performed and no marker is placed.
4. Start the count according to the *Taxonomy Options* setting.

#### Note

The keyboard Fn (function) keys F1 - F10 can be used as a shortcut for classification/class selection. Since there are 70 classes available, certain key combinations must be pressed in order to select/classify the right class:

<i>Key Combination</i>	<i>Class Number</i>
<i>Fn</i>	<i>n</i>
<i>Ctrl + Fn</i>	<i>10 + n</i>
<i>Alt + Fn</i>	<i>20 + n</i>
<i>Shift + Fn</i>	<i>30 + n</i>
<i>Ctrl + Alt + Fn</i>	<i>40 + n</i>
<i>Ctrl + Shift + Fn</i>	<i>50 + n</i>
<i>Shift + Alt + Fn</i>	<i>60 + n</i>

5. Perform the measurement until all classes are counted.

## 8.3 Automated Measurement

(requires: MEAS (Automatic Measurement))

Automated measurement is the most powerful of image analysis features. In combination with user macros, *NIS-Elements* can become a semi- or fully-automatic image analysis tool. There are several key procedures which a proper automated measurement should involve:

1. Optical system calibration (see the 4.1.4 Optical Configurations chapter)
2. Image acquisition (see the 5. Image Acquisition chapter)
3. Defining threshold, creating the binary layer (see the 7.2.1 Thresholding chapter)
4. Performing the automated measurement
5. Results presentation

### **Image Layers Involved in Measurement**

The binary layer and also the color image can get involved in the automated measurement:

- *Binary layer* - is typically used for shape and size measurements (area, perimeter, etc).
- *Color layer* - intensity or hue measurements are carried out on this layer. The area covered with the binary layer is used as the source data.

### **Object/Field Measurement**

*NIS-Elements* distinguishes two types of automated measurements: object and field.

**Object Measurement** Groups of neighboring pixels of the binary layer are called objects. Results of the object measurement are usually the object data such as Length, Area, CentreX/Y (the X,Y position of the central pixel of an object), etc.

**Field Measurement** Field measurements produce information regarding the whole measurement area (measurement frame, ROI). Such information can be Area Fraction, Mean Brightness, Density Variation, etc.

Please see the complete list of measurement features in the 8.4 Measurement Features chapter.

### **Restricting the Area of Measurement**


The area of measurement can be restricted by the measurement frame or by a user defined region of interest (ROI).

**Region of Interest - ROI** ROI is a user defined area of the image. Whenever the ROI is displayed (by the *View > Layers > Measurement ROI* command), it is also applied to restrict the area of measurement.

**Measurement Frame** The measurement frame is a re-sizable rectangular frame which serves for bounding the area accepted for measurement. Turn the measurement frame ON via the *Measure > Use Measurement Frame* command.

Binary objects touching edges of ROI or the measurement frame can be treated differently (excluded from the measurement or included). This behavior can be specified within the *Measure > Options* window.

### **Statistics and Data Presentation**

*NIS-Elements* calculates basic statistics - mean value, standard deviation and distribution of all measured features. The features to be measured can be selected in the *Measure > Object Features* and the *Measure > Field and ROI Features* windows. Results of all automated measurements are presented and can be exported from the *View > Analysis Controls > Automated Measurement Results*  control panel.

## 8.4 Measurement Features

(requires: MEAS (Automatic Measurement))

The following list contains all features which can be measured within *NIS-Elements* Measurements. There are the following. Different measurement tools measure different features. If not stated otherwise, all distances and areas are measured in the current (image) calibration units. If the image is not calibrated, pixels are used instead.

**MAN** Features available in manual measurement.

**OBJ** Object features - measured under binary objects (layers).

**FLD** Field features - measured over the whole image area.

**ROI** ROI features - measured within ROIs.

**EDF** EDF features - can be measured only after EDF focused Image was created.

**AcqTime** **OBJ**, **ROI**, **FLD**

Acquisition time of the current image in seconds. If the image is a part of an image sequence, the time is measured from the beginning of the sequence. Otherwise it is the time elapsed since the software was started.


**Angle** **MAN**

Angle between two user-defined lines. See 8.2.3 Measurement Tools.

**Area** **OBJ**, **MAN**

Area is a basic quantity for measuring object size.

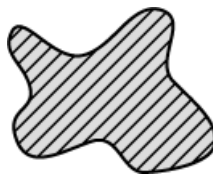Figure 8.16. Area



**Axis A, Axis B, Axis A/B** **MAN**

Lengths of the major (A) and minor (B) axes of an ellipse and their ratio. See 8.2.3 Measurement Tools.

**BinaryArea** **FLD**

Sum of areas (*Area*) of all binary objects.

**BinaryAreaFraction** **FLD**

$$\text{BinaryAreaFraction} = \frac{\text{BinaryArea}}{\text{MeasuredArea}}$$

See *BinaryArea* and *MeasuredArea*.


### Blue **MAN**

Arithmetic mean of pixel intensities of the blue component within the measured area. See 8.2.3 Measurement Tools. It equals 0 on other than RGB images.

### **BoundsAbsLeft, BoundsAbsRight, BoundsAbsTop, BoundsAbsBottom** **OBJ**

These are absolute distances (X/Y coordinates) of the left, right, top, and bottom edges of the object bounding rectangle within a motorized stage range. This feature can be measured only if a motorized stage is connected.

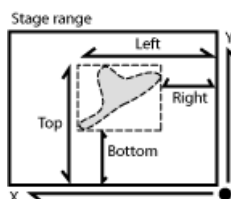
Figure 8.18. *BoundsAbs*



### **BoundsLeft, BoundsRight, BoundsTop, BoundsBottom** **OBJ**

These are distances (X/Y coordinates) of the left, right, top, and bottom edges of the object bounding rectangle.

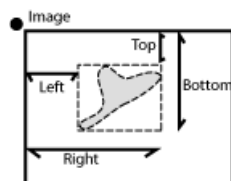Figure 8.19. *Bounds*



### **BoundsPxLeft, BoundsPxRight, BoundsPxTop, BoundsPxBottom** **OBJ**

These are distances (X/Y coordinates) of the left, right, top, and bottom edges of the object bounding rectangle measured in pixels.

### **BrightVariation** **OBJ**, **ROI**, **FLD**

It is the standard deviation of brightness values within the object.

### **CentreX, CentreY** **OBJ**, **ROI**, **FLD**, **MAN**

Coordinates of the center of gravity of the object (ROI, field).

### **CentreXabs, CentreYabs** **OBJ**, **ROI**, **FLD**

Absolute coordinates of the center of gravity in the scope of the stage XY range. These features can be measured only when a motorized stage is connected. See also *BoundsAbsLeft*.

### **CentreXpx, CentreYpx** **OBJ**, **ROI**, **FLD**

---

Coordinates of the center of gravity measured in pixels.

**Channel** **OBJ** , **ROI** , **FLD** , **MAN**

Comma-separated list of channel names on which the measurement was performed.

**Circularity** **OBJ**

Circularity of a circle equals 1. All other shapes are characterized by circularity smaller than 1. It is a derived shape measure, calculated from *Area* and *Perimeter*. It is useful for examining shape characteristics.

$$Circularity = \frac{4 \times \pi \times Area}{Perimeter^2}$$

**Comment** **OBJ** , **ROI** , **FLD** , **MAN**

This feature adds an editable column to the results table.

**Convexity** **OBJ**

Indicates convexity of the object edges.

$$Convexity = \frac{Area}{Convex\ Hull\ Area}$$

*Convex Hull Area* is the area of objects after running the *Binary > Convex Hull* command.

**DensityVariation** **OBJ** , **ROI** , **FLD**

It is the usual standard deviation of the density values.

**Diameter** **MAN**

Diameter of a circle measured manually. See 8.2.3 Measurement Tools.

**EdfRoughness** **OBJ** , **ROI** , **EDF** , **FLD** , **MAN**

Based on the 3D EDF surface model, it indicates how much the surface is rough. Its value for a flat surface is 1.

$$Edf\ Roughness = \frac{Edf\ Surface}{Area}$$

See also *EdfSurface* and *Area*.

**EdfRoughnessRA** **OBJ** , **ROI** , **EDF** , **FLD** , **MAN**

Based on the 3D EDF surface model, it indicates how much the surface is rough. It is a sum of absolute values of differences between Z value of each pixel and mean of all Z values divided by *Area*. Its value for a flat surface is 0. It is the most widely used one-dimensional roughness parameter



$$EdfRoughnessRA = \frac{1}{n} \sum_{i \in \Omega} |Z_i - EZ|$$

$$Area = |\Omega| = n \quad EZ = \frac{1}{n} \sum_{i \in \Omega} Z_i$$

**EZ** The mean height of the Z-map (mean of all Z coordinates of the EDFSurface).


**n, Area,  $\Omega$**  Number of pixels in the measured area (image/object/ROI).

**$Z_i$**  EDFSurface Z coordinate of a particular pixel.

**EdfSurface** **OBJ** , **ROI** , **EDF** , **FLD** , **MAN**

Based on the 3D surface model, it is an approximation of the surface area. For a flat surface, *EdfSurface* = *Area*.

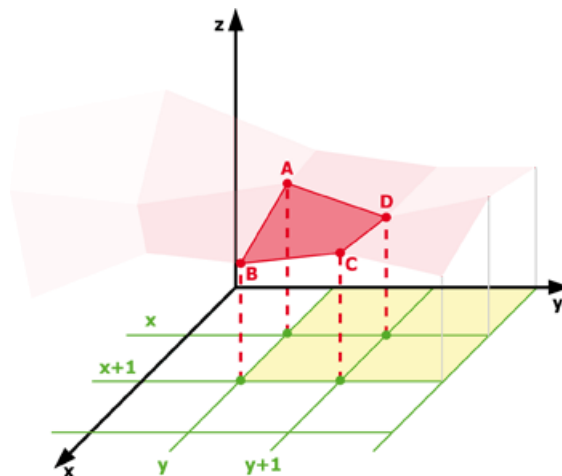Figure 8.24. EdfSurface highlighted red, Area of four Edf surfaces highlighted yellow



$$A(x, y, z1)$$

$$B(x + 1, y, z2)$$

$$C(x + 1, y + 1, z3)$$

$$D(x, y + 1, z4)$$

$$a = |AB|$$

$$b = |BC|$$

$$c = |CD|$$

$$d = |DA|$$

---

$$EdfSurface = \frac{a \times b}{2} + \frac{c \times d}{2}$$

$$EdfSurface \geq Area$$

**Elongation** **OBJ**

Characterizes the object shape. Ratio of *MaxFerret* and *MinFerret* features.


$$Elongation = \frac{MaxFerret}{MinFerret}$$

**EqDiameter** **OBJ** , **MAN**

A size feature derived from the area. It determines the diameter of a circle with the same area as the measured object:

$$EqDiameter_{2D} = \sqrt{\frac{4 \times Area}{\pi}}$$

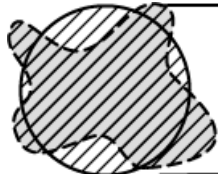Figure 8.31. EqDiameter


**FieldID** **OBJ** , **FLD** , **ROI**

Automatically generated field ID. Each measurement per field generates a new ID.

**FillArea** **OBJ**

If an object contains holes, *FillArea* remains the same while *Area* is reduced by the area of the holes. If an object does not contain holes, then *FillArea* = *Area*.

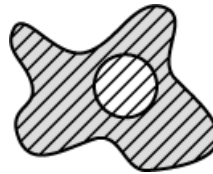Figure 8.32. FillArea

**FillRatio** **OBJ**

*FillRatio* is the ratio of *Area* and *FillArea*:

$$FillRatio = \frac{Area}{FillArea}$$

If an object does not contain holes, *FillRatio* = 1. If an object contains holes, *FillRatio* is less than 1. This feature helps distinguish objects with and without holes.

**Green** **MAN**

Arithmetic mean of pixel intensities of the Green component within the measured area. See 8.2.3 Measurement Tools. It equals 0 on non-RGB images.

**Height** **MAN**

Projection of 3D length measurement to the Z axis. See 8.2.3 Measurement Tools.

**HueTypical** **OBJ** , **ROI** , **FLD**

HueTypical (H) is the hue value with maximum frequency in the hue-values histogram. This feature describes the most frequent hue (color) in an object or field.

$$\alpha = \frac{1}{2} (2R - G - B)$$

$$\beta = \frac{\sqrt{3}}{2} (G - B)$$

$$H = atan2(\beta, \alpha)$$

**HueVariation** **OBJ** , **ROI** , **FLD**

HueVariation is the usual standard deviation of hue values. This feature describes hue (color) distribution of inner structure of an object or a field.

$$\alpha = \frac{1}{2} (2R - G - B)$$

$$\beta = \frac{\sqrt{3}}{2} (G - B)$$

$$H = \text{atan2}(\beta, \alpha)$$

**IntensityVariation** **OBJ**, **ROI**, **FLD**


IntensityVariation is derived from the intensity histogram. It is an usual standard deviation of intensity values. This feature describes the inner structure of an object or a field.

**Length** **OBJ**, **MAN**

Length is a derived feature appropriate for elongated or thin structures. Since it is based on the rod model, it is useful for calculating length of medial axis of thin rods.

$$\text{Length} = \frac{\text{Perimeter} + \sqrt{(\text{Perimeter}^2 - 16 \times \text{Area})}}{4}$$

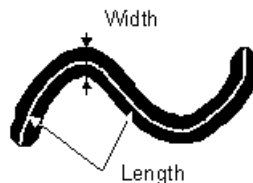Figure 8.37. Length



**Length xyz, Length xy** **MAN**

3D length measured manually. *Length xyz* is the actual length measured on *Slices View* or *Volume View*. *Length xy* is length of a projection of the measured line to the XY plane. See 8.2.3 Measurement Tools.

**LineLength** **OBJ**


It is the length of an object with elongated shape.

**MaxFerret, MaxFerret90** **OBJ**, **MAN**

It is the maximal value of the set of Feret's diameters. Generally (for convex objects), Feret's diameter at angle  $\alpha$  equals the projected length of the object at angle  $\alpha$ ,  $\alpha (0,180)$ ; *NIS-Elements* calculates Feret's diameters for  $\alpha = 0, 10, 20, 30, \dots, 180$ .

The MaxFerret90 is a length projected across the MaxFerret diameter.

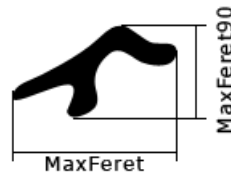Figure 8.38. MaxFeret



**MaxIntensity** **OBJ** , **ROI** , **MAN** , **FLD**

It is derived from the intensity histogram as the maximal of all pixel intensity values (I).

Figure 8.39. Intensity

$$I = \frac{1}{3}(R + G + B)$$

**Mean\*ChannelName\***, **MeanRed**, **MeanGreen**, **MeanBlue** **OBJ** , **ROI** , **FLD**

Arithmetic mean of pixel intensities of one image component. See also *MeanIntensity*.

**MeanBrightness** **OBJ** , **ROI** , **MAN** , **FLD**

Arithmetic mean of brightness values of pixels.

**MeanChord** **OBJ** , **ROI** , **FLD**

The mean value of secants in the 0, 45, 90 and 135 degrees directions. It is a derived feature and is calculated from the *Area* and mean projection according to the following formula.

$$MeanChord = \frac{4 \times Area}{(Pr_0 + Pr_{45} + Pr_{90} + Pr_{135})}$$

**MeanDensity** **OBJ** , **ROI** , **MAN** , **FLD**

Arithmetic mean of density (O.D.) values of pixels.

$$O.D. = -\log \frac{PixelIntensityValue}{MaxIntensityValue}$$

**MeanIntensity** **OBJ** , **ROI** , **MAN** , **FLD**

It is derived p from the intensity histogram as the arithmetic mean of pixel intensities (I).

Figure 8.42. Intensity

$$I = \frac{1}{3}(R + G + B)$$

$$I = \frac{1}{3}(R + G + B)$$

**MeanRatio, MeanCa2+, MeanCorrFRET, MeanFRETeff, MeanTitration** **OBJ**, **ROI**, **FLD**

Arithmetic mean of intensities of the *Ratio*, *Ca2+*, *FRET / Corr FRET / FRET Eff / Ph* channels. This feature is activated upon creation of the channel view.

**MeanSaturation** **OBJ**, **ROI**, **FLD**

Arithmetic mean of saturation values of pixels (S).

Figure 8.43. Saturation

$$S_{HSI} = \begin{cases} 0, & \text{if } C = 0 \\ 1 - \frac{m}{I}, & \text{otherwise} \end{cases}$$

$$M = \max(R, G, B)$$

$$m = \min(R, G, B)$$

$$C = M - m$$

**MeasuredArea** **OBJ**, **ROI**, **FLD**


The area of the measurement frame or a ROI within the measurement frame (if the ROI is on).

**MinDistanceTo** **OBJ**

**MinFerret** **OBJ**, **MAN**

The minimal value of the set of Feret's diameters. Generally (for convex objects), Feret's diameter at angle  $\alpha$  equals the projected length of object at angle  $\alpha$ ,  $\alpha$  (0,180); *NIS-Elements* calculates Feret's diameter for  $\alpha = 0, 10, 20, 30, \dots, 180$ .

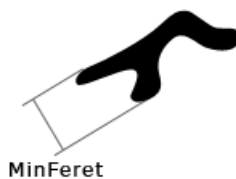Figure 8.45. MinFerret



**MinIntensity** **OBJ** , **ROI** , **MAN** , **FLD**

Its value is derived from the intensity histogram. It is the minimum of intensity values of pixels.

Figure 8.46. Calculation of Intensity ( $I$ )

$$I = \frac{1}{3}(R + G + B)$$

**ND.M** **FLD** , **ROI** , **MAN**

Frame index within the Multipoint dimension.

**ND.T** **FLD** , **ROI** , **MAN**

Frame index within the time dimension.

**NearestObjDist** **OBJ**

The smallest distance to another object (measured between centers of gravity).

Note

If ROIs are used during the measurement, NearestObjDist is calculated only between the objects within the specific ROI (objects having the same ROI ID).

**NumberObjects** **FLD** , **ROI**

The number of objects inside the measured area. Exclusion rules for counting objects are taken into account (see 8.1.4 Measurement Options).

**NumberObjectsRestricted** **FLD**

Number of objects left after applying restrictions.

See also *View > Analysis Controls > Restrictions* .

**Numbers** **MAN**

Number of points clicked within one *Count* measurement. See 8.2.3.7 Count and Taxonomy.

**Obj ID** **OBJ**

Binary object ID. Each binary object has a unique ID

**ObjectAreaFraction** **ROI**

Ratio of *BinaryArea* and *ROIArea*.

$$\text{ObjectAreaFraction} = \frac{\text{BinaryArea}}{\text{ROIArea}}$$


**Objects Area** **ROI**

Area of all binary objects placed within *ROIArea*.

**Orientation** **OBJ**

The angle at which Feret's diameters have their maximum. The diameters are calculated with 5 degrees angle increment.

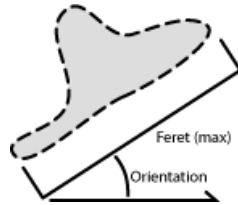Figure 8.47. Orientation



**OuterPerimeter** **OBJ**

Length of the outer perimeter (unlike in *Perimeter*, holes inside the object do not count).

**Perimeter** **OBJ**, **ROI**, **MAN**, **FLD**

Perimeter is the total boundary length. It includes both the outer and inner boundaries (if there are holes inside an object). The perimeter is calculated from four projections in the directions 0, 45, 90 and 135 degrees using Crofton's formula

$$Perimeter = \frac{\pi \times (Pr_0 + Pr_{45} + Pr_{90} + Pr_{135})}{4}$$

Figure 8.48. Perimeter



**PerimeterContour** **OBJ**

Total boundary length computed from object contours. Compared to *Perimeter*, it is more accurate (especially on elongated objects) but slower to get the result.

**Radius** **MAN**

Radius of a circle measured manually. See 8.2.3 Measurement Tools.

**Red** **MAN**

Arithmetic mean of pixel intensities of the Red component within the measured area. See 8.2.3 Measurement Tools. It equals 0 on non-RGB images.

**Roi ID** **OBJ**, **ROI**

ROI ID. Each ROI has a unique ID.

**ROIArea** **ROI**



Area of each ROI.

### **Roughness** **OBJ**, **MAN**

This feature indicates how much the object is rough. “1” means the object roughness is minimal (it is circular). *Roughness* values range is <0;1>.

$$Roughness = \frac{Convex\ hull\ perimeter}{Perimeter}$$

### **RoughnessInf** **OBJ**

This feature indicates how much the object is rough. “1” means the object roughness is minimal (the object is circular). *RoughnessInf* values range is <1;inf>.


$$RoughnessInf = \frac{1}{Roughness}$$

### **ShapeFactor** **OBJ**, **MAN**

This feature is to define whether the object is rough or not.

$$ShapeFactor = \frac{4 \times \pi \times Area}{Convex\ hull\ perimeter^2}$$

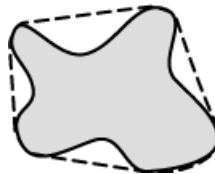Figure 8.52. ShapeFactor




### **Source** **OBJ**, **FLD**, **ROI**, **MAN**

Name of the measured image file.

### **StartX, StartY** **OBJ**

When you scan the image from the origin in left-to-right direction, the first pixel of the object you hit is the one with *StartX* and *StartY* coordinates.

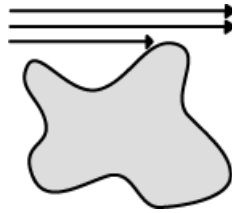Figure 8.53. StartX, StartY




**StartXpx, StartYpx** **OBJ**

StartX and StartY coordinates measured in pixels (see StartX).

**StgPosX, StgPosY** **FLD** , **ROI**

Coordinates of the absolute position of the measured field. It is available only for systems equipped with a motorized XY stage.

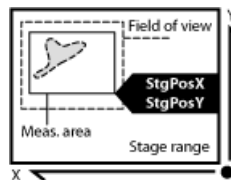Figure 8.54. StgPos



**Sum\*ChannelName\*** **OBJ** , **ROI** , **FLD**

The sum of intensity of all pixels of the object, ROI or field.

**SumBrightness** **OBJ** , **ROI** , **MAN**

A sum of brightness in every pixel of the object.

**SumDensity** **OBJ** , **ROI** , **MAN** , **FLD**

A sum of individual optical densities (O.D.) of each pixel in the measured area. This feature describes, for instance, the amount of a substance in biological sections. Optical density is evaluated according to the following formula:

$$O.D. = -\log \frac{PixelIntensityValue}{MaxIntensityValue}$$

**SumIntensity** **OBJ** , **ROI** , **FLD** , **MAN**

Sum of intensity in every pixel/voxel of the object.

**SurfVolumeRatio** **ROI** , **FLD**

SurfVolumeRatio is a feature with a strong stereological interpretation: if you measure on fields that are sampled systematically and independently of the content of the sections, then the feature is an unbiased estimator of the surface area of objects (inner structure) per volume of the whole sample.

$$\text{SurfVolumeRatio} = \frac{\frac{4}{\pi} \times \text{Perimeter}}{\text{MeasuredArea}}$$


See also *Perimeter* and *MeasuredArea*.

### Time **MAN**

Time interval measured on a 1DT document. See also *Velocity*.

### Velocity **MAN**

Speed measured between two points defined manually. This feature can be measured on a 1DT image using length measurement tools. 1DT image can be created from a timelapse ND file.

1. Open a timelapse ND2 file and run *View > Image > ND View > Slices View* .
2. Right-click one of the side views and select *Create New Document from this View*. A new image will be created.
3. Perform length measurement on the new image.

### VolumeEqCylinder **OBJ**


This parameter is based on the rod model. *Length* is interpreted as height and *Width* is interpreted as the base diameter of a cylinder.

$$\text{VolumeEqCylinder} = \frac{(\pi \times d^2)(l - d)}{4} + \frac{\pi \times d^3}{6}$$

$$l = \max(\text{MaxFeret}, \text{Length})$$

$$d = \min(\text{MinFeret}, \text{Width})$$

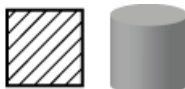Figure 8.58. VolumeEqCylinder




### VolumeEqSphere **OBJ**

Volume of a sphere whose diameter is equal to *EqDiameter* of the measured 2D object.

$$\text{VolumeEqSphere} = \frac{\pi \times \text{EqDiameter}^3}{6}$$

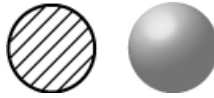Figure 8.59. VolumeEqSphere




**Width OBJ**

Width is a derived feature appropriate for elongated or thin structures. It is based on the rod model and is calculated according to the following formula:


$$Width = \frac{Area}{Length}$$

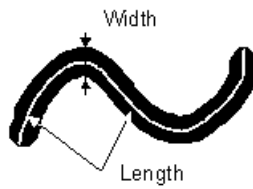
Figure 8.61. Width



## 9. Creating Reports

### 9.1 Report Generator

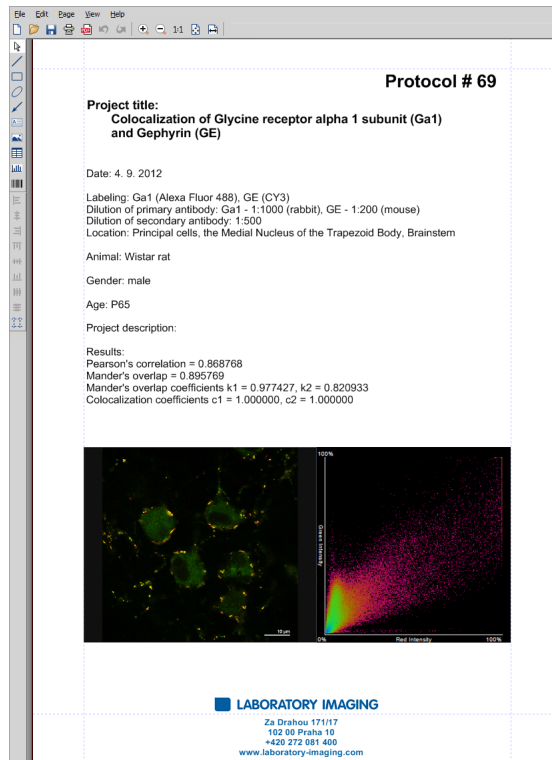Report Generator allows the user to create a print-ready document which can contain measured data, images, graphs, etc. Once you start it by the *File > Report*  *> New Blank Report* command, a graphical editor appears.



Creating of a report in *Report Generator* is very similar to creating any document in a common text editor. You can insert text fields, simple shapes, images, tables or graphs. As well as in text editors, you can create templates containing objects which do not change such as headers and footers, but also dynamic objects to be filled with data (e.g.: image name, author, etc) filled upon creation of the report. This comes especially handy in connection with the *Database* module where database fields are loaded to the report(s) automatically.

### 9.2 Report Objects

A report is composed from report objects. In the left toolbar, there are tools for inserting different objects.

1. Select one of the tools

2. Click and move the mouse inside the report to determine the area reserved for the object. Release the mouse button.
3. Double-click the report object to open a window where its properties can be adjusted.
4. Confirm the changes by **OK**.

Tip


Vector objects (report objects, annotations, interactive measurement objects, ROIs) can be copied by “drag and drop” while holding the **Shift** key down.

1. Select the object(s).
2. Press **Shift** and drag the object somewhere.
3. A copy of the object is placed to where you release the mouse button.

## 9.2.1 Aligning Objects


### **Align Objects**

You can align two objects to the same horizontal or vertical level.

1. Select more objects (e.g. by holding down **CTRL** and clicking with the primary mouse button).
2. Right-click one of the selected objects and select the *Align or Distribute > Align* command from the context menu.
3. The objects will be aligned as indicated on the command icon.
4. If the *Relative to Page* (  ) option is turned on, the objects will be aligned to the edges or the center of the page.

### **Distribute Objects**

Objects of similar size can be distributed uniformly in the horizontal/vertical direction.

1. Select three or more objects.
2. Right click one of the objects and select the *Align or Distribute > Distribute Horizontally* command from the context menu.
3. Distances between the objects will be adjusted to equal.
4. If the *Relative to Page* option (  ) is turned ON, the outer objects will be moved to the edges of the page.

### **Match Object Sizes**

Sizes of objects can be unified too.

1. Select two or more objects.
2. Right-click the “master” object to the size of which you want to resize the other objects.
3. Select one of the *Resize Objects* sub-menu commands.

## 9.2.2 Dynamic Data

Results of automatic-interactive measurement, graphs, or the current image can be inserted to reports.

1. Insert an object which can contain dynamic data (text, picture, table, or graph).
2. Right-click the object and select *Insert Dynamic Data* or *Insert Dynamic Picture* from the context menu.
3. A window appears.
4. Select one of the available sources, and click *Next*.
5. Finish the source definition and click . The data appear on the report page.

### **List of Available Dynamic Data Sources:**



- *Data inserted by user (available for: image, text, table)* - the system will ask you to type a text or browse for an image to insert during the report creation. When creating a dynamic object of this type, a query text which will be used to prompt you for the data can be defined.
- *System data (available for: text, table)* - enables you to insert some general data such as date, the name of the user account currently logged in, the page number, or the page count.
- *Macro (available for: text, table)* - enables you to insert expressions, values, or results of a macro.
- *Measurement (available for: text, table, graph)* - enables you to insert results of automatic/interactive measurement.
- *Database (available for: image, text, table)* - this data source is displayed only when creating a database report template (see below). It enables you to insert a link to any of the database records.

The real purpose of the dynamic data turns up when creating reports using report templates.

## 9.3 Report Templates


A report template is a layout defining the appearance of future reports, which is ready to be filled with data. The data can be inserted by the user “on demand” or automatically (dynamic data).

### **To Create a Report Template**

1. Run the *File > Report*  *> New Blank Report* command. An empty report appears.
2. In the Report Generator, select *File > Change to Template*.
3. Edit the report template in the same way as a common report.
4. Insert the dynamic data where it is appropriate - measurement results to text boxes or tables, the current image to the image field, etc.).
5. Save the report template (\*.rtt) via the *File > Save*  command.

---

### **To Create a Report FROM Template**

1. If the template is opened inside the Report Generator, run the *File > Create Report* command. Otherwise, use the *File > Report  > New Report from Template > Browse* command inside the main application window to open a template from hard disk.
2. A report opens and the dynamic data is inserted automatically.
3. Save the report, print it, or export the page(s) to PDF using the commands from the *File* menu.




# 10. Macros

## 10.1 Creating Macros

A macro - an executable sequence of commands - can make the work very effective. *NIS-Elements* provides a C-like programming language utilizing its internal set of functions. The sequence of functions can be created either by recording the performed actions, by writing the functions within the macro editor, or by modifying the command history (the history is recorded automatically during the work). The macro can be saved to an external (\*.mac) file for later use.

### **Recording a Macro**

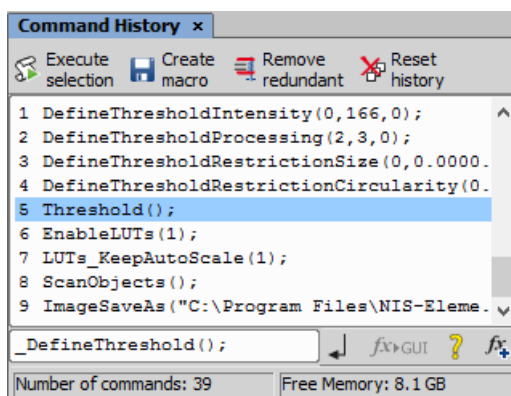
The fastest way to create a macro is to record it.

- Start the *NIS-Elements* macro recorder by selecting the *Macro > Record* command.
- Perform the series of actions you would like to record.
- Finish the recording by the same command again (its name changes to “Stop Recording”)
- It is recommended to check the macro in the macro editor before saving. Run the *Macro > Edit*  command to display it.
- Save the macro to a file via the *Macro > Save As* command.

### **Creating a Macro from History**


You can create a macro using the list of recently performed commands.

- 1) Run the *Macro > Command History* command to display the *Command History* control panel:



- 2) Push the *Create Macro* button.

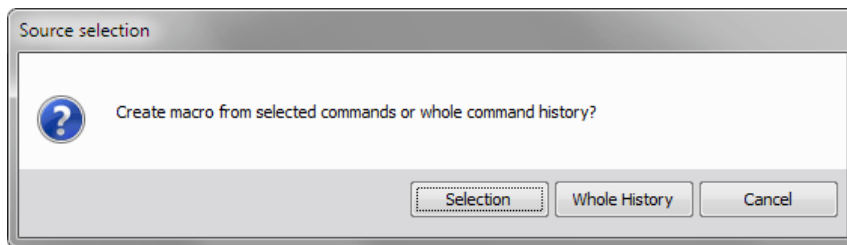- 3) Select what portion of the command history will be used in the macro:



**Selection** This option will use only commands which you have previously selected by mouse within the *Command History* window. Group selection (holding **Shift** or **Ctrl**) is available.


**Whole History** All commands listed in the *Command History* window will be used for the new macro.


- 4) After you selected one of the options, the macro editor appears containing the new macro source code. Use the editor to fine-tune it.

Note


*The Remove Redundant command is automatically performed before the macro is opened in the editor.*




### **Writing/Editing a Macro**

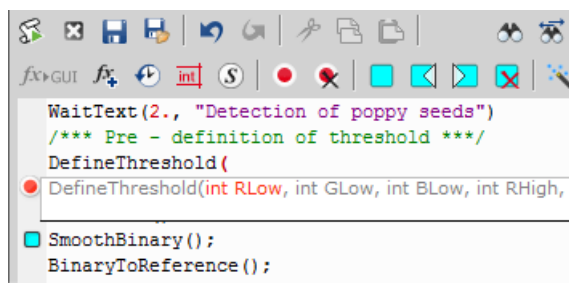If you possess programmers skills, a macro can be written by hand in the built-in macro editor. Run the *Macro > Edit*  command to display it.






The macro editor provides:

- Undo/Redo functionality.
- Command insertion from the list of available commands.
- Interactive command names list. Press **Ctrl+Space** to display a simplified list of all commands while typing the macro.
- Syntax hints (parameter types and names) appear as you type a command name.
- Bookmarks  can be placed to the code so you can easily roll to the important parts of the macro.

- Breakpoints  can be placed to the code. A breakpoint forces the macro to stop the execution at a certain point so you can check the state of variables and therefore inspect the macro functionality thoroughly.
- Syntax highlighting.
- Help on commands with detailed description.
- Use the *Save Macro*  and the *Save Macro As...*  commands in the macro editor to save the created macro. *NIS-Elements* automatically remembers the folder from where the macros were run the last time and will open this folder when user clicks on Open/Save command in Macro menu.

## 10.2 Running a Macro

*NIS-Elements* provides several ways to run a macro.

- You can run the current macro loaded to *NIS-Elements* by choosing the *Macro > Run*  command or by pressing F4.
- A *hotkey* combination (Ctrl+Alt+1,2,3 ... 9) can be assigned to it in the *Macro > Options*  window.
- You can run a macro at the beginning of the *NIS-Elements* session, by assigning the *StartUp* flag to the macro in the *Macro > Options*  window.
- You can run a macro saved to a disk by using the *Macro > Run From* command.
- A toolbar button can be assigned to run a macro. See 3.3 Modifying Toolbars.

### **Breaking the Macro**

The macro execution can whenever be stopped by pressing the Ctrl+Break key shortcut.

## 10.3 Interactive Advanced Macro (API)

(requires: *INTADV (Interactive Advanced Macro - API)*)

If the API module is installed, it dramatically extends the macro programming capabilities. It adds the following functionality to the main application:

**Loops and conditions** Loops and conditions are enabled (if, for, while...)

**User interaction functions** Contains functions, which enable interaction with user like: type value, text, combo boxes, list boxes and many other Windows standard controls with easy and intuitive code handling.

**Support of importing functions from external DLLs** Functions from external DLLs can be imported at the start of the macro and used during the macro execution.

See 10.4 Macro Language Syntax.

---

## 10.4 Macro Language Syntax

Specifies the *NIS-Elements* Macro Language features.

### 10.4.1 Variable types

The following data types are implemented:

**char** text

**char8** <-128, +127>

**byte** <0, +255>

**int** <-32.768, +32.767>

**int64** <-9.223.372.036.854.775.808, +9.223.372.036.854.775.807>

**word** <0, +65.535>

**long** <-2.147.483.648, +2.147.483.647>

**dword** <0, +4.294.967.295>

**double** <1.7E +/- 308 (15 digits)>

### 10.4.2 Structures and Unions

Structures and unions are not supported.

### 10.4.3 Arrays

One and two dimensional arrays are supported.

### 10.4.4 Local and Global Variables

You should declare local variables at the beginning of macro or function only. You should declare global variables only at the beginning of macro. You can run two nested macros declaring the same global variables, but they must be of the same type. You can declare global variables by prefixing the declaration *global*. Eg.:

```
global int Number_Rows;  
global char buffer[200];
```

The “global” keyword in front of the variable definition assigns the variable to the global scope. Such variable is then accessible from all function scopes within the macro interpreter.

## 10.4.5 Statements

### **Supported Statements**

**for** (requires: *INTADV (Interactive Advanced Macro - API)*)

The *for* statement lets you repeat a statement a specified number of times.

Syntax

```
for([init-expr]; [cond-expr]; [loop-expr])
    statement
```

First, the initialization (*init-expr*) is evaluated. Then, while the conditional expression (*cond-expr*) evaluates to a non zero value, Statement is executed and the loop expression (*loop-expr*) is evaluated. When *cond-expr* becomes 0, control passes to the statement the following the *for* loop.

**while** (requires: *INTADV (Interactive Advanced Macro - API)*)

The *while* statement lets you repeat a statement until a specified expression becomes false.

Syntax

```
while(expression) statement
```

First, the expression is evaluated. If expression is initially false, the body of the *while* statement is never executed, and control passes from the *while* statement to the next statement in the program. If expression is true (nonzero), the body of the statement is executed and the process is repeated.

**if, else** (requires: *INTADV (Interactive Advanced Macro - API)*)

Conditionally executes a statement or group of statements, depending on the value of an expression.

Syntax

```
if(expression) statement1
    [else statement2]
```

The *if* keyword executes *statement1* if expression is true (nonzero); if *else* is present and expression is false (zero), it executes *statement2*. After executing *statement1* or *statement2*, control passes to the next statement.

**goto** Transfers control of the program execution.

Syntax

```
goto name;
    . . .
    name: statement
```

You cannot use *goto* to jump inside the block from outside. E.g. following is not allowed:

```
goto label;
    if(k>5)
    {
    label:
        DilateBinary(3, 5);
```

```
FillHoles();
}
```

You cannot use `goto` to jump out from the block more than 2 block levels down. E.g. the following is not allowed:

```
for(i=0; i<64; i=i+1)
{
  for(j=0; j<64; j=j+1)
  {
    if(a[i] > b[i])
    {
      value = i;
      goto end; // crossing 3 right brackets
    }
  }
}
end
```

**break** Terminates the execution of the nearest enclosing statement.

Syntax

```
break;
```

The *break* keyword terminates the execution of the smallest enclosing *for* or *while* statement in which it appears. Control passes to the statement that follows the terminated statement.

**continue** Passes control to the next iteration of the statement in which it appears.

Syntax

```
continue;
```

The *continue* keyword passes control to the next iteration of the *for* or *while* statement in which it appears. Within a *while* statement, the next iteration starts by reevaluating the expression of the *while* statement. Within a *for* statement, the first expression of the *for* statement is evaluated. Then the compiler reevaluates the conditional expression and, depending on the result, either terminates or iterates the statement body.

### **Statements which are not supported**

**do, switch, case, default, typedef** These statements are not supported.

## 10.4.6 Directives

The following directives are supported by the system:

**define** The *#define* directive assigns a meaningful name to a constant in a program.

```
//Syntax
```

```
#define identifier token-string
```

The directive substitutes token-string for all subsequent occurrences of the identifier in the source file. Token-string can be a value or a string (only for 32-bit version of *NIS-Elements*).

```
//Example
```

```
#define ERROR_SPRINTF 0
#define MAINDIR "c:\Images"

int main()
{
char buf[256];
int retval;

retval = sprintf(buf, "%s", "MAINDIR");
if(retval == ERROR_SPRINTF)
Beep();
else WaitText(0., buf);
return TRUE;
} \
```

**include** Specifies the name of the file to be included.

```
//Syntax
```

```
#include filename
```

The *#include* directive includes the contents a file with a specified filename in the source program at the point where the directive appears.

Example

// if you do not specify the full path, NIS-Elements assumes, that it is a \ relative path to a main directory

```
#include "macros\my_macro.h"
#include "c:\NIS-Elements\macros\my_macro1.h"
```

**import** (requires: *INTADV* (*Interactive Advanced Macro - API*))

The *#import* directive is used to incorporate information from an external library.

```
//Syntax
```

```
#import("DLLname");
#import function_declaration
```

*NIS-Elements* can call functions from external DLL's. You should import the DLL, where the functions resides and then make a declaration of the functions. You should not import the following system DLL's: kernel32.dll, user32.dll, gdi32.dll, com32.dll, comdlg32.dll. This feature is available only for 32 bit version of *NIS-Elements*.

//Example

```
#import("luc_13.dll");
#import int RTF_ReplaceVariables(LPSTR destfile, LPSTR sourfile);
#import int RTF_FindQuestion(LPSTR sourfile, LPSTR question, long *length, \
LPSTR defvar);
#import int RTF_ReplaceQuestion(LPSTR destfile, LPSTR sourfile, LPSTR \
replacement);
```

**\_\_underC** (requires: *INTADV (Interactive Advanced Macro - API)*)

The **\_\_underC** directive assigns a function to be interpreted by the UnderC engine instead of the standard interpreter.

```
//Syntax
//works only on 32bit operating systems

__underC int inter_sharpen(int cols, int rows) { }
```

**#importUC** (requires: *INTADV (Interactive Advanced Macro - API)*)

The **#importUC** directive imports an API function to the UnderC engine so that it can be used from within there.

```
//Syntax
//works only on 32bit operating systems

#importUC DisplayCurrentPicture;
```

## 10.4.7 Operators

The following operators are supported by the system.

Caution

---

*Precedence of operators in NIS-Elements is different than might be expected. See 10.4.8 Priority of Operators (!).*

### **arithmetic operators**

+ Addition

- Subtraction

\* Multiplication

/ Division



% Modulo

### **assignment operators**

The assignment operator assigns the value of the right operand to the left operand.

= Addition

### **bitwise operators**

The bitwise operators compare each bit of the first operand to the corresponding bit of the second operand. The following bitwise operators are supported.

& Bitwise AND

| Bitwise OR

~ One's complement

### **pointer operators**

& Address of ?

\* Indirection

### **relational operators**

< Less than

<= Less than or equal to

> Greater than

>= Greater than or equal

== Equal

!= Not equal

### **logical operators**

The logical operators perform logical operations on expressions. The following logical operators are supported:

&& Logical AND

|| Logical OR

! Logical NOT

Caution

*Precedence of operators in NIS-Elements is different than might be expected. See 10.4.8 Priority of Operators (!).*

## 10.4.8 Priority of Operators (!)

If an expression contains more than one operator:

- Logical operators are evaluated *from right to left!!!*

- It is recommended to always use brackets to define different evaluation order:

*Example 10.1. Priority of logical operators in NIS-Elements*

```
if( !a && !b )
```

is evaluated by *NIS-Elements* as

```
!(a && !(b)) \
```

To make it work as expected it should be written as:

```
if( (!a) && (!b) )
```

- Priority of arithmetic operators is standard, division (/), multiplication (\*) and modulo (%) have higher priority than other operators.

## 10.4.9 C-like functions

The system can interpret your own C-like functions. An entry point to program is the `main()` function in your macro. If `main()` is not presented, for backward compatibility, the body of the macro is also considered as "main function". The new macros should use `main()` as an entry point.

The general C-like functions (also called interpreted functions, as opposed to basic system functions from *NIS-Elements*) have the following syntax.

Syntax

```
int MyFunction(int a, LPSTR str, double d)
{
int retval;
. . .
return retval;
}
```

**Return Value** The return value can be any basic data types (char, int, word, dword, int, double or pointer).

**Parameters** Parameters can be any basic data types (char int, word, dword, int, double or pointer).

Note

---

*Int and double parameters are automatically converted to string type, when they are assigned to text variable, or are one of the parameters of a function.*

Example

```
int main()
{
char buf[256];
my_function1(buf);
WaitText(0., buf);
return TRUE;
}
```

```

int my_function1(char *buf)
{
    strcpy(buf, "This function has a pointer to char array as a parameter");
    return TRUE;
}

```

## 10.5 Controlling Cameras by Macro



Every camera can be controlled from within a macro by changing its properties. There is a `CameraGet_*` and `CameraSet_*` macro function available for each camera property (e.g. `CameraGet_ExposureTime()`). The number of functions (beginning with `CameraSet_` and `CameraGet_`) and their actual names depend on the type of camera(s) you have currently connected to *NIS-Elements*.

**CameraGet\_** The *CameraGet\_...* functions retrieve current values of the properties determined by the function name postfix. So the `CameraGet_Exposure(int Mode, double *Exposure);` function retrieves the current value of Exposure of the specified Mode.

**CameraSet** The *CameraSet\_...* functions allows you to adjust camera properties. So the `CameraSet_Exposure(int Mode, double *Exposure);` function sets a new Exposure value for the specified Mode.

### How To Set Camera Properties

Instead of describing each property function, it is better to show you a universal way to handle a camera by macro:

- 1) Display the *View > Acquisition Controls > [Camera name] Settings*  control panel and find the property you would like to control.
- 2) Make a change of the property from within the control panel. For example, change the exposure time.
- 3) Display the *View > Macro Controls > Command History*  window and see the last function which has been called and its parameters (e.g. `CameraSet_ExposureTime(1, 500);`). This is the function you will control the exposure time with.
- 4) Experiment with the camera control panel in order to determine the right parameter values of the functions to be called within your macro.

## 10.6 Macro Preferences

This window enables you to configure key shortcuts to macros and set a macro to run automatically on startup.

Run the *Edit > Options*  command and switch to the *Macro* tab.

**Macros** Lists of macros that can be executed using hot keys or automatically at the beginning of the *NIS-Elements* session.

**Start Up** To launch a macro automatically after starting the *NIS-Elements* program, select it and press the `StartUp` button. StartUp field of the selected macro in the list box is filled.

---

**Hot Key** To assign a hot key to a macro, select the macro in the macro list box and press the **Hotkey** button. Select one of the predefined key combination and press **OK** button.

**Full Path** Shows the full pathname of the selected macro.

**Add** Adds a macro to the macro list box. Select Macro dialog box appears and you can search the disk for the macro.

**Remove** Removes a selected macro from the macro list box.

**Edit** Enables to edits a selected macro from the macro list box.

**Filename substitution** These lines display the current values of substitution strings available in macro commands.

**#** Letter representing a drive e.g. "C"

**##** A customizable path e.g. "C:\Images"

**###** Path to a directory according to the macro function used. It leads to the folder where *NIS-Elements* EXE file is placed and to a subdirectory - "Images" for functions concerning image files or "macros" for functions concerning macros. E.g.:

```
ImageOpen ("###\agnor.tif");           //leads to the IMAGES subdir  
RunMacro ("###\macro_001.mac");       //leads to the MACROS subdir
```

**####, #####** The current filename including/excluding the path. These substitutions can be used with the **Sequences** or **SequencesEx** macro commands.

**Information Shown in Caption** Optionally, the name of the current macro and an arbitrary user text can be displayed in the *NIS-Elements* caption.

**Defaults for this Page** Restores the default setting for Macro Options.

# 11. Movies

## 11.1 Capturing AVI Movie

1. Display the AVI Acquisition control panel by the *Acquire > AVI/MP4 Acquisition* command.
2. Adjust the advanced settings, especially define the file name and the destination folder.
3. Click the *Record* button. The live image appears and recording starts according to the settings.
4. Stop recording by the *Stop* button, or wait until the time defined in the *Duration* field passes.
5. When stopped, the AVI file remains opened in a new image window.

## 11.2 Save ND2 as AVI

You can easily create an AVI movie from an ND2 file. Run the *File > Save As* command and select the AVI format in the *Save As Type:* pull-down menu. The *Save As* window is extended by the following options:

**Use Default Video Compression** Either leave this box checked or select which video compression will be applied to the movie file. See 11.4 About Video Compression.

**XY Loop** When the ND document contains XY points, you can decide how these are saved.

**Current XY** Only the currently selected XY loop will be saved as a movie.

**All XY in separate files** A movie will be saved for each XY loop present in the ND2 file. An automatic number will be appended to the file name.

**Output Timing Parameters** Timing of frames of the resulting AVI movie may equal the Original Acquisition Time. If this does not suit your needs, define a custom interval in *ms*.

**Experiment Loop** Select the dimension, frames of which will be used to create the AVI movie.

**Use Time Loop** The T dimension frames of the current Z stack position will be used.

**Use Z-stack Loop** The Z dimension frames of the current Time position will be used.

**Use Z-stack and Time Loop** All frames of the ND2 file will be used in the AVI movie.

### Note



*If you have the QuickTime Player installed on your PC, a QuickTime movie (\*.mov) can be created as well.*

## 11.3 Creating a Movie from EDF Surface View

*(requires: EDF Module (Extended Depth of Focus))*

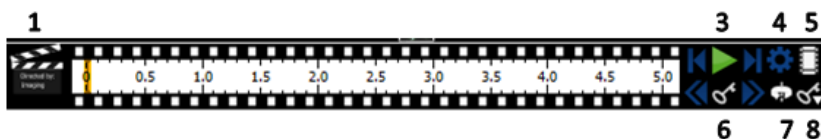
### 11.3.1 Creating a Movie Step by Step

1. Open a Z-Stack image and use the *Applications > EDF > Create Focused Image* .

2. Display the surface view by the *Applications > EDF > Show Surface View*  command.
3. Run the Movie Maker by the  button.
4. Adjust the movie settings by the *Settings* button a window appears where the movie duration and frame rate can be defined.
5. You can use one of the predefined presets or import some previously defined key frames by the *Export / Import* button.
6. Either enable or disable the Director's mode by the clapper button. See 11.3.3 Directors Mode.
7. Go to the beginning of the time line and define the initial zoom, position, rotation or component appearance of the object. This step is not necessary, but very much recommended. Press the *Add/Update key frame* button to fix the initial appearance.
8. Browse the time line, modify the object position each time, and set key frames by clicking the *Add/Update key frame* button. Press the *Play Movie* button for a preview.
9. When all the key frames are properly set, press the *Create Movie* button. The movie is created and the AVI file will open in a new image window.
- 10 Save the new file by the standard *File > Save As* command.

### 11.3.2 Movie Maker Controls

When you start the movie maker, the following bar appears at the bottom of the image view. The user is supposed to create a set of key-frames. The rest of the movie will be created by interpolation between these key frames. This gives the user absolute freedom in how the movie will look like.



1. The *Clapper* button. By this button, the director's mode can be turned on/off. See 11.3.3 Directors Mode.
2. Time line. By clicking to the time line you can browse the movie frames. When a key frame is added, a bold line appears in the time line to indicate its position.
3. *First Frame*, *Play*, *Last Frame* buttons.
4. *Settings* - by this button, you can define the movie length and the frame rate.
5. *Create Movie*
6. The key frame buttons. The existing key frames may be browsed by pressing the *Previous frame / Next Frame* buttons. The *Add / Update Key Frame* button inserts a key frame to the current position on the time line.
7. *Preset*s - a predefined rotation can be loaded. When you press the button and select the rotation type, the existing key frames settings will be overwritten.

8. *Import/Export Key Frames* - this enables you to save and load the key frame settings. A file with the \*.KEY extension will be created during the export.

### 11.3.3 Directors Mode

In the directors mode it is not needed to press the *Add/Update key frame* button every time the position of the object is changed. The director's mode detects all changes made to the object position and adds a key frame to the time line whenever such change is detected. A typical work-flow would look like this:

1. Turn ON the directors mode by the *Clapper* button. The time line gets highlighted with red color.
2. Click to the time line where a certain position of the object shall occur.
3. Move the object by mouse or adjust it in some way (zoom, crop, etc.) - at this moment, the key frame will be added to the time line automatically.
4. Click to the time line where another position shall occur, and so on...

The director's mode shall be used when you know precisely how the movie shall look.

## 11.4 About Video Compression

It is common to encode (compress) movies in order to save some disk capacity. This can be done using various types of codecs. A video codec is a software tool which can code (compress) and decode (decompress) video files. Remember that the codec you use to save the video will be required for playback as well (when playing the video on other computers).

Look at the list of codecs coming with *NIS-Elements* by default. If other codecs are installed on the computer, they become also available to *NIS-Elements*.

#### Caution

*32bit operating systems provide better choice of codecs because a 64bit codec must be used on a 64bit operating system, but there are not many of those in the meantime.*

#### *NIS-Elements* compression options

We performed a test by converting a 1 GB / 4000 frames ND2 file to AVI using different codecs. The results are displayed in the list:

**No compression** Original quality, file size: 787 MB.

**DV Video Encoder** This codec produces only videos with 640 x 480 resolution. Videos with other resolutions are stretched to fit, therefore we do not recommend to use this codec.

**MJPEG Compressor** Very good quality, file size: 240 MB.

**Cinepack Codec by Radius** Average quality, file size: 70 MB.

**Intel IYUV codec** Very good quality, file size: 390 MB. Provides the best quality / compression rate.

**Microsoft RLE** Very poor quality, file size: 90 MB.

**Microsoft Video 1** Very poor quality, file size: 6 MB.





---

# Index

## Symbols

3D, 140

[Camera name] Settings, 53, 173

## A

A Basic Way to Threshold an Image, 110

About ND Acquisition, 73

About Organizer, 101

About Video Compression, 177

Accepted Region, 132

AcqTime, 144

Acquisition and Analysis Modules, 11

Additional Module/Device Installation, 9

Additional Modules Available in NIS-Elements D, 11

Adjusting Program Preferences, 19

Advanced Layout, 25

Advanced options, 76

Aligning Objects, 160

Angle, 138, 144

Annotations and Measurements, 37, 130, 135, 136, 138, 141

Appearance Options, 36

Applying Restrictions, 126

Area, 138, 144, 146, 147, 149, 151

Arranging User Interface, 30

Arrays, 166

Assigning Colors to Channels, 89

Assigning Objectives to Nosepiece Positions, 45

Attaching Binary Layers to Channels, 115

Auto Capture Folder, 83

Auto Scale Settings, 98

Auto-Detect Tool, 121

Automated Measurement, 130, 142

Automated Measurement Results, 143

Automatic Image Sequence Recognition, 84

AVI/MP4 Acquisition, 175

AWB, 99

Axis A, 144

Axis A/B, 144

Axis B, 144

## B

Basic Workflow, 136

Basic Workflows, 41, 84, 123

Behavior, 37

Binary Editor, 113, 115

Binary Layer Color and Transparency, 116

Binary Layers, 113, 116

Binary Layers Control Panel, 113

Binary Operations, 114

Binary Toolbar, 116

BinaryArea, 144, 145, 153

BinaryAreaFraction, 144

Blue, 145

BoundsAbsBottom, 145

BoundsAbsLeft, 145

BoundsAbsRight, 145

BoundsAbsTop, 145

BoundsBottom, 145

BoundsLeft, 145

BoundsPxBottom, 145

BoundsPxLeft, 145

BoundsPxRight, 145

BoundsPxTop, 145

BoundsRight, 145

BoundsTop, 145

BrightVariation, 145

## C

C-like functions, 172

Calibrating an (Uncalibrated) Image, 123

Calibration, 123

Camera Control Panel, 38

Camera ROI, 68

Camera Settings, 53, 67

Capture, 67, 70, 130

Capture Multipoint, 76

Capture Timelapse, 36, 75

Capture Z-Series, 79

Capturing AVI Movie, 175

Cast Iron, 12

CentreX, 145

CentreXabs, 145

CentreXpx, 145

CentreY, 145

CentreYabs, 145

CentreYpx, 145

Change to Template, 161

Channel, 146

Channel Coloring, 88

Check Version Update, 22

Circularity, 146

Close, 85

Close All, 85

Close All but Current, 85

Closing Another Instance of NIS-Elements, 14

Closing Images, 85

Color Modes, 111

Command History, 163, 173

---

- Command Line Installation Options, 9
- Command Line Startup Options, 1
- Comment, 146
- Common ND Experiment Options, 73
- Connecting a Device to NIS-Elements, 49
- Connecting Cameras, 61
- Connecting Devices, 62
- Connectivity, 117
- Contrast, 34
- Control Bar, 90
- Controlling Cameras by Macro, 173
- Controlling Illumination Devices, 59
- Convex Hull, 146
- Convexity, 146
- Copying Channels by Drag and Drop, 89
- Count and Taxonomy, 142
- Create Custom LUTs, 100
- Create Focused Image, 141, 175
- Create Full View Snapshot (8bit RGB), 125
- Create Report, 162
- Creating a Movie from EDF Surface View, 175
- Creating a Movie Step by Step, 175
- Creating a Shared Layout, 14
- Creating Macros, 163
- Creating New Optical Configuration, 42
- Custom Metadata, 74

## D

- Data Export Options, 135
- DensityVariation, 146
- Device Manager, 41, 50, 52
- Device Modules, 12
- Device Updates, 10
- Diameter, 146
- Different Ways to Open an Image File, 83
- Directives, 168
- Directors Mode, 177
- Docking Panes, 26
- Drawing tools, 115
- Dual, Triple, Quad Camera Modes, 41
- Dynamic Data, 161

## E

- EDF Z-Profile, 29
- EdfRoughness, 146
- EdfRoughnessRA, 146
- EdfSurface, 146, 147
- Edit, 163, 164
- Elongation, 148
- Enable Mouse Joystick Z in Live, 53
- EqDiameter, 128, 148, 157
- Erasing Single objects, 116

- Export, 108
- Export Destinations, 134
- Exporting Results, 134

## F

- Field and ROI Features, 143
- FieldID, 148
- FillArea, 148, 149
- FillRatio, 149
- Filters and Shutters, 51
- Fixed Grabber Startup, 2
- Fixes, 11
- Floating Point Images, 86

## G

- General, 19
- Grain Size, 12
- Graph Memorizing, 108
- Graticules Properties, 124
- Green, 149
- Groups Tab Options, 17

## H

- Hardware licences, 21
- HASP Info, 10
- Height, 149
- Histogram, 107
- Histogram Options, 109
- Histogram Scaling, 108
- Homotopic Transformations, 117
- How Do I Display the Binary Layer?, 113
- How do I find out the content of my USB key?, 22
- How Do I Get a Binary Layer?, 113
- HueTypical, 149
- HueVariation, 149

## I

- Image Filter, 102
- Image Layers, 86
- Image processing, 109
- Image Properties, 20, 74
- Image Segmentation, 110
- Image Types, 87
- Image Window, 28
- Industrial Modules, 11
- Installation, 5
- Installation and Updates, 5
- Installing the Database Module on 64-bit Systems, 11
- Intensity Profile, 29
- IntensityVariation, 150
- Interaction with ROIs, 120

Interactive Advanced Macro (API), 165  
Introduction, 123  
Introduction to Binary Layers, 113  
Introduction to Image Acquisition, 67  
Introduction to Image Layers, 86  
Introduction to Look-Up Tables (LUTs), 96  
Introduction to ND Acquisition, 73  
Introduction to Optical Configurations, 42  
Introduction to ROIs, 120  
Is there an update for my software?, 22

## K

Keeps Z position and centers Piezo Z, 53

## L

Large Image (Covering) Multi-Point, 77  
Large Images, 94  
Layout Manager, 14, 31, 33, 35, 36  
Legacy Cameras and Devices, 65  
Length, 139, 150, 157  
Length xy, 150  
Length xyz, 150  
Lightpath Scheme, 63  
LineLength, 150  
Live - Fast, 67  
Live - Quality, 67  
Local and Global Variables, 166  
Locating XY Positions Between Images, 95  
LUTs - Non-destructive Image Enhancement, 96  
LUTs on Monochromatic Images, 99  
LUTs on Multichannel Images, 101  
LUTs on RGB Images, 97

## M

Macro Language Syntax, 166  
Macro Preferences, 173  
Main Menu, 25  
Main Window Components, 25  
Managing Binary Layers, 113  
Managing Objectives, 44  
Managing Optical Configurations, 43  
Manual Calibration, 46  
Manual Measurement, 136  
Mathematical Morphology Basics, 116  
Mathematical Morphology Examples, 118  
MaxFeret, 148, 150  
MaxFeret90, 150  
MaxIntensity, 151  
Mean\*ChannelName\*, 151  
MeanBlue, 151  
MeanBrightness, 151

MeanCa2+, 152  
MeanChord, 151  
MeanCorrFRET, 152  
MeanDensity, 151  
MeanFRETEff, 152  
MeanGreen, 151  
MeanIntensity, 151  
MeanRatio, 152  
MeanRed, 151  
MeanSaturation, 152  
MeanTitration, 152  
Measure Layers, 12  
MeasuredArea, 145, 152, 157  
Measurement Features, 144  
Measurement on Graph, 131  
Measurement Options, 130  
Measurement Tools, 138  
Methods Used for Stitching Large Images, 95  
Microscope Control Pad, 45, 51  
MinDistanceTo, 152  
MinFeret, 148, 152  
MinIntensity, 153  
Modifying Layouts, 31  
Modifying Menus, 35  
Modifying Toolbars, 33  
Monitor Calibration, 49  
Mouse Joystick and Auto Focus Z, 53  
Move Piezo Z to Home Position, 53  
Movie Maker Controls, 176  
Multi-point Acquisition, 76  
Multiple Binary Layers, 116

## N

Navigation in ND2 Files, 90  
ND Dimensions, 90  
ND Sequence Options, 74  
ND Views, 92  
ND.M, 153  
ND.T, 153  
ND2 Information, 93  
NearestObjDist, 153  
New Optical Configuration, 42  
NIS-Elements authentication, 14  
NIS-Elements Installation Steps, 5  
NIS-Elements Preferences, 19  
NumberObjects, 153  
NumberObjectsRestricted, 153  
Numbers, 153

## O

Obj ID, 153  
Object Count, 125, 128

- Object Count Procedure Example, 128
- Object Features, 143
- ObjectAreaFraction, 153
- Objective Calibration, 45
- Objective Calibration Using XY stage, 47
- Objective Clearance, 52
- Objectives, 44, 45, 123
- Objects Area, 153
- OC Panel, 59
- Open, 20, 83
- Open First, 83
- Open Last, 83
- Open Next, 19, 83
- Open Previous, 83
- Open, Save, Close Images, 83
- Opening Files in Progressive Mode, 94
- Opening Image Files, 83
- Operations with Binary Layers, 114
- Operations with Images within Organizer, 103
- Operations with Optical Configurations, 43
- Operators, 170
- Optical Configurations, 42, 43
- Options, 19, 36, 92, 129, 130, 143, 165, 173
- Organizer, 101
- Organizer Layout, 83, 101
- Orientation, 153
- Other, 13
- OuterPerimeter, 154
- Overrun Indication, 108

## **P**

- Perimeter, 146, 154, 157
- PerimeterContour, 154
- Pixel Classifier, 132, 133
- Point by Point (Manual) Multi-Point, 76
- Preparation, 107
- Preview Control Panel, 39
- Print, 38
- Priority of Operators (!), 171
- Privileges Tab Options, 18
- Processing On: Intensity/RGB/Channels, 109

## **R**

- Radius, 141, 154
- Random Multi-Point, 77
- Recalibrate Document, 123
- Recent Files, 83
- Record, 163
- Red, 154
- Region Settings, 132
- Regions of Interest - ROIs, 120

- Removing Dust from Camera Chip Using Shading Correction, 70
- Repetition issues, 118
- Report, 159, 161, 162
- Report Generator, 159
- Report Objects, 159
- Report Templates, 161
- Reporting, 105
- Resizing the Organizer Panes, 105
- Restrictions, 153
- Results, 127
- Roi ID, 154
- ROIArea, 153, 154
- Rough Measurement, 124
- Roughness, 155
- RoughnessInf, 155
- Run, 165
- Run From, 165
- Running a Macro, 165
- Running a Macro Upon Layout Change, 36

## **S**

- Sample Database Installation, 10
- Save, 84, 161
- Save As, 20, 84, 85, 163, 175, 176
- Save Control Panel, 38
- Save ND2 as AVI, 175
- Save Next, 19, 84
- Saving Image Files, 84
- Saving Images with UAC, 84
- Select [Camera Name], 1
- Select HW Configuration, 41
- Semixaxes, 141
- Shading Correction, 68
- Shading Correction Panel, 69, 70, 71
- Shading Correction per Optical Configuration, 70
- ShapeFactor, 155
- Show Lightpath Scheme, 64
- Show Physical Zoom, 49
- Show Surface View, 29, 176
- Show Z-Map Image, 29
- Simple Control, 12, 37
- Simple ROI Editor, 121
- Simple Shading Correction, 69
- Simplified User Interface, 37
- Slices View, 140, 157
- Software Copy Protection, 10
- Source, 155
- Source Data, 107
- Special Options, 75, 78
- Split Image, 94, 95
- Splitting Large Images, 94

---

Startup Hardware Configuration Selection, 41  
StartX, 155, 156  
StartXpx, 156  
StartY, 155  
StartYpx, 156  
Statements, 167  
Status Bar, 26  
StgPosX, 156  
StgPosY, 156  
Structures and Unions, 166  
Structuring Element = Kernel = Matrix, 118  
Sum\*ChannelName\*, 156  
SumBrightness, 156  
SumDensity, 156  
SumIntensity, 156  
Supported Image Formats, 85  
SurfVolumeRatio, 156  
Switching Between Loaded Images, 83

## T

The Installation DVD-ROM Content, 5  
Thresholding, 110, 113  
Thresholding the Image, 125  
Time, 157  
Time-lapse Acquisition, 75  
Tips and Tricks, 111  
To Change a Single Z Coordinate, 78  
Toolbar Buttons, 38  
Toolbars, 25  
Types of ROI, 120

## U

Units, 123  
Use Measurement Frame, 130, 143  
User Rights, 13  
User Rights Options, 15  
Users Tab Options, 16  
Using a Multiview Device, 41  
Using Auto Exposure while Switching between Optical Configurations, 44  
Using ROIs for analysis, 122

## V

Variable types, 166  
Velocity, 157  
Visual Device Manager, 59  
VolumeEqCylinder, 157  
VolumeEqSphere, 157

## W

Well Plate (Rectangular) Multi-Point, 77

What are Logical Devices?, 50  
What is a hardware licence?, 21  
What is SUA?, 22  
Width, 157, 158  
Working with the Measurement and Annotation Objects, 137

## X

XY Stages and Z Drives Tips, 52

## Z

Z-series Acquisition, 79  
Zoom, 37  
Zoom Configuration, 52





